

WTW Product Families...


WTW Product Families ...



Laboratory Meters **inoLab®** p. 6

720/730/735/740/750

- pH
- ISE
- D.O.
- Cond
- Multi

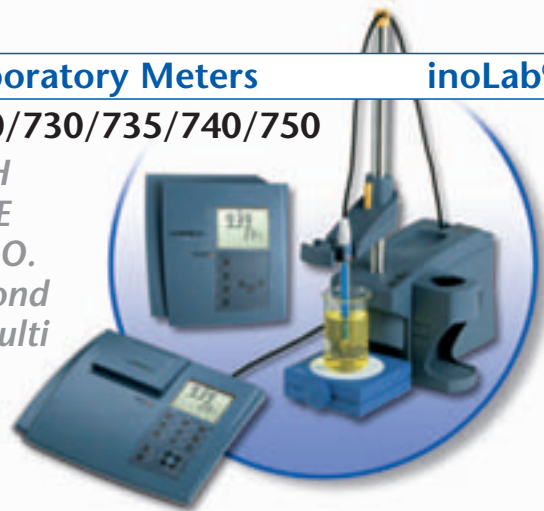

Field Meters p. 8

ProfiLine 197i

- pH
- D.O.
- Cond
- Multi



Handheld Meters p. 9

315i/330i/340i/350i

- pH
- ISE
- D.O.
- Cond
- Multi



VARIO p. 11

VARIO pH/C_{ond}

- pH
- Cond



Photometers p. 80

photoLab® S6/S12/Spektral
pHotoFlex

- Photometrical Tests

wide range of tests available



Turbidity Meters p. 104

Turb 430 / 355 / 550 / 555
NTU



Respirometer OxiTop® p. 62


OxiTop®/OxiTop® Control

- BOD
- Respiration

soil respiration, biodegradation
OECD, Biogas determination, respiration rate




Typical Applications



Environmental Monitoring
e.g. Photometric Measurements with pHotoFlex
see pp. 86




Swimming Pools
e.g. pH Control Measurement with pH 315i
see pp. 17




Pharmaceutical Water
e.g. Conductivity Flow Measurement with inoLab® Cond 730
see pp. 40




Food and Beverage Industry
e.g. Turbidity Measurement with Turb 430 + LabStation
see pp. 108




Chemical Water
e.g. pH/Conductivity/ISE Measurement with inoLab® pH/ION/Cond 750
see pp. 54




Cosmetics/Detergents
e.g. pH Measurement with VARIO pH
see pp. 18

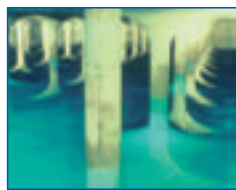


Groundwater
e.g. D.O./pH/Conductivity Measurement with Multi 197i
see pp. 56




Semi-Conductor Industry
e.g. pH/Conductivity Measurement with pH/Cond 340i
see pp. 60




Surface Water
e.g. D.O./pH/Conductivity Measurement with Multi 350i
see pp. 57




Process Technology
e.g. D.O./pH/Conductivity Measurement with Multi 340i
see pp. 60




Depths Profiles, Limnology
e.g. D.O./pH-/Conductivity Profiles with Multi 197i + Depths Armatures
see pp. 56




Laboratory Measurement
e.g. pH Routine Measurement with inoLab® pH 720
see pp. 14




Biotechnology (not autoclavable)
e.g. D.O./pH/Conductivity Measurement with Multi 350i
see pp. 57




Waste Water Treatment Plant: Aeration Basin
e.g. D.O. Control Measurement with Oxi 330i
see pp. 35




BOD Measurement
e.g. Dilution BOD with Analysis Program: inoLab® BSB/BOD 740; Self-Check Measurement with OxiTop®
see pp. 64



Biodegradation according to OECD
OECD 301: Determination with OxiTop® Control
see pp. 76



Fish Farming
e.g. D.O. Measurement with Oxi 315i and DurOx® incl. Protection Cap
see pp. 35



Wastewater: COD
e.g. COD Determination in Wastewater with photoLab® S12
see pp. 84

Contents

Contents

	page
New Instruments	2 - 5
General Description of Meters	6 - 11
Laboratory Meters inoLab®	6 - 7
Portable Meters ProfiLine	8
Handheld Meters	9 - 10
VARIO	11
Parameter	12 - 114
pH	12 - 21
Meters	14 - 18
Electrodes/Accessories	19 - 21
ORP	22 - 23
ISE	24 - 29
Oxygen (D.O.)	30 - 37
Meters	32 - 36
Sensors/Accessories	37
Conductivity	38 - 49
Meters	40 - 45
Conductivity Cells/Accessories	46 - 49
Multi-Parameter	50 - 61
BOD/Respiration	62 - 79
Photometry	80 - 103
Laboratory Photometers photoLab®	84 - 85
Portable Photometers pHotoFlex	86 - 89
Thermoreactors	90 - 91
Turbidity	104 - 109
Colony Counter	110 - 111
Software, Printers	112 - 114
General Information	115 - 128
Customer Service	115 - 117
On-line Instrumentation	118 - 119
WTW – About us	120 - 123
Index	127 - 128

Publisher



Wissenschaftlich-Technische Werkstätten GmbH

Dr.-Karl-Slevogt-Strasse 1

D-82362 Weilheim

Germany

Tel: +49 (0) 881 183-0
+49 (0) 881 183-100

Fax: +49 (0) 881 183-420

E-Mail: Info@WTW.com

Internet: <http://www.WTW.com>



New New Instruments

pHotoFlex Series and Turb 430

both units available in a convenient field case!

NEW

- A Complete Laboratory in the Field
- – Photometry
 - pH
 - Turbidity
- Convenient



These new complete SETs include everything you need for accurate field analysis – even an integrated tray for reagents and accessories.

Sets include:

- pH electrode SenTix® 41 (pHotoFlex models)
- 1 adjustable 5 ml pipette (pHotoFlex models)
- Calibration standards (pHotoFlex Turb and Turb 430 IR/T)
- Many useful accessories: empty cuvettes, buffer solutions pH 4,01 + 7,00, PC cable AK 540 B, cleaning tissues, stand, screw driver
- Storage space for accessories

In combination with the LabStation and its new software package, this is a complete solution for lab and field use as well.



New Instruments

The right test for every application!

Whether you require highly accurate tests with certificates of analysis or convenient tests for environmental surveys, WTW continues to offer an increasing number of test sets.

WTW now offers:

- Cuvette and reagent tests with barcode and certificate
- The cuvette tests for field use, without barcode and certificate
- The economical "powder pillow" test sets for field use and monitoring tasks

NEW



LabStation and LSdata

for the pHOtoFlex Series and Turb 430

Data management the smart way!



NEW

- Direct Excel Compatibility
- Conforms to GLP
- Complete Calibration Data

The new LSdata software package for all pHOtoFlex and Turb 430 models allows the user to process measured data collected in the field conveniently and according to GLP on your PC. The software is supplied with the LabStation (Order No. 251301). It makes your work considerably easier by providing:

- Data export from instrument to PC, password protected
- Excel processing, e.g. clear documentation of individual collection points
- Easy management of user defined methods and calibration data



New Instruments

Turb 430 T

Portable Turbidity Meter with Tungsten Lamp

NEW

- According to US EPA 180.1
- Scattered light characteristics according to Pharmacopoeia 5.0
- Suited for drinking water



Available in a convenient field case, too: With a tray, fittings for instrument and cuvettes, calibration set and many useful accessories.



Depending on the application, conditions, turbidity and user-defined standards: There is no measurement method with a wider range of uses than turbidity measurements!

Therefore we have developed the Turb 430T with tungsten lamp to complement the Turb 430IR with infrared light source. The T model is the ideal instrument for small particles in unstained solutions.

The convenient laboratory solution for up to 1100 NTU:

Turb 430 IR/T and LabStation

In combination with a LabStation the Turb 430 models are ideally suited for laboratory use:

The Turb 430 models have 1000 data point sets capability with data filter and interface. The data including calibration interval and protocol can be transferred to the PC according to GLP for further documentation via the LabStation's software.





New "Specialty Electrodes"

NEW

for acids, bases and everything in between!

WTW has broadened its selection of SenTix® pH sensors by 8 new sensors for special applications.

**Small sample size?
Measuring in narrow spaces?**

SenTix® Mic-D and SenTix® Mic-B will help you:

- Micro electrode
- Platinum diaphragm
- For smallest volumes

High concentrations or low ionic strength? What about temperature compensation?

No problem with SenTix® HWS:

- Electrode with SMEK head
- Ground joint diaphragm
- Integrated temperature sensor

Contaminated samples? Looking for a low-maintenance electrode?

The answer is SenTix® RJS:

- Polymer electrolyte
- Annular gap diaphragm
- Integrated temperature sensor

SenTix® L



pH measurement in large and in deep vessels?

Precise and with temperature sensor?


SenTix® L is the solution:

- Extra long electrode
- Platinum diaphragm
- Integrated temperature sensor

Difficult samples, need separate measuring and reference cells?

The glass pH cell SenTix® pH in combination with the reference electrode SenTix® R or the reference electrode SenTix® B with double electrode system will master these applications:

- Half cells
- Reference electrode with or without electrolyte bridge

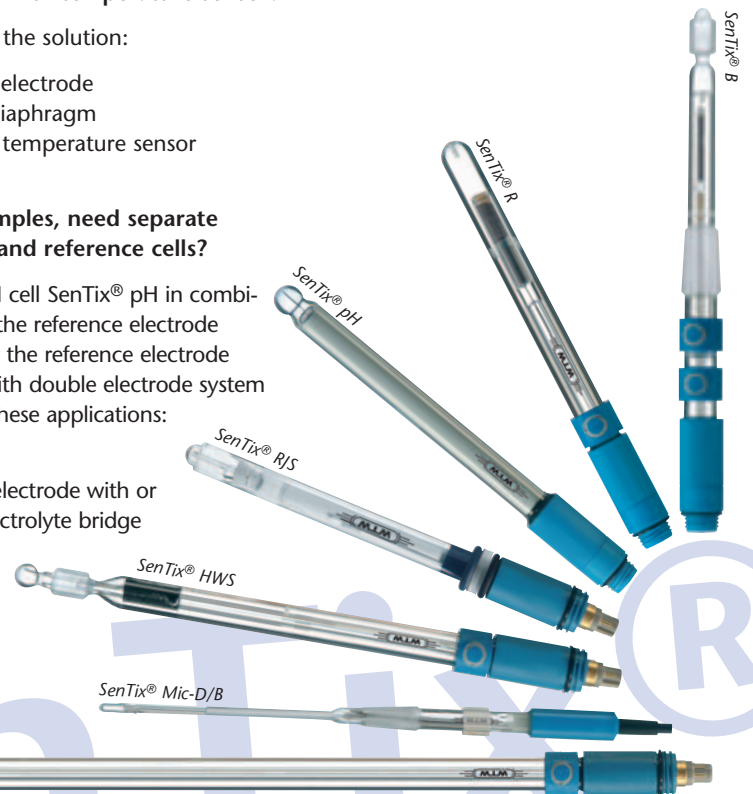

New ORP electrodes –

NEW

precious metals for relevant measurements!

Not just for Argentometry: **SenTix® Ag**

Alternative to Platinum: **SenTix® Au**

For use in contaminated samples: **SenTix® PtR**





inoLab[®]

inoLab[®] Laboratory Meter

inoLab[®] – innovative and progressive

The new inoLab[®] series: more powerful, more innovative. Modern design, new instrumentation. For reliable and precise laboratory measurements. With different function levels to meet all requirements concerning measuring safety, reliability, ease of use and flexibility.

inoLab[®] 720

- Large, bright display
- Easy-to-clean surface
- Simple Operation



Simple and reliable measurements. Ideal for measuring pH and ORP, oxygen, conductivity and temperature. Instrument for routine measurements at an excellent price.

Simple

All inoLab[®] function keys are well-laid-out. Automatic calibration and AutoRead functions ensure stable and reproducible results. The multi-function display for pH, oxygen, conductivity and temperature is large and easy to read.

Flexible

Battery or AC power for maximum mobility. Battery power means the instrument is immediately ready for use and is not influenced by interferences.

inoLab[®] 730

- Data transmission via bidirectional interface or optional built-in printer
- GLP/AQA
- Datalogger and memory for 800 entries



No matter whether pH, ORP, ISE, dissolved oxygen, conductivity, TDS or salinity: it is the perfect system for all standard measurements. Supports anyone who needs to perform documentation: AQS, GLP, according to ISO 9000.

Safe

The built-in printer (optional) with its high-quality thermal paper is guaranteed not to fade for 10 years. Functions such as a real-time clock, a serial number and a calibration protocol with sensor characteristics provide GLP-support. In addition to the bidirectional RS 232 interface the datalogger has memory for 800 data sets.

Comfortable

Modern and well-laid-out display for pH, oxygen, conductivity and temperature. The touch-sensitive keypad (with defined pressure point and fingertip guide recesses) is easy to clean and well-suited for biological laboratories.



inoLab® Laboratory Meter

inoLab® 740/750

Communicative

inoLab® 740 can be connected directly to a PC. The included "MultiLab® pilot" software package communicates with the instrument. Configuration, calibration, measurement: all functions are controlled by means of pull-down menus on a clearly arranged user interface. Similar to data base and export to PC standard formats. All measurement data are available as tables and graphics, both online or offline.

- Control via PC or terminal
- Comprehensive data acquisition
- Measurement functions for demanding laboratory analysis

Intelligent

The active multifunction box contains all measuring functions, numerous extension possibilities, connections, plug contacts and an electrically isolated interface for control via PC or terminal. The measuring procedure on the inoLab® 740 is initiated by pressing either the Run/Enter or AutoRead key. The built-in electronics process all incoming measuring signals immediately and make them available for communication with other peripheral devices.



inoLab® 750

Precise

With its outstanding technical features, the inoLab® 750 provides the utmost of measuring comfort for scientific demands. It offers easy operating for routine measuring, but its real capability shows up with the most difficult measuring tasks. A multitude of measuring combinations can easily be selected with a clearly structured menu. The instrument features password-protected operator levels, multi-level menu functions and complies with requirements of the pharmaceutical industry (CFR 21 Part 11, USP 28 etc.).





ProfiLine Portable Meters

ProfiLine 197i

- Robust, waterproof
- Precise, versatile
- For measurements at depths down to 100 m

Indestructible and waterproof, reliable and safe:

Both the housing and the technical specifications of **ProfiLine 197i** portable instruments are top class – suitable for rough field operation as well as for use in laboratories.

These portable field meters are characterized by their extreme robustness and shock-resistance. The instruments are hose-proof (IP 66) and submersible (IP 67). Special probes with cable lengths of up to 330 ft (100 m) are available for depth measurements.

The new built-in, environmentally-friendly NiMH rechargeable battery with an operation time of 600 hrs or mains supply, a carrying strap, carrying handle as well as the built-in electrode storage sleeve make the **ProfiLine 197i** a multi-talent for field or laboratory use.

ProfiLine 197i has excellent technical specifications, data transmission by means of the "MultiLab® pilot" software and can be ordered with or without probes.

No matter whether pH, dissolved oxygen, conductivity or multiparameter: there is a suitable ProfiLine available for every measuring task.



e.g. Measurement of D.O. depth profiles

Oxygen depth armature TA 197 Oxi with built-in temperature sensor, up to 100 m cable with waterproof plug (IP 67), pressure-resistant steel armoring and a screw-off protective hood, fits into 2" boreholes.

Use with battery-powered stirrer BR 325 for depth measurement.





Handheld Meters

Portable and Handheld Meters

The 315i, 330i and 340i series of handheld meters provide reliable measurements under difficult conditions both in the Lab and in the field. Lightweight and compact, the Series 3xxi meters are impact-resistant, hose-proof to IP 66 and meet the requirements for IP 67.

Model 315i

Simple

Simple operation means that measuring errors are avoided. All the function keys of the 315i can be operated even when wearing gloves; automatic calibration and AutoRead functions ensure stable and reproducible results. The multifunctional display for pH, oxygen, conductivity and temperature is easy to read.

Model 330i

Proven

The 330i, for field use with proven reliability, has a datalogger for 800 data records, multifunctional display, manual input for data, and 3000 hours of continuous operation.

Model 340i

All-rounder

In addition to all of the advanced features of all 3xxi series units, the model 340i can operate either with line or battery power. The model 340i also has a bi-directional interface for output to a PC or printer.



- For routine measurements
- Simple operation
- Reliable, Accurate

- Routine and standard measurements
- Wide Range of Sensors Available
- Datalogger

- Bidirectional interface RS 232
- Remote Control





The System

The key word is MobiLab

Our professional case can be used as a mobile laboratory. It contains a work platform with measuring beakers, a plug-in stand, all the necessary calibration and cleaning solutions as well as waterproof instructions – simply Ready to go.



Handheld Meter Multi 350i

- Simultaneous measurement of 3 parameters + temperature
- Back-lit graphic display
- Menu-driven

Flexible and robust. Compact precision without compromises. For simultaneous measurement of up to 4 parameters. Graphics display with user-friendly menu navigation as well as the highest level of accuracy and resolution make this instrument top class.

Exclusively for the Multi 350i – the new multiparameter Probes from WTW:

ConOx: combined conductivity and oxygen sensor for laboratory and field applications

MPP 350: outdoor multiparameter probe for pH, oxygen and conductivity

More on page 58.



VARIO

Handheld Meters and VARIO

General description of Meters

inoLab®

ProfiLine

Handheld Meters




VARIO

Simple Measurements at your Fingertip!

You notice it straight away: VARIO has no keys! Instead it has an innovative touch screen: all functions can easily be called up and set – with one hand. A light fingertip touch on the display is enough – and the VARIO switches on itself.

Placing the electrode in solution starts the measurement automatically. Even very small volumes can be measured with the optimized electrode.

The VARIO also functions as a laboratory clock, with stopwatch and timer.

-  Touch screen
-  Up to 1000 h continuous operation
-  Laboratory clock with timer function

Simply clever...



Handheld Instrument with added Value!


Due to their ruggedness and ergonomic form, the VARIO instruments are particularly suited for service applications. Therefore, VARIO is supplied with a comfortable carrying case assuring a trouble-free storage of the sensors.



pH Meters


pH Value

The water molecule has the property of dissociating into two ionic components in aqueous solutions.

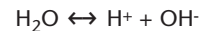

The H^+ ion is termed hydrogen ion or proton, the OH^- ion hydroxide ion.

The pH value describes the activity of hydrogen ions in aqueous solutions on a scale of -1 to 15. Based on this scale, liquids are characterized as being acidic, alkaline or neutral: a solution which is neither acidic or alkaline is neutral. This corresponds to a value of 7 on the scale. Acidity indicates a higher activity of hydrogen ions and a pH value lower than 7. Alkaline solutions are characterized by a lower hydrogen ion activity or higher hydroxide ion activity, respectively and a pH value above 7. The graph below uses examples to illustrate the pH scale.

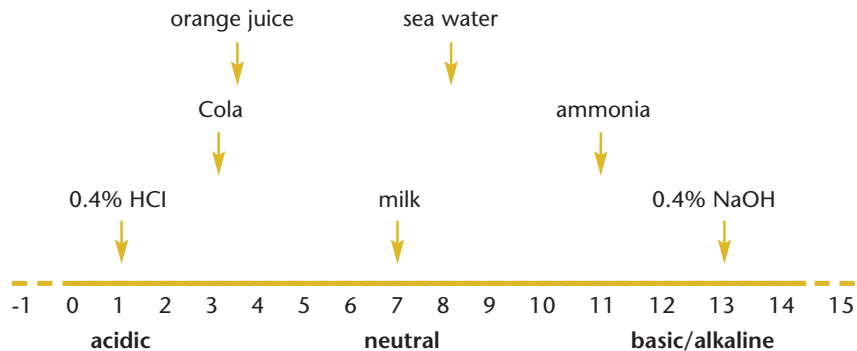

The pH scale is logarithmic. A difference of one pH unit represents a tenfold, or ten times increase or reduction of hydrogen ion activity in the solution. This explains how a solution's aggressiveness increases with the distance from the neutral point.

The pH value can be measured using electrochemical measuring systems, litmus paper, indicators and colorimeters. Of these methods, electrochemical sensors provide the most accurate results.

The pH electrode is an electrochemical sensor which consists of a measuring electrode and a reference electrode. The measuring electrode is made of special glass which, due to its surface properties, is particularly sensitive to hydrogen ions. It is filled with a buffer solution which has a pH value of 7. When placing the pH electrode into a test solution, the change in voltage is measured by the electrode by comparing the measured voltage to the stable reference electrode. This change is recorded by the meter and converted into the pH value displayed.



pH Meters

● recommended by WTW ○ conditionally applicable – not recommended

Application range	inoLab®					Profiline pH 197i	Handheld Meters				
	pH 720	pH 730	pH/ION 735	pH 740	pH/ION 740		VARIO pH	pH 315i	pH 330i	pH 340i	pH/ION 340i
Routine measurement	●	○	○	○	○	○	●	●	●	○	○
Routine measurement with documentation	-	●	●	●	●	●	-	-	-	●	●
AQA with documentation	-	●	●	●	●	●	-	-	-	●	●
R&D high resolution and precision	-	●	●	●	●	●	-	-	●	●	●
Control measurements	-	●	●	●	●	●	●	-	●	●	●
LIMS connection	-	●	●	●	●	●	-	-	-	○	○
Quality assurance	-	●	●	●	●	●	-	-	●	●	●
Training	●	●	●	●	○	○	●	●	●	○	○
Service	-	-	-	-	-	●	●	●	●	●	●
Laboratory measurements	●	●	●	●	●	●	●	-	-	○	○
Field measurements	-	-	-	-	-	●	-	●	●	●	●
Depth measurements	-	-	-	-	-	●	-	-	-	-	-
External control/PC connection/ PC control	-	●/○/-	-/○/-	●/○/●	●/○/●	●/○/-	-	-	-	●/○/-	●/○/-
pH/ION function	-	-	●	●	●	-	-	-	-	-	●
Ion-specific measurement programs	-	-	●	-	●	-	-	-	-	-	-

see page

for pH measurement with multi-parameter instruments see page 52





pH Meters

Laboratory pH Meters

Along with weighing and temperature measurements, pH is the most measured parameter in the laboratory. With inoLab® WTW offers a family of laboratory instruments which meet all measurement requirements from routine measurements to research applications.

inoLab® pH 720

simple and reliable

Easy to use routine laboratory pH/mV meter with large multifunctional display for pH and temperature, automatic temperature compensation, MultiCal® calibration system; for battery or line power operation.



inoLab® pH 730

compact and precise

Precision pH/mV meter with large multifunctional display for pH and temperature, automatic temperature compensation, MultiCal® calibration system, built-in measurement storage with GLP-conforming documentation and digital interface. Shown with optional built-in printer.



- Routine meter for precise measurement values (0.001 pH)

- Large display

- Easy-to-clean membrane keypad

- Supports all GLP needs

- Built-in printer (optional)

- Datalogger with memory for 800 data sets



Laboratory pH Meters

inoLab® pH 740

flexible and powerful

High-performance pH/mV/ION meter with graphic display and digital recorder function for pH, temperature and ion-selective measurement, automatic temperature compensation, high resolution (0.001 pH), MultiCal® calibration system, built-in measurement storage with GLP-conform documentation and digital interface. PC keyboard interface for connecting an external keyboard or a barcode reader. Includes software for direct control by PC. Built-in printer option available.

additional features

- 5-point calibration
- Selectable buffers
- Real-Time Graphic Display
- Built-in digital recorder
- Connection for bar-code reader or PC keyboard
- User Selectable Languages
- Multi-Level GLP Functions (password-protected operator levels)
- Free-of-charge downloads for MultiLab® pilot or terminal
- Firmware Updates

Computer-controlled precision meter

EMC-stabilized

Upgradeable firmware/software



IP 43



cETLus

3 Year Warranty

Technical Data

Model	pH 720	pH 730	pH 740
Range/Resolution	pH: -2.000 ... +19.999 pH; -2.00 ... +19.99 pH mV: -999.9 ... +999.9 mV; -1999 ... +1999 mV Temp.: -5.0 ... +105.0 °C (23 ... 221 °F)		-2.000 pH ... +20.000 pH, -2.00 ... +2.00 pH -999.9 ... +999.9 mV, -2000 ... +2000 mV -5.0 °C ... +105.0 °C
Accuracy (±1 digit)	pH: ±0.005 pH ±0.01 pH mV: ±0.3 mV, ±1 mV Temp.: ±0.1 K	±0.005 pH ±0.01 pH ±0.3 mV, ±1 mV ±0.1 K	±0.004 pH ±0.01 pH ±0.2 mV, ±1 mV ±0.1 K
Calibration	MultiCal® automatic calibration: AutoCal 2-/3-point AutoCal-Tec 2-/3-point ConCal® 1-/2-point ISECal -	2-/3-point 2-/3-point 1-/2-point -	2-/3-/4-/5-point 2-/3-/4-/5-point 1-/2-point 2- and 3-point

Ordering Information

□ with BNC plug ▲ with DIN plug

inoLab® Laboratory pH Meter SETs – with wide-range power supply 100-240 VAC (50/60 Hz) included	Order No.	Order No.
inoLab® pH 720 Simple and reliable pH meter, including SenTix® 42/41, without passive multifunction box and accessories	1A10-2117	1A10-1112
inoLab® pH 730 Compact precision pH meter with serial interface, including SenTix® 82/81, passive multifunction box and accessories	1A20-2119	1A20-1114
inoLab® pH 740P The intelligent pH measuring station, additionally equipped with built-in printer, incl. terminal, SenTix® 82/81 and accessories	1A31-2119	1A31-1114
Passive multifunction box (not included in pH 720 Set)	109 810	109 810
Other SETs or electrodes in SET see brochure "Product Details"		






pH Meters

Portable pH Meters

ProfiLine pH 197i

All WTW meters in the ProfiLine pH 197i series are both waterproof (IP 66) and submersible (IP 67). In addition, these units are able to float providing a high degree of comfort when used in field applications. With GLP memory functions, real-time clock, a display corresponding to the recorder output, 800 data records memory capacity, a carry handle, strap standard, the ProfiLine 197i is a complete pH measuring system. When used with the TA 197 Depth Armature, the ProfiLine 197i with its built-in preamplifier, is accurate to a depth of 330 ft (100m).

New: Powerful NiMH rechargeable batteries.

-  Robust, shockproof
-  Fully waterproof
-  Standard pH measurement and pH measurement down to depths of 330 ft (100 m)








Technical Data

Model	ProfiLine pH 197i
Range/Resolution	pH: -2.00 ... +19.99 pH, mV: -199.9 ... +199.9 mV; -1999 ... +1999 mV Temp.: -5.0 ... +105.0 °C (23 ... 221 °F)
Accuracy (± 1 digit)	pH: ±0.01 pH, mV: ±0.5 at +15 °C ... +35 °C (59 ... 95 °F), ±1 at +15 °C ... +35 °C (59 ... 95 °F) Temp.: ±0.1 K
Calibration	MultiCal® automatic calibration 1-, 2- and 3-point calibration, AutoCal, AutoCal-Tec and ConCal®

Ordering Information

Portable pH Meter – with wide-range power supply 100-240 VAC (50/60 Hz) included	Order No.
ProfiLine pH 197i	Robust, waterproof, submersible pH/mV meter 3A30-110
Depth armatures for measurements down to depths of 330 ft (100 m) see brochure "Product Details"	



pH Meters

Handheld pH Meters

pH Meters

pH 315i, pH 330i, pH 340i

WTW handheld pH meters are optimized for use on-site and field use, but can also be used in the laboratory. In particular the pH 340i, with its optional line power supply and serial interface, is suitable for applications in which precise measurements are required both in the laboratory and the field.

WTW handheld pH meters are available in three versions:

pH 315i:

Robust and waterproof battery-operated pH/mV meter. Measuring errors are avoided by the silicone keypad with only 5 keys and a simplified calibration method with automatic buffer recognition and display for standard buffers. AutoRead ensures stable and reproducible results.

pH 330i:

Robust and waterproof battery-operated pH/mV meter with built-in datalogger, real-time clock, GLP-supporting functions, display for calibration switchable between pH and mV. Configured with MultiCal[®] automatic calibration for buffer recognition, and automatic temperature compensation.

pH 340i:

Features of pH 330i, with additional analog and digital RS 232 output.




IP 66
IP 67

CE

cETLus

3 Year
Warranty

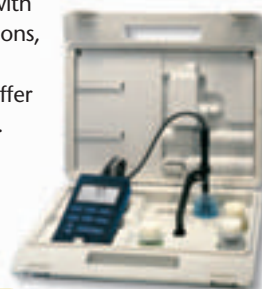

Robust

Waterproof (IP 67)

Large silicone keys

Technical Data

Model	pH 315i	pH 330i	pH 340i
Range/ Resolution	pH: -2.00 ... +16.00 pH mV: -1999 ... +1999 mV Temp.: -5.0 ... +105.0 °C (23 ... 221 °F)	-2.000 ... +19.999 pH; -2.00 ... +19.99 pH -999.9 ... +999.9 mV; -1999 ... +1999 mV -5.0 ... +105.0 °C (23 ... 221 °F)	
Accuracy (± 1 digit)	pH: ±0.01 pH mV: ±0.3 mV at +15 °C ... +35 °C (59 ... 95 °F) Temp.: ±0.1 K	±0.005 pH at +15 ... +35 °C (59 ... 95 °F) ±0.3 mV at +15 °C ... +35 °C (59 ... 95 °F) ±0.1 K	
Calibration	Simplified 1, 2 or 3-point calibration with automatic buffer recognition AutoCal automatic 3-point calibration with DIN buffers	MultiCal[®] automatic calibration: AutoCal automatic 1-, 2- or 3-point calibration with DIN buffers AutoCal-Tec automatic 1-, 2- or 3-point calibration with WTW technical buffers ConCal [®] conventional 2-point calibration with any buffers	

Ordering Information

Handheld pH Meter SETs	Order No.
pH 315i Robust and waterproof handheld pH meter, for battery operation, in mobile case set with SenTix [®] 41	2A10-1012
pH 330i Robust and waterproof handheld pH meter with datalogger, for battery operation, in mobile case set with SenTix [®] 41	2A20-1012
pH 340i Robust and waterproof handheld pH meter with datalogger and serial interface, in mobile case set with SenTix [®] 41	2A30-1012
Universal wide-range power supply 100 V - 240 V, 50-60 Hz; for 340i series	902 867
Other electrodes in SET see brochure "Product Details"	

Parameter

pH

ORP

ISE

Oxygen (D.O.)

Conductivity

Multi-parameter

BOD/Respiration

Photometers

Turbidity

Colony Counter

Software/Printers



NEW

- Compatible with most electrode types
- One-hand operation
- Twistable display

IP 65



3 Years Warranty



VARIO pH

You notice it immediately: apart from its ergonomic form the new VARIO has no keys, but has an innovative touch screen instead. In this way all functions can be called up and set – a simple touch is enough.

VARIO pH

Measuring in no time at all

Just a fingertip touch on the display – and VARIO is ready for use. Immersion in the solution starts the measurement automatically. The stable measurement can be read from the large display together with the temperature and can also be “frozen”. The memory has space for up to 50 measured values for later evaluation.

If the VARIO is not being used for pH measurements then it can be used as a laboratory clock or timer.

Light, handy, robust – it finds a place in every laboratory coat without dripping or leaving nasty stains, as it can be stored without KCl.

The VARIO can operate continuously for 1500 hours on one easily replaceable AA (1.5 V) battery. In addition, the VARIO can operate on line power.

VARIO comes standard with an intelligent glass electrode that is protected by a plastic casing. The conical protective cap does not require KCl which prevents the electrode from dripping and protects it from drying out.



The VARIO can do even more.

The adapter contained in the VARIO Set makes the instrument compatible with commercially available precision electrodes. The VARIO measures as accurately and reliably as a handheld instrument.

You can twist and turn as much as you want, the VARIO is an essential aid whenever speed is required in the laboratory or in production.

Technical Data

	VARIO pH
pH range	-2.00 ... 16.00
pH accuracy	±0.01 pH
Temperature range	-5.0 ... 100.0°C (23 ... 212 °F)
Automatic buffer recognition	TEC/NIST
Calibration points	3 (MultiCal®)

Ordering Information

VARIO	Order No.
VARIO Set	VARIO in the mobile case set, incl. short electrode with built-in temperature sensor and technical buffer 4 and 7
	2V00-001V
Other electrodes see brochure "Product Details"	



pH Meters

pH Electrodes and Accessories



SenTix® PLUS Electrodes	SenTix® 20 103 630			SenTix® 21 103 631			SenTix® 21-3 103 632			SenTix® 22 103 633			SenTix® 41 103 635			SenTix® 41-3 103 636			SenTix® 42 103 637			SenTix® 51 103 651			SenTix® 52 103 652			SenTix® 60 103 639			SenTix® 61 103 640			SenTix® 62 103 641			SenTix® 81 103 642			SenTix® 82 103 643			SenTix® 91 103 695			SenTix® 92 103 696								
	Measuring range pH	0 ... 14 pH									0 ... 14 pH									0 ... 14 pH									0 ... 14 pH									0 ... 14 pH									0 ... 14 pH							
Operating range °C (°F)	0...80 °C (32...176 °F)									0...80 °C (32...176 °F)									0...80 °C (32...176 °F)									0...100 °C (32...212 °F)									0...100 °C (32...212 °F)									0...100 °C (32...212 °F)								
Reference electrolyte	Gel									Gel									KCl 3 mol/l, Ag ⁺ -free									KCl 3 mol/l, Ag ⁺ -free									KCl 3 mol/l, Ag ⁺ -free									KCl 3 mol/l, Ag ⁺ -free								
Membrane shape	cylindrical									cylindrical									cylindrical									conical									conical									spherical								
Membrane resistance	<1 GΩ at 25 °C (77 °F)									<1 GΩ at 25 °C (77 °F)									<1 GΩ at 25 °C (77 °F)									<600 MΩ at 25 °C (77 °F)									<600 MΩ at 25 °C (77 °F)									<600 MΩ at 25 °C (77 °F)								
Diaphragm	Fiber									Fiber									ceramics									Platinum									Platinum									Platinum								
Shaft material	Noryl									Noryl									ceramics									Glass									Glass									Glass								
Shaft length (±0.08 in./±2 mm)	4.72 in. (120 mm)									4.72 in. (120 mm)									4.72 in. (120 mm)									4.72 in. (120 mm)									4.72 in. (120 mm)									6.79 in. (170 mm)								
Shaft Ø (±0.02 in./±0.5 mm)	0.47 in. (12 mm)									0.47 in. (12 mm)									0.47 in. (12 mm)									0.47 in. (12 mm)									0.47 in. (12 mm)									0.47 in. (12 mm)								
Temperature Sensor	-									built-in NTC (30 KΩ)									built-in NTC (30 KΩ)									-									built-in NTC (30 KΩ)									built-in NTC (30 KΩ)								
Connection	(1)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(1)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)	(2)																	
Electrode cable	(3)	(4)	(5)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(3)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)	(4)																	
Electrode plug	(6)/(7)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)/(7)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)	(6)																



SenTix® PLUS Electrodes	SenTix® L 103 655		SenTix® Mic-D 103 660		Mic-B 103 661	SenTix® HWS 103 662	SenTix® RJS 103 663	SenTix® pH 103 667	SenTix® R 103 668	SenTix® B 103 669
	Measuring range pH	0 ... 14 pH		0 ... 14 pH		0 ... 14 pH	0 ... 14 pH	2 ... 13 pH	0 ... 14 pH	-
Operating range °C (°F)	10...100 °C (50...212 °F)		-5...100 °C (23...212 °F)		-5...100 °C (23...212 °F)	-5...100 °C (23...212 °F)	0...80 °C (32...176 °F)	0...80 °C (32...176 °F)	-5...100 °C (23...212 °F)	-5...100 °C (23...212 °F)
Reference electrolyte	KCl 3 mol/l		KCl 3 mol/l		KCl 3 mol/l	KCl 3 mol/l	Referid®	-	KCl 3 mol/l	Double electrolyte system
Membrane shape	Spherical		Cylindrical		Spherical	Spherical	Calotte	Spherical	-	-
Membrane resistance	<600 MΩ at 25 °C (77 °F)		<1 GΩ at 25 °C (77 °F)		<600 MΩ at 25 °C (77 °F)	<600 MΩ at 25 °C (77 °F)	<600 MΩ at 25 °C (77 °F)	<600 MΩ at 25 °C (77 °F)	-	-
Diaphragm	Platinum		Platinum		Cut	Cut	Split ring	-	Platinum	Cut
Shaft material	Glass		Glass		Glass	Glass	Glass	Glass	Glass	Glass
Shaft length (±0.08 in./±2 mm)	16.73 in. (425 mm)		3.78 in. (96 mm)		6.69 in. (170 mm)	4.72 in. (120 mm)	4.72 in. (120 mm)	4.72 in. (120 mm)	4.72 in. (120 mm)	4.06 in. (103 mm)
Shaft-Ø (±0.02 in./±0.5 mm)	0.47 in. (12 mm)		0.12 in. (3 mm)		0.47 in. (12 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)
Temperature sensor	built-in NTC (30 KΩ)		-		built-in NTC (30 KΩ)	built-in NTC (30 KΩ)	built-in NTC (30 KΩ)	-	-	-
Connection	(9)		(6), (7)		(9)	(9)	(9)	(3)	(10)	(10)

(1): Plug head, (2): Fixed cable, (3): AS/DIN, AS/DIN-3 or AS/BNC, (4): Cable length 3 ft (1 m), (5): Cable length 9 ft (3 m), (6): DIN plug, (7): BNC plug, (8): Banana plug, (9) AS S/D1 or AS S/D3 or AS S/B1 or AS S/B3, (10) AS S/R



SenTix® PLUS Special Electrodes

	SenTix® H	SenTix® HW	SenTix® SP	SenTix® Sur	SenTix® Mic	SenTix® V	SenTix® FET-D	.../-B
Order No.	103 644	103 650	103 645	103 646	103 647	103 690	103 700	103 702
Measuring range pH	0 ... 14 pH	0 ... 14 pH	2 ... 13 pH	2 ... 13 pH	0 ... 14 pH	0 ... 14 pH	0 ... 14 pH	
Operating range °C (32 ... 176 °F)	0 ... 80 °C (32 ... 176 °F)	0 ... 60 °C (32 ... 140 °F)	0 ... 80 °C (32 ... 176 °F)	0 ... 50 °C (32 ... 122 °F)	0 ... 100 °C (32 ... 212 °F)	0 ... 80 °C (32 ... 176 °F)	0 ... 60 °C	
Reference electrolyte	KCl 3 mol/l, Ag ⁺ -free	KCl 3 mol/l, Ag ⁺ -free	Referid®	Referid®	KCl 3 mol/l, Ag ⁺ -free	Gel	KCl 3,3 mol/l, Ag ⁺ -free	
Membrane shape	Cylindrical	Cylindrical	Spear	Flat	Cylindrical	Flat	ISFET	
Membrane resistance	<2 GΩ at 25 °C (77 °F)	<800 MΩ at 25 °C (77 °F)	<400 MΩ at 25 °C (77 °F)	<1 GΩ at 25 °C (77 °F)	<700 MΩ at 25 °C (77 °F)	<500 MΩ at 25 °C (77 °F)		
Diaphragm	Cut	Cut	Hole	Split ring	Ceramic	Fiber	fritted polyethylene	
Shaft material	Glass	Glass	Glass	Glass	Glass	Noryl	ABS	
Shaft length (±0.08 in./±2 mm)	6.69 in. (170 mm)	6.69 in. (170 mm)	2.56/0.98 in. (65/25 mm)	4.72 in. (120 mm)	1.57/3.15 in. (40/80 mm)	1.22/0.79 in. (31/20 mm)	3.39 in. (86 mm)	
Shaft Ø (±0.02 in./±0.5 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)	0.59/0.02 in. (15/5 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)	0.67/0.75 in. (17/19 mm)	0.67 ... 0.51 in. (17 ... 13 mm)	
Connection	Plug head	Plug head	Plug head	Plug head	Plug head	—	DIN	BNC
Electrode cable* or AS/BNC	AS/DIN, AS/DIN-3, or AS/BNC	AS/DIN, AS/DIN-3, or AS/BNC	AS/DIN, AS/DIN-3, or AS/BNC	AS/DIN, AS/DIN-3, or AS/BNC	AS/DIN, AS/DIN-3, or AS/BNC	—	fixed cable (39.37 in./1 m)	
Electrode plug as selected	DIN plug or BNC, as selected	DIN plug or BNC, as selected	DIN plug or BNC, as selected	DIN plug or BNC, as selected	DIN plug or BNC, as selected	—		
Temperature sensor	—	—	—	—	—	NTC (30 KΩ)	NTC (30 KΩ)	

* not included

Calibration and Maintenance Supplies

The new buffer bottles from WTW

- easy to dispense
- easy to use
- reliable calibration



All WTW Technical buffers are certified accurate and are NIST/DIN traceable (see page 115, Services).

Applicable buffers	PL 4/7/9 DIN/NIST	APL 4/7/9 STAPL 4/7/9 DIN/NIST	TEP 4/7 Trace	TEP 10 Trace	TEP 10 Tec	TPL 4/7 Trace	TPL 10 Trace	TPL 10 Tec
inoLab® 7xx/197i/ Multi 350i	●	●	●	●	-	●	●	-
VARIO pH	●	●	●	●	-	●	●	-
pH 315i, 330i, 340i, pH/ION 340i	●	●	●	●	-	●	●	-
pH/Cond 340i, pH/Oxi 340i, Multi 340i	● not Multi 340i	● not Multi 340i	●	●	-	●	●	-
inoLab® Level 1,2,3/ pH 197	●	●	●	-	●	●	-	●
pH 330, 340, pH/ION 340	●	●	●	-	●	●	-	●
MultiLine P3/P4	-	-	●	-	●	●	-	●

Ordering informations for calibration and maintenance supplies see brochure "Product Details".



pH Electrodes and Accessories

Applications for SenTix® PLUS Electrodes

	SenTix® V	SenTix® 20 21-..., 22	SenTix® 41, 41-3, 42, RJS	SenTix® 51, 52	SenTix® 60, 61 62	SenTix® 81, 82	SenTix® 91, 92, L	SenTix® H	SenTix® HW, HWS	SenTix® Sp	SenTix® Sur	SenTix® Mic, MIC-D, MIC-B	SenTix® FET	SenTix® ORP**, PtR, Ag, Au Au, ORP*
Acids				○	●	●	●	○	○					
Ammonia				○	○	○	○	●						
Aquarium water	●	●	●	●	○	○	○							ORP, PtR*
Beer				●	●	●	●	○						
Beverages				●	●	●	●	○	○				○	
Bleach solution				○	○	○	○	●						
Boiler feedwater					○	○	○		●					
Bread										●				●
Cheese										●				●
Coffee extract				○	●	●	●	○	●					●
Condensate									●					
Cosmetics	○							○	●					●
Deminerlized water									●					
Developer				○	○	○	○	●	○					
Dispersion colors	○		RJS*					○	●					
Distilled water									●					
Drinking water	○	○	○	●	●	●	●	○	○					
Electroplating baths	○		RJS*	●	●	●	●	○	○					
Electroplating wastewater	●	●	●	○	○	○	○	○	○					○
Extracts				○	○	○	○	○	○	●				
Fixing baths			RJS*	○	○	○	○	●	●					ORP, PtR*
Fruit										●				●
Fruit juice	○			●	●	●	●	○	○					○
Groundwater		○	○	○	○	○	○	○	○					PtR*
Household cleansers	○	○	○	○	●	●	●	●	○					
Juice	○			●	●	●	●	○	○					○
Leather	○										●			
Lemonade				●	●	●	●	○	○					○
Lyes								●						
Margarine										●				●
Meat										●				○
Milk								○	●					○
Mineral water	○	○	○	●	●	●	●	○	○					○
Non-aqueous liquids				○	○	○	○		○					
Oil/water emulsions			RJS*					○	●					
Paint, water-soluble	○		RJS*					○	●					●
Paper	○										●			
Paper extract				○	●	●	●							
Protein-containing liquids					●	●	●	○	●			MIC-D/-B*		
Rainwater				○	○	○	○		●					
Saliva	●										●	○	●	
Salt solutions	○	○	○	○	●	●	●	○	○					
Sausage										●				●
Seawater				○	○	○	○	●	●					
Shampoo	○								●					●
Skin	○										●			
Soil extract	●	●	●	●	○	○	○	○						
Solids (penetration)										●				
Solids (surface)	○										●			
Sulfide-containing liquids			RJS*					○	●					PtR*
Surface water	○	○	○	●	●	●	●	○	○					
Suspensions			RJS*					●	●					
Swimming pool water	●	●	●	●	○	○	○							
Tapwater	○	○	○	●	●	●	●	○	○					
Tris buffer solutions							●		●					
Vegetable juice	○	○	○	●	●	●	●	○	○					○
Vegetables										●				●
Wastewater	○	●	●	●	○	○	○							PtR*
Wine				●	●	●	●		●					
Yogurt	○	○	○	○	●	●	●			●				●

● recommended by WTW ○ suitable for this application

* recommended for the mentioned type only

** for ORP Measurement see page 22

Parameter
pH
ORP
ISE
Oxygen (D.O.)
Conductivity
Multi-parameter
BOD/Respiration
Photometers
Turbidity
Colony Counter
Software/Printers



ORP measurements

ORP measurements

Reduction and oxidation are two central chemical terms that describe the ability of chemical agents to accept (reduction) or donate electrons (oxidation). In aqueous solutions, the ORP (Oxidation-Reduction Potential) voltage can be measured using a standard hydrogen electrode as reference. The reducing or oxidizing properties of a solution first are a matter of the reactants. By using an ORP electrode this change in potentials would be recorded as a positive or negative voltage.

ORP measurements monitor chemical reactions such as checking the denitrification of wastewater and disinfectant effect of detergents or the strength of plating baths.

Measurement of ORP voltage is carried out with ORP combination electrodes. Similar to pH electrodes, these consist

of a measuring electrode and a reference electrode. A metal electrode (normally a precious metal like gold, silver or platinum) is used in ORP combination electrodes in place of a glass membrane for carrying out the measuring function. The tendency for the chemical agents to accept or donate electrons determines the potential of the metal and thus the electrical potential of the combination electrode. ORP combination electrodes in use today contain a silver/silver chloride reference electrode, i. e. the indicated potential refers to this potential. Conversion to the standard hydrogen electrode system (U_H) and that of the silver/silver chloride reference electrode is easily possible.

$$U_H = U_{Meas} + U_{Ref}$$

ORP measurements can be carried out with all WTW pH/mV meters.

SenTix® ORP reference electrode potential against the standard hydrogen electrode

Temperature in °F (°C)	Potential in mV
32 (0)	+ 224
41 (5)	+ 221
50 (10)	+ 217
59 (15)	+ 214
68 (20)	+ 210
77 (25)	+ 207
86 (30)	+ 203
95 (35)	+ 200
104 (40)	+ 196
113 (45)	+ 192
122 (50)	+ 188
131 (55)	+ 184
140 (60)	+ 180
149 (65)	+ 176
158 (70)	+ 172



ORP measurements

NEW



SenTix® PLUS Electrodes	SenTix® ORP 103 648	SenTix® Ag 103 664	SenTix® Au 103 665	SenTix® PtR 103 666
Working range °C (°F)	0...100 °C (32...212 °F)	-5...100 °C (23...212 °F)	-5...100 °C (23...212 °F)	-5...100 °C (23...212 °F)
Reference electrolyt	KCl 3 mol/l	ELY/ORP/Ag	KCl 3 mol/l	Gel
Sensor	Platinum	Silver	Gold	Platinum
Sensor shape	Round 0.16 in. (4 mm)	Cylindrical cap	Cylindrical cap	Round 0.24 in. (6 mm)
Diaphragm	Ceramic	Ceramic	Ceramic	Split ring
Shaft material	Glass	Glass	Glass	Glass
Shaft length (±0,08 in./±2 mm)	4.72 in. (120 mm)	4.72 in. (120 mm)	4.72 in. (120 mm)	4.72 in. (120 mm)
Shaft-Ø (±0,02 in./±0,5 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)	0.47 in. (12 mm)
Temperature sensor	-	-	-	-
Connection	AS DIN/AS DIN-3, AS BNC			

Ordering Informations for accessories

Testing and maintenance supplies for ORP measurements	Order No.
SORT/RH Reagents for regenerating ORP electrodes consisting of activation powder (10 g) and clorina powder (30 g)	109 730
RH 28 ORP buffer solution 1 bottle of 250 ml: pH 7, U _H = 427 mV	109 740
ELY/ORP/AG Electrolyte with 2 mol/l KNO ₃ +0.001 mol/l KCl for combined ORP electrode with silver electrode	109 735

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.

Parameter

pH

ORP

ISE

Oxygen (D.O.)

Conductivity

Multi-parameter

BOD/Respiration

Photometers

Turbidity

Colony Counter

Software/Printers



Ionmeters

ISE

Ion-Selective Measurements

Determination of	Application
Lead (Pb ²⁺)	Soil samples
Bromide (Br ⁻)	Wine, plants
Cadmium (Cd ²⁺)	Soil samples
Calcium (Ca ²⁺)	Dairy products
Chloride (Cl ⁻)	Drinking water, foodstuffs
Cyanide (CN ⁻)	Electroplating baths
Fluoride (F ⁻)	Toothpaste, cement
Iodide (I ⁻)	Salt water
Potassium (K ⁺)	Wine, fertilizer
Copper (Cu ²⁺)	Electroplating baths
Sodium (Na ⁺)	Wine, boiler feed water
Nitrate (NO ³⁻)	Baby food, fertilizer, wastewater
Silver (Ag ⁺)	Electroplating baths
Sulfide (S ²⁻)	Proteins, sediments

Ion-selective measurement is a method for determining the concentration of dissolved ions. Potassium ions, sodium ions, fluoride or chloride are examples of such cations and anions which are directly measured in solutions. Indirect methods such as –titration allow the determination of aluminum, nickel ions, or sulfate.

Measurement with ISE's, like the measurement of pH, is a potentiometric method. ISE's are in two configurations:

1. separate ion-selective electrode and reference electrode
2. combined ion-selective electrode with built-in reference electrode

The ion-selective membrane of the electrode consists of a sparingly soluble salt of the ion to be measured (solid state electrodes), a PVC-membrane, modified by an ion exchanger or ion carrier (matrix electrodes), glass (glass electrode) or a gas-permeable plastic (gas-sensitive electrodes).

The activity of the ions to be measured determines the electrode current. With increasing activity of the anions the voltage turns more negative, with increasing activity of cations, more positive. A pH/ionmeter uses the electrode signal to calculate the concentration of the sample.

The wide range of possible applications includes, for example, the measurement of fluoride concentration according to DIN 38 405. Chloride content determination in concrete samples or nitrate concentration determination in fruit juices are further examples of the ways in which ion-selective measurement technology can be applied. An introduction to ion-selective measurement technology as well as application reports are available on our CD-ROM entitled "Principles of measurement technology".

	inoLab®					Handheld Meters
	pH/ION 735	pH/ION 740	pH 740, pH/Cond 740, Multi 740	pH/ION/Cond 750	pH/ION 340i, Multi 350i	
Application Range						
Occasional, simple ISE measurement	○	○	●	○	●	
Routine and standard measurement	●	●	○	●	○	
Advanced methods and procedures	●	●	–	●	–	
<i>see page</i>	26	26	15, 50	54	29, 57	



NEW

Ion-Selective Electrodes

WTW offers a complete range of ion selective electrodes and measuring instruments; you can choose between two types: **Half cells of the 500 series** which require a reference electrode from the R 503 type (an exception is the gas sensitive NH 500/2 which has a built-in reference electrode), or the **combination electrodes from the 800 series**.

These combination electrodes with built-in reference are excellent in handling and additionally offer the possibility to measure in small volumina. More over this they have an outstanding price performance ratio.



Type 800

Ion-Selective and Gas-sensitive Electrodes

Electrode type	Membrane®	Determinable ions	ISE Type 500 (Half cell, reference electrode necessary)	Reference electrode	combination ISE Type 800 (built-in reference electrode)	Measuring range	Bridge electrolyte	Ionic strength adjustment solution	Standard-solution (conc. 10 g/l)	pH range
Ammonium (NH ₄ ⁺)		Ammonium	NH 500/2	—	—	0,02...900 mg/l 10 ⁻⁶ ...5 x 10 ⁻² mol/l	—	MZ/NH ₃ /CN	ES/NH ₄	4-12
Bromide (Br ⁻)	S	Bromide	Br 500	↑ for all ion selective electrodes from the 500 series: R 503/P (0.08 in./ 2 mm pin plug) or R 503/D (0.16 in./ 4 mm banana plug) ↓ for all ion selective electrodes from the 500 series:	Br 800	0,4...79000 mg/l 5 x 10 ⁻⁶ ...1 mol/l	ELY/BR/503	ISA/FK	ES/Br	1-12
Cadmium (Cd ²⁺)	S	Cadmium	Cd 500		Cd 800	0,01...11000 mg/l 10 ⁻⁷ ...10 ⁻¹ mol/l	ELY/BR/503	ISA/FK	—	2-8
Calcium (Ca ²⁺)	L	Calcium, magnesium ^③	Ca 500 ^①		Ca 800 ^①	0,02...40000 mg/l 5 x 10 ⁻⁷ ...1 mol/l	ELY/BR/503	ISA/Ca	ES/Ca	2,5-11
Chloride (Cl ⁻)	S	Chloride	Cl 500		Cl 800	2...35000 mg/l 5 x 10 ⁻⁵ ...1 mol/l	ELY/BR/503	ISA/FK	ES/Cl	2-12
Copper (Cu ²⁺)	S	Copper, nickel ^③	Cu 500		Cu 800	0,0006...6400 mg/l 10 ⁻⁸ ...10 ⁻¹ mol/l	ELY/BR/503	ISA/FK	ES/Cu	2-6
Cyanide (CN ⁻) ^②	S	Cyanide	CN 500		CN 800	0,2...260 mg/l 8 x 10 ⁻⁶ ...10 ⁻² mol/l	ELY/BR/503	MZ/NH ₃ /CN	—	0-14
Fluoride (F ⁻)	S	Fluoride, aluminum phosphat ^③ , lithium ^③	F 500		F 800	0,02...gesätt. mg/l 10 ⁻⁶ ...gesätt mol/l	ELY/BR/503	TISAB	ES/F	5-7
Iodide (I ⁻)	S	Iodide, thiosulfate mercury	I 500		I 800	0,006...127000 mg/l 10 x 10 ⁻⁸ ...1 ⁻¹ mol/l	ELY/BR/503	ISA/FK	ES/I	0-14
Lead (Pb ²⁺)	S	Lead	Pb 500		Pb 800	0,2...20000 mg/l 10 ⁻⁶ ...10 ⁻¹ mol/l	ELY/BR/503	ISA/FK	ES/Pb	4-7
Nitrate (NO ₃ ⁻) ^②	L	Nitrate	NO 500 ^①		NO 800 ^①	0,4...62000 mg/l 7 x 10 ⁻⁶ ...1 mol/l	ELY/BR/503/N	TISAB/NO ₃	ES/NO ₃	2,5-11
Potassium (K ⁺) ^②	L	Potassium	K 500 ^①		K 800 ^①	0,04...39000 mg/l 10 ⁻⁶ ...1 mol/l	ELY/BR/503/K	ISA/K	ES/K	2-12
Sodium (Na ⁺) ^②	G	Sodium	DX 223 NA			0,05...23000 mg/l 2 x 10 ⁻⁶ ...1 mol/l	—	ISA/Na	ES/Na	>10
Silver (Ag ⁺) ^②	S	Silver	Ag/S 500		Ag/S 800	0,01...108000 mg/l 10 ⁻⁷ ...1 mol/l	ELY/BR/503	ISA/FK	—	2-12
Sulfide (S ²⁻) ^②	S	Sulfide	Ag/S 500		Ag/S 800	0,003...32000 mg/l 10 ⁻⁷ ...1 mol/l	ELY/BR/503	④	—	2-12

① Exchange measuring head

② S = solid state electrode, L = matrix electrode, G = glass electrode

③ Titration

④ Use according to operating instructions

⑤ Formulations for additionally required solutions are given in the application steps and operating instructions

Ordering Information for ISE electrodes and accessories see brochure "Product Details".

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.



Laboratory Ionmeters

inoLab® pH/ION 735



pH, mV and Concentration Measurements with a single Instrument

Whether routine measurements or demanding analysis: the pH/ION 735 is the ideal precision instrument for all uses. A graphical user interface makes high-resolution pH and ion measurement easy and comfortable. 5-point calibration for pH and up to 7 calibration points for ion measurements guarantees a high-precision measurement by calculating non-linear calibration curves. The Model pH/ION 735 has user defined method capability as well as preprogrammed incremental techniques.

For those who need to document their results: a datalogger with storage for 4,500 entries, bi-directional RS 232 interface, real-time clock, and GLP-supporting calibration protocols, as well as date, time and selectable sample identification number identify all data sets.

This instrument is also available with a built-in printer.

Incremental Methods

Advanced Incremental Methods

Menu-driven user interface

IP 43



CE TLus

3 Years Warranty

Advanced increment methods

Free software downloads

Comprehensive documentation options

inoLab® pH/ION 740

flexible and powerful

High-performance pH/mV/ion meter with graphic display and digital recorder function for pH, temperature and ion-selective measurement, automatic temperature compensation, high resolution (0.001 pH), MultiCal® calibration system, built-in measurement storage with GLP-conform documentation and digital interface. PC keyboard interface for connecting an external keyboard or barcode reader, and software for direct control by PC is included. With a built-in printer option available.

Features

- 5-point calibration by linear regression
- Selectable buffer sets
- Graphic evaluation possible
- Built-in digital recorder
- Connection for bar-code reader or PC keyboard
- User Selectable Languages
- Multi-Level GLP Functions (password-protected operator levels)
- Free-of-charge downloads for MultiLab® pilot or terminal
- Four to seven point ISE calibration with a modified nikolski algorithm
- Known addition, double-known addition and known subtraction
- Sample addition/subtraction
- Blank value addition



IP 43



CE TLus

3 Years Warranty



Laboratory Ionmeters

inoLab[®] pH/ION/Cond 750

Premium class from WTW:

Two galvanically isolated inputs allow independent measurements of pH-value, ORP or ion concentration. For accurate measurements along the characteristic curve of the electrode it is possible to carry out calibrations with up to seven standard solutions. The calculation of the calibration curve by using a modified Nikolski algorithm also takes the non-linear parts of the curve in to account. The following methods are used to evaluate the ionic concentration:

- Known addition/known subtraction
- Sample addition/sample subtraction
- Double known addition
- Blank value correction
- Known addition with blank value correction
- Reference measurement

A further highlight of this instrument is the possibility of conductivity measurement. Not only can specific resistance, salinity, and TDS be determined but also sample specific temperature coefficients. A wide range of additional functions like data administration, PC-operation by MultiLab[®] pilot, GLP-compliant calibration and data recording allow an easy integration into modern laboratory environment.

For further details see page 54.

- Two galvanically isolated pH/mV/ISE inputs
- Menu-operated with back-lit graphic display
- One pH and one ISE calibration record per each input





Technical Data inoLab® pH/ION 735 and 740

Model	pH/ION 735	pH/ION 740
Range/Resolution	pH: -2.000 ... +20.000 pH mV: -999.9 ... +999.9 mV -2000 ... +2000 mV Temperature: 23 ... 221 °F (-5 ... +105 °C/0.1 °C) Conc.: 0.000 ... 10.000 mg/l 0.00 ... 100.00 mg/l 0.0 ... 1000.0 mg/l 0 ... 2000 mg/l	-2.000 ... +20.000 pH -999.9 ... +999.9 mV - 2000 ... +2000 mV 23 ... 221 °F (-5 ... +105 °C/0.1 °C) Measuring range 1 (Resolution): 0.000 ... 9.999 (0.001) mg/l Measuring range 2: 0.00 ... 99.9 (0.01) mg/l Measuring range 3: 0.0 ... 999.9 (0.1) mg/l Measuring range 4: 0 ... 1999 mg/l
Accuracy (±1 digit)	±0.004 pH ±0.01 pH ±0.2 mV, ±1 mV ±0.1 K	±0.004 pH ±0.01 pH ±0.2 mV, ±1 mV ±0.1 K
Calibration	MultiCal® automatic calibration: AutoCal 2-/3-/4-/5-point AutoCal-Tec 2-/3-/4-/5-point ConCal® 1-/2-point ISECal 2- to 7-point Special functions: Known addition (single) Known subtraction, Sample addition, Sample subtraction, Blank value addition, Blank value correction	2-/3-/4-/5-point 2-/3-/4-/5-point 1-/2-point 2- to 7-point Special functions: Known addition (single and double)

Technical Data inoLab® pH/ION/Cond 750

Model	pH/ION/Cond 750	
Range/Resolution	pH: -2 ... 20.000 pH -2.00 ... 20.00 pH mV: -999.9 ... +999.9 mV -2000 ... +2000 mV Conc.: (mg/l) 0.000 ... 10.000 0.00 ... 100.00 0.0 ... 1000.0 0 ... 2000 Temperature: 23 ... 221 °F (-5 ... +105 °C)	
Accuracy (±1 digit)	pH: ±0.004 pH ±0.01 pH mV: ±0.2 mV, ±1 mV	
Temperature compensation	Automatic: 23 ... 221 °F (-5 ... +105 °C) 23.0 ... 212 °F (-5.0 ... 100 °C) Manual: -4 ... 266 °F (-20 ... +130 °C) NTC: 30 KOhm: ±0.1 Pt 1000: ±0.1 K	
Calibration	MultiCal® automatic calibration: AutoCal 2-/3-/4-/5-point AutoCal-Tec 2-/3-/4-/5-point ConCal® 1-/2-point ISECal 2- to 7-point Special functions: Known addition (single and double) Known subtraction, Sample addition, Sample subtraction, Blank value addition, Blank value correction	

Ordering Information

inoLab® Laboratory Ionmeters – with wide-range power supply 100-240 VAC (50/60 Hz) included		<input type="checkbox"/> with BNC plug	<input checked="" type="checkbox"/> with DIN plug
		<input type="checkbox"/> Order No.	<input checked="" type="checkbox"/> Order No.
pH/ION 735P	inoLab® pH/ION 735P with built-in printer for GLP-conform documentation	1G21-210	1G21-110
pH/ION 740P	inoLab® pH/ION 740P with built-in printer for GLP-conform documentation; extended measuring and storage options	1G31-210	1G31-110
pH/ION/Cond 750	Flexible and powerful precision bench-top pH/mV/Ion/conductivity-meter with two inputs, single instrument	1K30-210	1K30-110



Handheld Ionmeter ISE

Laboratory and Handheld Ionmeters

pH/ION 340i

pH, mV and Concentration Measurements in one Hand

The pH/mV and ionmeter pH/ION 340i offers the highest degree of flexibility possible. For pH measurements the instrument can be calibrated manually or automatically and offers simultaneous display of pH and temperature. For measurements with ion-selective electrodes the pH/ION 340i offers concentration display in mg/l. Direct display in mV to ± 999.9 mV in 0.1 mV steps; and to ± 1999 mV in 1 mV steps.

Of course, even in these higher ranges the concentration is calculated from a mV resolution of 0.1 mV. Calibration is carried out with up to three standards (selected from 16 standards in the range of 0.01 to 1000 mg/l).

Operating on either line power or rechargeable battery for up to 1500 hours with "LoBat" warning means that the instrument can be used in the Lab or in the field.

Lightweight and compact, these impact meters are IP 66 (hose proof) and IP 67 (submersible).

The built-in datalogger for up to 500 measurements together with GLP calibration protocol offers a comprehensive documentation of the results. With analog or digital data transfer (RS 232), automatic recognition of stable measurements (AutoRead), electrode evaluation and calibration interval monitoring functions ensure reproducible and comprehensible measurements.



Technical Data

Model	pH/ION 340i	
Range/Resolution	pH:	-2.000 ... +19.999 pH
	mV:	-999.9 ... +999.9 mV
		-1999 ... +1999 mV
Temperature:	23 ... 221 °F (-5 ... +105 °C/0.1 °C)	
Conc.:	0.01 ... 1999 mg/l	
Accuracy (±1 digit)	±0.005 pH ±0.01 pH ±0.3 mV, ±1 mV ±0.1 K	
Calibration	MultiCal® automatic calibration:	
AutoCal	2-point	
AutoCal-Tec	2-point	
ConCal®	1-/2-point	
ISECal	2-/3-point	

- Handy, waterproof
- Up to 1500 hours continuous operation
- GLP

Ordering Information

Handheld Ionmeter		Order No.
pH/ION 340i	Robust and waterproof handheld ionmeter with datalogger and serial interface	2G30-100
Universal wide-range power supply	100 V - 240 V, 50-60 Hz; for 340i series	902 867

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.

Parameter

pH

ORP

ISE

Oxygen (D.O.)

Conductivity

Multi-parameter

BOD/Respiration

Photometers

Turbidity

Colony Counter

Software/Printers



DO Meters

Oxygen Meters

Dissolved Oxygen

Some oxygen is dissolved in practically every liquid. For example, at a temperature of 68 °F (20 °C) and an atmospheric pressure of 1013 mbar saturated water contains about 9 mg/l oxygen. Ethanol can contain up to 40 mg/l whereas glycerol only has about 2 mg/l.

Each liquid takes up oxygen until the partial pressure of oxygen in the liquid is in equilibrium with the air or gas phase in contact with it. The actual oxygen concentration depends on a number of factors, such as temperature, air pressure, oxygen consumption by microorganisms in a biodegradation process or oxygen production by algae, etc.

The oxygen concentration is important for the:

- Living conditions for fish and microorganisms in waters
- Degradation processes in wastewater treatment
- Corrosion processes in pipelines
- Shelf life of beverages, etc.

The determination of the oxygen concentration was formerly carried out by the WINKLER titration method. Today electrochemical measurement is a recognized method in numerous standard procedures.

In its simplest form an oxygen sensor contains a working electrode and a counter-electrode. Both electrodes are located in an electrolyte system which is separated from the sample by a gas-permeable membrane. The working electrode reduces the oxygen molecules to hydroxide ions. In this electrochemical reaction a current flows from the counterelectrode to the working electrode. The more oxygen present in the sample, the larger the current signal. The oxygen meter calculates the oxygen concentration in the sample from this signal.



Oxygen Meters

● recommended by WTW ○ conditionally applicable – not recommended

Application Range	inoLab®			Profi-Line	Handheld Meters		
	Oxi 730	Oxi 740	BSB/BOD 740	Oxi 197i	Oxi 315i	Oxi 330i	Oxi 340i
Routine measurement	-	-	-	-	●	●	-
Routine measurement with documentation	●	●	●	●	-	-	●
AQA with documentation	●	●	●	●	-	-	●
R&D high precision	●	●	●	●	-	●	●
Control measurements	●	●	●	●	○	●	●
LIMS connection	●	●	●	●	-	-	○
Quality assurance	●	●	●	●	-	●	●
Training	●	●	●	○	●	●	○
Service	-	-	-	●	●	●	●
Laboratory measurements	●	●	●	●	-	-	○
Field measurements	-	-	-	●	●	●	●
Depth measurements	-	-	-	●	-	-	-
External control/PC connection/ PC control	●/●/-	●/●/●	●/●/●	●/●/-	-	-	●/●/-
BOD measurements with self-stirring sensor	●	●	●	●	-	-	○ with adaptor
BOD measurements with assessment program	-	-	●	-	-	-	-

see page

32

33

64

34

36

36

36

For oxygen measurement with multi-parameter instruments see page 50 ff

Application Range Sensors	ConOx	DurOx®	CellOx® 325	StirOx® G	TA 197 Oxi
BOD measurements	-	-	○	●	-
Fish farming	●	●	○	-	-
Surface waters	●	○	●	-	-
Ground water	○	-	○	-	●
Control measurements	●	●	●	○	-
Depth measurements	-	-	-	-	●
Laboratory measurements	○	-	●	○	-
Pharmaceutics	○	○	●	-	-
Biotechnology (non-autoclavable)	○	○	●	-	-
Wastewater treatment plant: aeration basin	○	●	○	-	-

applicable instruments:


Multi 350i

Oxi 3xxi
Multi 350i

all

inoLab®, 197i

197i

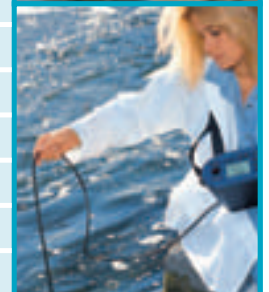

DO Meters

Laboratory Oxygen Meters

Laboratory Oxygen Meters

Oxygen is one of the most frequently measured parameters in the laboratory. It plays a large role in the degradation of substances and the growth of microorganisms, both in environmental technology and in biotechnology.

The inoLab® Oxi 740 instrument is particularly suitable for monitoring such processes. For routine measurements the inoLab® Oxi 730 is equipped with all the necessary functions which are required for documentation according to GLP.

inoLab® Oxi 730

compact and communicative

Standard laboratory oxygen meter with large multifunctional display for routine use in the laboratory. Automated functions such as air pressure correction, temperature compensation and OxiCal® rapid calibration make work easier. Datalogger, analog and digital output RS 232 as well as the optional built-in printer guarantee QA-conform documentation.

Features

- Simplified operator convenience
- Easy-to-clean touch-sensitive keypad
- Optional built-in printer



IP 43



cETLus

3 Year Warranty

- GLP documentation
- Comprehensive datalogging (800 data sets)
- Interface for self-stirring oxygen sensor StirrOx® G

The StirrOx® G can be operated by the inoLab® Oxi 730 for BOD₅ determination according to DIN EN 1899-1 resp. DIN EN 1899-2 with manual start/stop functions.



Laboratory Oxygen Meters

inoLab® Oxi 740

inoLab® Oxi 740 with Terminal or PC Software:
flexible and powerful

High-performance oxygen meter with graphic display and digital recorder function for oxygen measurement in the laboratory. Automated functions such as air pressure correction, temperature compensation by the IMT method and OxiCal® rapid calibration make work easier. The built-in measurement store, the digital output RS 232 and the optional built-in printer (paper width 4.4" resp. 112 mm) guarantee QA-conform documentation.

Features

- Built-in digital recorder
- Graphic evaluation possible
- Connection of bar-code reader or PC keyboard
- Automatic memory function if bar-code reader is used
- User Selectable Languages
- Multi-Level GLP Functions (password-protected operator levels)
- Limit input with acoustic alarm
- Free-of-charge software downloads for MultiLab® pilot or terminal

- Precision instrument
- Firmware/Softwareupdate
- External control (MultiLab® pilot)



IP 43



GETLUS

3 Year Warranty

Technical Data

Model	Oxi 730	Oxi 740
Range/ Resolution	O ₂ concentration: 0.00 ... 19.99 mg/l, 0.0 ... 90.0 mg/l* O ₂ saturation: 0.0 ... 199.9%, 0 ... 600%* O ₂ part. pressure: 0.0 ... 199.9 mbar, 0 ... 1250 mbar Temperature: 23 ... 122 °F (-5.0 ... +50.0 °C) * = depends on oxygen sensor and medium	O ₂ concentration: 0.00 ... 20.00 mg/l, 0.0 ... 90.0 mg/l* O ₂ saturation: 0.0 ... 200.0%, 0 ... 600%* O ₂ part. pressure: 0.0 ... 200.0 mbar, 0 ... 1250 mbar Temperature: 23 ... 122 °F (-5.0 ... +50.0 °C) * = depends on oxygen sensor and medium
Accuracy (± 1 digit)	O ₂ concentration: ±0.5% of value O ₂ saturation: ±0.5% of value Temperature: ±0.1 K	
Temperature compensation	32 ... 122 °F (0 ... +50 °C) automatic via IMT compensation	
Salinity correction	Automatic from 0.0 ... 70.0 adjustable via display	
Calibration	OxiCal® rapid calibration in OxiCal®-SL	

Ordering Information

inoLab® Laboratory Oxygen Meter SETs – with wide-range power supply 100-240 VAC (50/60 Hz) included		Order No.
inoLab® Oxi 730P	With built-in printer, including CellOx® 325, passive multifunction box and accessories	1B21-0111
inoLab® Oxi 740	Including terminal, PC software, CellOx® 325 and accessories	1B30-0111

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.

Parameter

pH

ORP

ISE

Oxygen
(D.O.)

Conductivity

Multi-
parameter

BOD/
Respiration

Photometers

Turbidity

Colony
Counter

Software/
Printers

DO Meters

Portable Oxygen Meters

ProfiLine Oxi 197i



The WTW oxygen meter of the ProfiLine Oxi 197i series is both hose-proof (IP 66) and submersible (IP 67). This easy to use meter conforms to GLP and has an 800 data point datalogger and accurate recorder output. External control via PC with MultiLab® pilot. They are equipped with a carrying/support handle and carrying strap as standard. The Oxi 197i is suitable for depth measurements down to 330 ft (100 m) in combination with the TA 197 Oxi depth armature.

New: Powerful NiMH rechargeable batteries.

- High-precision, indestructible, waterproof
- Recorder output corresponding to display
- Depths measurement down to 330 ft (100 m)



IP 66 IP 67 CE cETLus 3 Year Warranty

Oxygen depth armature **TA 197 Oxi** with built-in temperature sensor, up to 100 m cable with waterproof plug (IP 67), pressure-resistant steel armoring with screw-off protective hood, fits into small boreholes (2" dia.).

Battery-powered stirrer **BR 325** for profile and depth measurements.

Technical Data

Model	ProfiLine Oxi 197i
Range/Resolution	O ₂ conc.: 0.00 ... 19.99 mg/l (19.9 mg/l*), 0.0 ... 90.0 mg/l (90 mg/l*) O ₂ saturation: 0.0 ... 199.9% (199%*), 0 ... 600% <i>*depends on oxygen sensor and medium</i>
Accuracy (± 1 digit)	O ₂ conc.: ±0.5% of value O ₂ saturation: ±0.5% of value Temperature: ±0.1 K
Air pressure compen.	Automatic with built-in pressure sensor (500 ... 1100 hPa)
Temperature compen.	<2% at 32 ... 104 °F (0 ... +40 °C)
Salinity correction	Automatic from 0.0 ... 70.0, adjustable via display
Calibration	OxiCal® rapid calibration in OxiCal®-SL or OxiCal®-D

Ordering Information

Portable Oxygen Meters		Order No.
ProfiLine Oxi 197i	Robust, waterproof, submersible oxygen meter	3B30-010
Depth armatures for measurements down to depths of 330 ft (100 m) see brochure "Product Details"		



DO Meters

Handheld Oxygen Meters

Oxi 315i, Oxi 330i, Oxi 340i

Oxygen Meters



Oxi 340i with field armoring
FM 325 (optional)

Oxi 330i with protective armoring
SM 325 (optional)

- Robust, shockproof, waterproof
- Keypad with large, easy-to-operate silicone keys
- Extremely Long Battery Life



complete as SET



D.O. measurement with Oxi 315i and DurOx® incl. protection cap in fish farming

WTW handheld oxygen meters are suitable for checking both oxygen values in the field and for laboratory measurements. In combination with galvanic oxygen sensors DurOx® 325 or CelloX® 325 (no polarization time required) the measuring system is ready for immediate use anywhere.



DO Meters

Handheld Oxygen Meters

WTW handheld oxygen meters are available in three versions:

Oxi 315i

Simple, robust handheld meter with simplified keypad, without memory function and with no manual input facilities; in combination with DurOx® 325 with its protective hood it is particularly suitable for measurements in the fish-farming sector.

Oxi 330i

Proven robust battery-operated handheld meter with manual input functions, built-in datalogger (800 data records), GLP-supporting functions, automatic temperature and air pressure compensation and salinity correction...

Oxi 340i

As Oxi 330i, but with additional analog and digital RS 232 output and optional line adaptor. External control via MultiLab® pilot software.



Oxi 315i

Technical Data

Model		Oxi 315i	Oxi 330i	Oxi 340i
Range/ Resolution	O ₂ concentration:	0.00 ... 19.99 mg/l (19.9 mg/l*), 0.0 ... 90.0 mg/l (90 mg/l*)		
	O ₂ saturation:	0.0 ... 199.9% (199%*), 0 ... 600%		
	O ₂ part. pressure:		0.0 ... 199.9 mbar (199 mbar*), 0 ... 1250 mbar	
	Temperature:	23 ... 122 °F (-5.0 ... +50.0 °C)		23 ... 122 °F (-5.0 ... +50.0 °C)
		<i>*if DurOx® oxygen sensor is used</i>		
Accuracy (±1 digit)	O ₂ concentration:	±0.5 % of value		
	O ₂ saturation:	±0.5 % of value		
	O ₂ part. pressure:		±0.5 % of value	
	Temperature:	±0.1 K		
Temperature compensation		2% at 32 ... 104 (0 ... +40 °C)		
Air pressure compensation		Automatic with built-in pressure sensor (500 ... 1100 hPa)		
Salinity correction		0 or 30 solid SAL	Automatic from 0.0 ... 70.0, adjustable via display	
Calibration		OxiCal® rapid calibration in OxiCal®-SL or OxiCal®-D		

Ordering Information

Handheld Oxygen Meter SETs	Order No.
Oxi 315i Handheld oxygen meter, including DurOx® 325-3, professional case and accessories	2B10-0017
Oxi 330i Handheld oxygen meter, including CellOx® 325, professional case and accessories	2B20-0011
Oxi 340i Handheld oxygen meter, including CellOx® 325, professional case and accessories	2B30-0011
Universal wide-range power supply 100 V - 240 V 50-60 Hz; for 340i series	902 867
Other sensors in SET see brochure "Product Details"	

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.



DO Meters

Oxygen Sensors and Accessories

Oxygen Meters



CelloX® 325

Membrane covered galvanic oxygen sensor

- Temperature compensation
- Up to 6 months usage of membrane
- High signal resolution
- Rapid response
- Waterproof sensor (IP 68 – 2 bar)
- Incl. calibration vessel OxiCal®-SL
- Membrane leakage monitoring

Galvanic sensors: immediately ready for measurement

Equipped with protective hood for rough conditions

Simple air calibration using calibration vessel



DurOx® 325

only for handheld and field meters Oxi 3xxi/Oxi 197i and Multi 350i

Membrane covered galvanic oxygen sensor

- Temperature compensation
- Long working life – up to 6 months with one electrolyte filling
- Low approach flow
- Waterproof sensor (IP 68 – 2 bar)
- Incl. calibration vessel OxiCal®-D
- With protective hood SK-D as standard

StirrOx® G

for inoLab® Oxi 730, inoLab® Oxi 740 and ProfilLine Oxi 197i

Self-stirring oxygen sensor – simultaneous stirring and measurement

- Single-handed operation for rapid series measurements
- Constant approach flow for high reproducibility
- Extremely low self-consumption of oxygen – only 0.008 µg h⁻¹ (mg/l)⁻¹
- Incl. calibration vessel OxiCal®-ST
- Temperature compensation
- Membrane leakage monitoring

Calibration and storage vessels

Various calibration and storage vessels are available for oxygen sensors – see brochure "Product Details".

Other sensors and accessories see brochure "Product Details".

Ordering Information

Oxygen sensors	(The sensor prices include the corresponding accessory case with spare parts and maintenance supplies)	Order No.
StirrOx® G	Self-stirring oxygen sensor for oxygen determination in Karlsruhe bottles and Winkler bottles, with OxiCal®-ST calibration and storage vessel	201 425
CelloX® 325	Galvanic oxygen sensor with OxiCal®-SL calibration and storage vessel, waterproof plug, , cable length 1.5 m	201 533
DurOx® 325-3	Galvanic oxygen sensor with OxiCal®-D calibration vessel, water proof plug, cable length 3 m	201 570




Conductivity Meters

Conductivity

Conductivity is a parameter used to measure the ionic concentration and activity of a solution. The more salt, acid or alkali in a solution, the greater its conductivity. The unit of conductivity is S/m , often also S/cm . The scale for aqueous solutions begins with pure water at a conductivity of $0.05 \mu S/cm$ ($77^\circ F/25^\circ C$). Naturally occurring waters such as drinking water or surface water have a conductivity in the range $100 - 1000 \mu S/cm$. At the upper end of the chart some acids and alkalines can be found.

Conductivity measurements are used for applications such as in the production of ultrapure water or determining the salinity of sea water.

Conductivity is measured by making a measurement of the electrical resistance. The simplest kind of measuring cell used consists of two similar electrodes. An alternating voltage applied to one of the electrodes causes the ions in the solution to migrate towards the electrodes. The more ions in the solution, the greater the current which flows between the electrodes. The instrument measures the current and uses Ohm's law to calculate first the conductance of the solution and then – by taken the cell data into account – the conductivity.

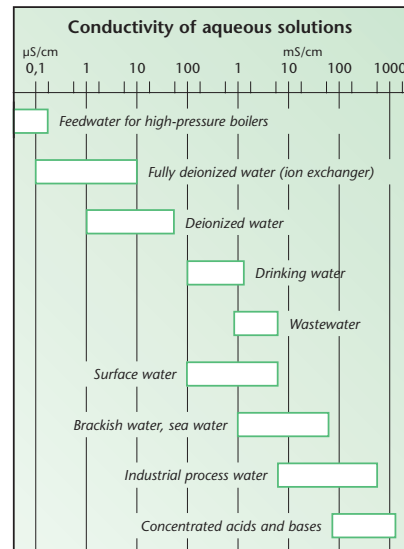



● recommended by WTW ○ conditionally applicable – not recommended

Application Range	inoLab®			Profi-Line	VARIO	Handheld Meters		
	Cond 720	Cond 730	Cond 740	Cond 197i	Cond	Cond 315i	Cond 330i	Cond 340i
Routine measurement	●	-	-	-	●	●	●	-
Routine measurement with documentation	-	●	●	●	-	-	-	●
AQA with documentation	-	●	●	●	-	-	-	●
R&D high precision	-	●	●	●	-	-	●	●
Control measurements	-	●	●	●	●	-	●	●
LIMS connection	-	●	●	●	-	-	-	○
Quality assurance	-	●	●	●	-	-	●	●
Training	●	●	●	○	●	●	●	○
Service	-	-	-	●	●	●	●	●
Laboratory measurements	●	●	●	●	●	-	-	○
Field measurements	-	-	-	●	-	●	●	●
Depth measurements	-	-	-	●	-	-	-	-
External control/PC connection/ PC control	-	●/●/-	●/●/●	●/●/-	-	-	-	●/●/-
Salinity/TDS measurement	●	●	●	●	●	only SAL	●	●
Specific resistance	●	●	●	-	-	●	●	●
Suitable for USP 28	●	●	●	●	-	-	●	●
Measurement of ultrapure water	●	●	●	●	●	●	●	●
Trace conductivity	●	●	●	●	-	-	●	●

see page 40 40 41 42 45 44 44 44
for conductivity measurements with multi-parameter instruments see page 50

Application Range Sensors	TetraCon® 325	TetraCon® 325/S	LR 325/01	LR325/001	TetraCon® DU/T	TA 197 LF
USP 28	-	-	●	●	-	-
Pharmaceutical water	○	-	●	●	-	-
Chemical water	○	-	-	-	●	-
Ground water	●	-	○	-	-	●
Surface waters	●	-	-	-	-	-
Depth measurements (barrages)	○	-	-	-	-	●
Laboratory measurements	●	-	●	●	-	-
Foods industry (juices)	●	-	-	-	○	-
Swimming pools	●	-	-	-	○	-
Pharmaceutics	●	-	●	○	○	-
Cosmetics/Detergents	○	●	-	-	-	-
Semi-conductor industry	-	-	●	●	-	-
Paint/Varnish (soluble)	●	○	-	-	-	-
Electroplating	●	-	-	-	-	-

applicable instruments: all/except VARIO 40 all/except VARIO/315i 40 all/except VARIO 41 all/except VARIO/315i 42 all/except VARIO/315i 45 Cond 197i 44

Parameter
pH
ORP
ISE
Oxygen (D.O.)
Conductivity
Multi-parameter
BOD/Respiration
Photometers
Turbidity
Colony Counter
Software/Printers



Conductivity Laboratory Conductivity Meters

Conductivity is an important parameter that is often used for monitoring water quality. In the laboratory sector this parameter has also increased in importance since the introduction of USP 28 Guideline for pharmaceutical water. WTW inoLab® Cond laboratory instruments meet all the requirements for measurements according to this standard.

- Application specific displays
- Touch-sensitive keypad with pressure point and tactile response
- Battery or line power operation

inoLab® Cond 720

Simple and reliable

Routine laboratory conductivity meter with large multifunctional display, parallel temperature display and automatic temperature compensation. 68 °F or 77 °F (20 °C or 25 °C) can be set as reference temperature. Both TDS and salinity measurements are possible as well as conductivity and temperature. The ability to set different cell constants means that other special conductivity cells can be connected as well as the TetraCon® 325 4-electrode conductivity cell and LR 325/01 ultrapure water conductivity cell.



- Simplified operator convenience
- GLP documentation via PC or optional built-in printer
- Meets all the requirements of USP 28

inoLab® Cond 730

Compact and precise

Standard laboratory conductivity meter with large multifunctional display, parallel temperature display and automatic temperature compensation. The datalogging capability and the real-time clock allows for all GLP functions. Data output can take place via the optional built-in printer on thermal paper accepted for use for official documents or via the built-in RS 232 digital interface via a PC or external printer.

68 °F or 77 °F (20 °C or 25 °C) can be set as reference temperature. Both TDS and salinity measurements are possible as well as conductivity and temperature. The ability to set different cell constants means that other special conductivity cells can be connected as well as the TetraCon® 325 4-electrode conductivity cell and LR 325/01 ultrapure water conductivity cell.



IP 43



cETLus

3 Year
Warranty



inoLab® Cond 740

**inoLab® Cond 740 with Terminal or PC software:
flexible and powerful**

High-performance laboratory conductivity meter with graphic display and digital recorder function, parallel display of temperature and automatic temperature compensation. A built-in datalogger and a real-time clock allows for all QA conforming functions. The optional built-in printer allows data printout on thermal paper accepted for use in official documents.

68 °F or 77 °F (20 °C or 25 °C) can be set as reference temperature. Both TDS and salinity measurements are possible as well as conductivity and temperature. The ability to set different cell constants means that other special conductivity cells can be connected as well as the TetraCon® 325 4-electrode conductivity cell and LR 325/01 ultrapure water conductivity cell.

A PC keyboard interface allows an external keyboard or a barcode reader to be connected.

additional features

- Built-in digital recorder
- Real-Time Graphic Display
- User Selectable Languages
- Multi-Level GLP Functions (password-protected operator levels)
- Limit input with acoustic alarm
- Meets all the requirements of USP 28
- Free-of-charge software downloads for MultiLab® pilot or terminal

IP 43



GETLUS

3 Year Warranty



Conductivity Meters

- Meets all the requirements of USP 28
- TDS and salinity measurement
- External control with MultiLab® pilot via PC

Technical Data

Model	Cond 720 and Cond 730	Cond 740
Range/Resolution	Conductivity: 0.0 µS/cm ... 500 mS/cm in 5 measuring ranges or AutoRange additionally for K = 0.1 cm ⁻¹ : 0.00 µS/cm ... 19.99 µS/cm K = 0.01 cm ⁻¹ : 0.000 µS/cm ... 1.999 µS/cm Temperature: 23 ... 211.82 °F (-5.0 ... +99.9 °C) Salinity: 0.0 ... 70.0 TDS: 0 ... 1999 mg/l Resistivity: 0.000 ... 1999 MΩcm	0.0 µS/cm ... 2000 mS/cm in 5 measuring ranges or AutoRange additionally for K = 0.1 cm ⁻¹ : 0.00 µS/cm ... 20.00 µS/cm K = 0.01 cm ⁻¹ : 0.000 µS/cm ... 2.000 µS/cm 23 ... 221 °F (-5.0 ... +105.0 °C) 0.0 ... 70.0 0 ... 2000 mg/l 0.000 ... 2000 MΩcm
Accuracy (± 1 digit)	Conductivity: ± 0.5 % of value Temperature: ± 0.1 K	
Reference temperature	68 ... 77 °F (20 °C or 25 °C) selectable	
Cell constants	0.475 cm ⁻¹ , 0.1 cm ⁻¹ and 0.01 cm ⁻¹ fixed; freely selectable 0.25 ... 2.5 cm ⁻¹ and 0.09 ... 0.11 cm ⁻¹	0.475 cm ⁻¹ , 0.1 cm ⁻¹ and 0.01 cm ⁻¹ fixed; freely selectable 0.25 ... 2.5 cm ⁻¹
Temperature compensation	Automatic or switched off	
Temperature coefficient	<ul style="list-style-type: none"> • Non-linear function for natural water to EN 27 888 • Linear compensation from 0.001 ... 2.999%/K • No compensation 	
Calibration	With 0.01 mol KCl	

Ordering Information

inoLab® Laboratory Conductivity Meter SETs – with wide-range power supply 100-240 VAC (50/60 Hz) included	Order No.
inoLab® Cond 720 Simple and reliable conductivity meter, including TetraCon® 325 and accessories	1C10-0111
inoLab® Cond 730 Compact precision conductivity meter, including TetraCon® 325, passive multifunction box and accessories	1C20-0111
inoLab® Cond 740P The intelligent conductivity measuring station with additional built-in printer, TetraCon® 325 and accessories	1C31-0111
Passive multifunction box (not included in Cond 720 Set)	109 810

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.



Conductivity Portable Conductivity Meters

ProfiLine Cond 197i



The WTW conductivity meter of the ProfiLine Cond 197i series is both hoseproof (IP 66) and submersible (IP 67). Along with an 800 data file datalogger, a real time clock and recorder output, the ProfiLine Cond 197i conforms to all GLP requirements. They are equipped with a carrying/support handle and carrying strap as standard.

New: Powerful NiMH rechargeable batteries.

The Cond 197i is suitable for depth measurements down to 330 ft (100 m) in combination with the TA 197 LF depth armature.

IP 66
IP 67



cETLus

3 Year
Warranty

Conductivity depth armature TA 197 LF with built-in temperature sensor, up to 100 m cable with waterproof plug (IP 67), pressure-resistant steel armoring (material VA 1.4571) with screw-off protective hood, pressure-resistant to max. 10 bar, fits into small boreholes (2" dia.).



Highly precise, indestructible, waterproof

Large, silicone keys for field use

Large, easy to read display

Measurement down to depths of 330 ft (100 m)

Technical Data

Model	ProfiLine Cond 197i	
Range/Resolution	LF:	0.0 µS/cm ... 500 mS/cm in 5 measuring ranges or AutoRange, 0.00 ... 19.99 µS/cm for K=0.1 cm ⁻¹ , 0.000 ... 1.999 µS/cm for K=0.01 cm ⁻¹
	Temp.:	23 ... 221 °F (-5.0 °C ... +105.0 °C)
	Salinity:	0.0 ... 70.0
	TDS:	0 ... 1999 mg/l
Accuracy (±1 digit)	LF:	±0.5 % of value
	Temp.:	±0.1 K
Reference temp.	68 ... 77 °F (20 °C or 25 °C), selectable	
Call constant	Calibratable 0.450...0.500 and 0.800...1.200 cm ⁻¹ , fixed: 0.01 cm ⁻¹ freely adjustable 0.25 ... 2.5 cm ⁻¹ and 0.09 ... 0.11 cm ⁻¹	
Temperature comp.	Automatic, can be switched off	
Temperature coefficient	<ul style="list-style-type: none"> • Non-linear function for natural waters to EN 27 888 and ultrapure water function • Linear compensation from 0.01 ... 2.99%/K • No compensation 	

Ordering Information

Portable Conductivity Meter – with wide-range power supply 100-240 VAC (50/60 Hz) included		Order No.
ProfiLine Cond 197i	Robust, waterproof, submersible conductivity meter	3C30-010
Depth armatures down to 330 ft (100 m) see brochure "Product Details"		



Conductivity Handheld Conductivity Meters

Cond 315i, Cond 330i, Cond 340i



Cond 340i in field armoring FM 325 (optional)

Cond 330i in protective armoring SM 325 (optional)

Conductivity Meters

- Robust, shockproof, water-proof
- Up to 1500 hours of continuous operation
- 4-electrode principle allows a wide range of applications



Complete as SET

Useable anywhere, WTW handheld instruments are robust, easy to operate and provide assured accurate readings. From the monitoring of on-line process systems to field studies, WTW's handheld Conductivity meters meet all your measurement needs.



For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.

Parameter

pH

ORP

ISE

Oxygen (D.O.)

Conductivity

Multi-parameter

BOD/Respiration

Photometers

Turbidity

Colony Counter

Software/Printers



Conductivity Handheld Conductivity Meters

WTW handheld conductivity meters are available in three versions:



Cond 315i

Cond 315i

The simple conductivity meter for meeting higher demands. With only 5 keys for error-free measurement with parallel temperature measurement, automatic temperature compensation, 4-electrode measuring technology.

Cond 330i

Robust and waterproof handheld conductivity meter for battery operation; with parallel temperature display, integrated datalogger, GLP supporting functions, automatic and manual temperature compensation with linear temperature function and non-linear function for ultrapure water and natural waters according to EN 27 888. The temperature compensation can be switched off; either 68 °F or 77 °F (20 °C or 25 °C) can be selected as the reference temperature.

Cond 340i

As Cond 330i, but with additional analog and digital RS 232 output and an optional power supply is available.

Technical Data

Model	Cond 315i	Cond 330i and Cond 340i
Range/ Resolution	LF: 0.0 µS/cm ... 500 mS/cm in 5 measuring ranges or AutoRange 0.00 µS/cm ... 19.99 µS/cm (for K=0.1 cm ⁻¹) Temp.: 23 ... 221 °F (-5.0 °C ... +105.0 °C) Salinity: 0.0 ... 70.0 TDS: - Resistivity: 0.000 ... 1999 MΩcm	0.000 µS/cm ... 1.999 µS/cm (for K=0.01 cm ⁻¹) 0 ... 1999 mg/l
Accuracy (± 1 digit)	LF: ±0.5 % of value Temp.: ±0.1 K	
Reference temp. T _{ref}	68 ... 77 °F (20 °C or 25 °C), selectable	
Cell constant	Fixed 0.475 cm ⁻¹ , 0.1 cm ⁻¹	0.475 cm ⁻¹ , 0.1 cm ⁻¹ and 0.01 cm ⁻¹ fixed and 0.25 ... 2.5 cm ⁻¹ and 0.09 ... 0.11 cm ⁻¹ adjustable
Temperature comp.	Automatic, can be switched off	
Temperature coefficient	<ul style="list-style-type: none"> • Non-linear function for natural waters (nLF) to EN 27 888 and ultrapure water function - • Linear compensation from 0.01 ... 2.99 %/K • No compensation 	

Ordering Information

Handheld Conductivity Meter SETs	Order No.
Cond 315i Robust and waterproof battery-operated handheld conductivity meter, including TetraCon® 325, professional case and accessories	2C10-0011
Cond 330i Robust and waterproof battery-operated handheld conductivity meter with datalogger, including TetraCon® 325, professional case and accessories	2C20-0011
Cond 340i Robust and waterproof battery-operated handheld conductivity meter with datalogger and serial interface, including TetraCon® 325, professional case and accessories	2C30-0011
Universal wide-range power supply 100 V - 240 V, 50-60 Hz; for 340i series	902 867
Other measuring cells in SET see brochure "Product Details"	



Conductivity Meters

VARIO C_{ond}

NEW

Simple measurement at a your fingertips – now also available for conductivity measurement!

- Touch screen
- Large operating range
- Plug-in cells – no cables

IP 65 CE 3 Years Warranty



VARIO C_{ond}

VARIO C_{ond} has a lot to offer – at an excellent price. This economical designed meter is ideal for use in process control monitoring or anywhere a small, accurate meter is needed. The VARIO is small, light, handy, waterproof and has a robust firm-grip rubber armoring.

Technical Data

Model	VARIO C _{ond}	
Range and Resolution	[µS/cm]	0.00 ... 19.99 (when using module LR01 V) 0.0 ... 199.9 0 ... 1999
	[mS/cm]	0.00 ... 19.99 0.0 ... 199.9
Resistivity [k*cm]		0.000 ... 1.999 0.00 ... 19.99 0.0 ... 199.9 0 ... 1999
Resistivity [M*cm]		0.00 ... 19.99 0.0 ... 199.9 0 ... 1999
	SAL	0.0 ... 70.0 according IOT
	TDS [mg/l]	0 ... 1999
	T [°F/°C]	41...221/5.0...+105.0

Miniature precision

The globally renowned measurement cell TetraCon® 325 was specially shortened and modified for the VARIO C_{ond}. With extra ultrapure water cell and flow vessel the Vario C_{ond} is uniquely suited for ultrapure water analysis.

Increased precision through the omission of cable connectors – the VARIO C_{ond} is an appropriate solution for servicing and maintaining water treatment equipment. No matter whether using it for pure water measurement in semi-conductor industry or in cell culture laboratories, the pure water conductivity cell with flow-through vessel always allows a rapid and easy control measurement.

Long-lasting power.

VARIO C_{ond} offers up to 500 hours of continuous operation – with just one standard battery. The low power technology shuts down the device after 10 minutes in standby. Changing the battery is easy and quick.



Conductivity Meters

Parameter

pH

ORP

ISE

Oxygen (D.O.)

Conductivity

Multi-parameter

BOD/Respiration

Photometers

Turbidity

Colony Counter

Software/Printers

Ordering Information

VARIO C _{ond}		Order No.
VARIO C _{ond} Set	VARIO C _{ond} in the mobile case set, incl. 4-electrode cell and KCl solution 0.01 mol/l	2X00-001A
VARIO C _{ond} Set	VARIO C _{ond} in the mobile case set, incl. ultra pure water cell and flow-through vessel	2X00-001B
Other accessories see brochure "Product Details"		



Conductivity Cells

Meters

TetraCon®

For more than 50 years WTW has been one of the leading manufacturers worldwide of precision conductivity meters and cells. The TetraCon® 4-electrode system is the result of our commitment to quality and sets a new standard for professional conductivity measurements. In comparison with conventional measuring cells with 2 electrodes, the TetraCon® conductivity cell offers numerous technical advantages:

- Highest degree of precision and linearity by optimized cell geometry
- Extremely large measuring range with just one cell
- Long-term cell constant stability with high-quality abrasion-resistant graphite electrodes
- With built-in temperature sensor as standard
- Smallest immersion depth possible
- No measuring errors even with very dirty electrode – contact resistance on the electrode surface is automatically compensated
- No measuring errors from cable influences
- No measuring errors from primary or secondary polarization effects
- No measuring errors due to contact with side walls or base of measuring vessels
- Robust, unbreakable epoxy body

Selection Guide

Measuring cell	VARIO C _{ond}	Cond 315i	LF 315	LF 318	LF 320/323/325	LF 330/340A	Cond 330i/340i	inoLab® Cond, pH/Cond, Multi	LF 537	LF 538	LF 539	LF 3000	MultiLab® 540	MultiLab® P4/P5	MultiLine® P4, Multi 340i, Multi 197i	MultiLine® P3 pH/LE, pH/Cond 340i	Multi 350i	LF 197	LF 597	Cond 197i
LTA 1			②	②	②	②	②	②	●	●	②		②			②	②			②
LTA 10									●		●	●	②				②			
LR 01/T									●		●	●		●						
KLE 1/T									●		●	●		●						
KLE 315			●																	
TetraCon® 96									●		●	●		●						
TetraCon® 96-1,5									●		●	●		●						
TetraCon® 325		●		●	●	●	●	●		●	●	●	●	●	●	●	●	●	●	●
TA 197 LF																		●		●
TetraCon® 325/Pt																				
TetraCon® DU/T					⑤	⑤	⑤	⑤	④	⑤	④	④	⑤	④			⑤	⑤	⑤	⑤
TetraCon® DU/TH					⑤	⑤	⑤	⑤	④	⑤	④	④	⑤	④			⑤	⑤	⑤	⑤
LR 325/01		●			●	●	●	●		●			●				●	●	●	●
LR 325/001						●	●	●					●				●	●	●	●
TetraCon® 325/S					●	●	●	●		●			●				●	●	●	●
ConOx																	●			
TetraCon® V	●																			
LR01 V	●																			

Adapter (possible conversion with cell constants) is required:
 ② Adapter cable K/LTA together with temperature sensor TFK 325 or TFK 150
 ④ Connection cable KKDU
 ⑤ Connection cable KKDU 325



Conductivity Cells



	Standard conductivity cell		Special conductivity cell TetraCon® 325/S	Ultrapure water conductivity cell		Trace conductivity cell LR 325/001	Conductivity flow-through cell TetraCon® DU/T
	TetraCon® 325	TetraCon® V		LR 325/01	LR01 V		
Order No.	301 960	301 990	301 602	301 961	301 992	301 962	301 252**
Electrode material	Graphite		Graphite	V4A steel		V4A steel	Graphite
Flow-thru vessel	-		-	-		V4A steel	-
Shaft material	Epoxy		Epoxy	V4A steel		V4A steel	Epoxy
Shaft length	4.72 in (120 mm)		4.72 in (120 mm)	4.72 in (120 mm)		4.72 in (120 mm)	6.10 in (155 mm)
Cell constant	K = 0.475 cm ⁻¹		K = 0.491 cm ⁻¹	K = 0.1 cm ⁻¹		K = 0.01 cm ⁻¹	K = 0.778 cm ⁻¹
Diameter	0.60 in (15,3 mm)		0.60 in (15,3 mm)	0.47 in (12 mm)		0.79 in (20 mm)	-
Cable length	4.9 ft (1.5 m)		4.9 ft (1.5 m)	4.9 ft (1.5 m)		4.9 ft (1.5 m)	3.3 ft (1 m) (only with KKDU 325)
Measuring range	1 µS/cm ... 2 S/cm*		1 µS/cm ... 2 S/cm*	0.001 µS/cm ... 200 µS/cm		0.0001 µS/cm ... 30 µS/cm	1 µS/cm ... 2 S/cm*
Temperature range	32 ... 212 °F (0 ... 100 °C)		32 ... 212 °F (0 ... 100 °C)	32 ... 212 °F (0 ... 100 °C)		32 ... 212 °F (0 ... 100 °C)	32 ... 140 °F (0 ... 60 °C)
Filling volume	-		-	17 ml (without sensor)		ca. 10 ml (without sensor)	7 ml
Min./max. immersion depth	36/120 mm	40 mm	40/120 mm	30/120 mm	40 mm	40/120 mm	-

* Measuring range depends on particular instrument, ** Adapter cable KKDU 325 (order no. 301 963), length 3.3 ft (1 m), is necessary for the connection

For additional special measuring cells or other cable lengths see brochure "Product Details"



Conductivity Meters

USP 28 and accessories

Calibration and testing agents



Kit for measuring conductivity according to USP 28

This kit contains LR 325/01 Ultrapure water cell, D01/T flow-through vessel made of glass (USP-KIT 1) or stainless steel (USP-KIT 2), NIST traceable 5 μ S standard with accuracy $\pm 2\%$ and 6R/SET/LabTesting set

Calibration standard 100 μ S/cm

Shelf life 2 years,
NIST traceable with accuracy $\pm 3\%$

Calibration standard 5 μ S/cm

Shelf life 1 year,
NIST traceable with accuracy $\pm 2\%$



Ordering-Information

Kit for measuring the conductivity according to USP 28		Order no.
USP Kit 1	Kit for measuring conductivity according to USP 28, consisting of LR 325/01 Ultrapure water cell, D01/T Glass flow-through vessel, NIST traceable 5 μ S standard with accuracy $\pm 2\%$ and 6R/SET/LabTesting set	300 569
USP Kit 2	As USP Kit 1, but flow-through vessel made of stainless steel instead of D01/T	300 568
Calibration agents		
KS 100 μ S	Calibration standard 100 μ S/cm, shelf life 2 years, NIST traceable with accuracy $\pm 3\%$ (300 ml)	300 578
KS 5 μ S	Calibration standard 5 μ S/cm, shelf life 1 year, NIST traceable with accuracy $\pm 2\%$ (300 ml)	300 580
EP/SET	Calibration and platinization set (6 x 50 ml bottles calibration and control standard, KCl 0.01 mol/l, 30 ml platinizing solution, 1 calibration vessel), only for platinized cells	300 570
E/SET	Calibration set (6 x 50 ml bottles calibration and control standard, KCl 0.01 mol/l)	300 572

USP 28 and accessories



Conductivity measuring kit according to USP 28, with stainless steel flow-through vessel for pharmaceutical water.

Flow-through vessels



Ordering Information

		Order no.
for LTA 1, LTA, LTA 01 and TFK 530		
D 530	Flow-through vessel of transparent PVC, suitable for conductivity cells and temperature sensors, I.D. 44 mm, V*=97 ml	108 060
for TetraCon® 325		
D 201	Flow-through vessel of transparent PVC, I.D. 18 mm, V*=13 ml	203 730
for TetraCon® 96, LTA 100 and KLE 1		
D 1/T	Flow-through vessel, glass I.D. 24 mm, V*=36 ml	302 730
for LR 01/T and LTA 01		
D 01/T	Flow-through vessel, glass I.D. 18 mm, V*=17 ml	302 750
V*: filling volume without sensor		



Multi-Meters

Multi-Parameter Instruments

pH, Oxygen and Conductivity Measurements

Multi-parameter Instruments

With simultaneous function for pH, ORP, Dissolved Oxygen, Conductivity, Salinity and Temperature, these small and compact meters are complete laboratories themselves. WTW multiparameter instruments solve the problem of multiple electrode interference by utilizing our advanced sensor technology. It is now possible to measure oxygen, conductivity and pH simultaneously in the same sample vessel or field location.

Long battery life or optional line adaptors allows WTW's Multiparameter instruments to be used either in the laboratory or in the field. The datalogging function insures accurate reporting. Our standard automatic documentation storage, user defined calibration procedures and GLP functions provide comprehensive quality assurance.





● recommended by WTW ○ conditionally applicable – not recommended

Application Range	inoLab®					Profi-Line	Handheld Meters			
	pH/Cond 720	Multi 720	pH/Cond 740	Multi 740	pH/ION/Cond 750	Multi 197i	pH/Cond 340i	pH/Oxi 340i	Multi 340i	Multi 350i
Routine measurement	●	●	○	○	○	-	○	○	-	○
Routine measurement with documentation	-	-	●	●	●	●	●	●	●	●
AQA with documentation	-	-	●	●	●	●	●	●	●	●
R&D high precision	-	-	●	●	●	-	-	-	-	●
Control measurements	●	●	●	●	●	●	●	●	●	●
LIMS connection	-	-	●	●	●	●	○	○	○	●
Quality assurance	●	●	●	●	●	●	●	●	●	●
Training	●	●	○	○	○	○	●	●	○	○
Service	-	-	-	-	-	●	●	●	●	●
Laboratory measurements	●	●	●	●	●	●	○	○	○	●
Field measurements	-	-	-	-	-	●	●	●	●	●
Depth measurements	-	-	-	-	-	●	-	-	-	up to 330 ft (100 m)
External control/PC connection/PC control	-/-/-	-/-/-	●/●/●	●/●/●	●/●/●	●/●/-	●/●/-	●/●/-	●/●/-	-/●/-
BOD measurements with self-stirring sensor	-	-	●	●	-	●	-	-	-	-
BOD measurements with assessment program	-	-	-	-	-	-	-	-	-	-
pH/ION function	-	-	●	●	●	●	-	-	-	●
Ion-specific measurement programs	-	-	-	-	●	-	-	-	-	-
Salinity/TDS measurement	●/●	●/●	●/●	●/●	●/●	●/-	●/-	-/-	●/-	●/●
Specific resistance	●	●	●	●	●	-	-	-	-	●
Suitable for USP 28	●	●	●	●	●	-	-	-	-	●
Measurement of ultrapure water	●	●	●	●	●	-	-	-	-	●
Trace conductivity	●	●	●	●	●	-	-	-	-	●

see page 52 52 52 52 54 56 60 60 60 57

Application Range Sensors	TetraCon® 325	LR 325/01	LR 325/001	DurOx®	CellOx® 325	ConOx	MPP	TA 197 pH	TA 197 LF	TA 197 Oxi
USP 28	-	●	●	-	-	-	-	-	-	-
Pharmaceutical water	○	●	●	-	-	-	-	-	-	-
Chemical water	○	-	-	○	○	○	○	-	-	-
Ground water	●	●	-	○	-	●	●	-	-	-
Surface waters	●	-	-	●	-	●	●	-	-	-
Depth measurements (barrages)	up to 66 ft (20 m)	-	-	-	up to 66 ft (20 m)	up to 66 ft (20 m)	up to 330 ft (100 m)	up to 330 ft (100 m)	up to 330 ft (100 m)	up to 330 ft (100 m)
Laboratory measurements	●	●	●	○	●	●	-	-	-	-
BOD measurements	-	-	-	-	●	-	-	-	-	-
Fish farming	-	-	-	●	○	●	●	-	-	-
Swimming pools	●	-	-	●	●	●	○	-	-	-
Foods industry (juices)	●	-	-	-	○	○	-	-	-	-
Pharmaceutics	●	●	○	○	●	●	-	-	-	-
Cosmetics/Detergents	○	-	-	-	-	-	-	-	-	-
Semi-conductor industry	-	●	●	-	○	○	-	-	-	-
Prozesstechnik	○	-	-	○	○	-	-	-	-	-
Biotechnology (non-autoclavable)	-	-	-	●	●	●	-	-	-	-
Wastewater treatment plant: aeration basin	-	-	-	●	○	○	●	-	-	-
Wastewater treatment plant: inflow	●	-	-	●	○	○	●	-	-	-

applicable instruments: all with Cond Meas. 350i inoLab® pH/Cond 7xx inoLab® Multi 7xx 350i all with D.O. Meas. 350i 350i Multi 197i Multi 197i Multi 197i

Parameter
pH
ORP
ISE
Oxygen (D.O.)
Conductivity
Multi-parameter
BOD/Respiration
Photometers
Turbidity
Colony Counter
Software/Printers



Multi-Meters

Laboratory Multi-Parameter Instruments



Compact

Versatile

Easy-to-use

IP 43



GETLUS

3 Year Warranty

With the growing number of measurements required by today's modern laboratory, instruments need to be able to perform multiple functions. The inoLab® multi-parameter instruments provide the most necessary features and functions at an outstanding price.

inoLab® 720

inoLab® pH/Cond 720 and inoLab® Multi 720

The routine instrument **inoLab® pH/Cond 720** is a favorably-priced model for initial processing of pH / conductivity measurements. The **inoLab® Multi 720** allows the measurement of pH and ORP, oxygen, conductivity and salinity as well as the temperature. Their large multifunction display means that they are suitable for standard measurements where GLP-documentation is not required.

additional features:

- Easy-to-clean touch-sensitive keypad
- Simplified operator convenience
- Bench and wall-mounted models
- Mains-independent
- Country-specific units
- Application-oriented displays

inoLab® 740

inoLab® pH/Cond 740 and inoLab® Multi 740

As high-end instruments the **inoLab® pH/Cond 740** and the **inoLab® Multi 740** have a resolution of 0.001 pH units. They also have all the functions necessary for making ion selective measurements. With the inoLab® Multi 740 oxygen is measured as %, mg/l or mbar. The pH/Cond 740 and the inoLab® Multi 740 can be used to measure the salinity and TDS (Total Dissolved Solids) as well as conductivity and specific resistance.

The optional built-in printer as well as the graphic display with operator guidance increase the measuring comfort and allow QA-conform documentation.

additional features:

- 5-point calibration by linear regression
- Freely selectable buffer sets
- Graphic evaluation possible
- Built-in digital plotter
- Connection for bar-code reader or PC keyboard
- User Selectable Languages
- Limit input with acoustic alarm
- Multi-Level GLP Functions (password-protected operator levels)
- Free-of-charge downloads for MultiLab® pilot or terminal

Multifunctional

Precise

Communicative

IP 43



GETLUS

3 Year Warranty





Laboratory Multi-Parameter Instruments

Technical Data inoLab®

	pH/Cond 720, Multi 720	Multi 720	pH/Cond 720, Multi 720
	pH measurement	Oxygen measurement	Conductivity measurement
Range/Resolution	pH: -2.00 ... +16.00 pH mV: -199.9 ... +199.9 mV -1999 ... +1999 mV Temp.: 23 ... 221 °F (-5.0 ... +105.0 °C)	O ₂ concentration: 0.00 ... 19.99 mg/l 0.0 ... 90 mg/l* O ₂ saturation: 0.0 ... 199.9% 0 ... 600%* O ₂ partial press.: 0.0 ... 199.9 mbar 0 ... 1250 mbar Temp.: 32 ... 122 °F (0.0 ... 50.0 °C) <i>*depends on oxygen sensor and medium</i>	0.0 µS/cm ... 500 mS/cm in 5 ranges or AutoRange, in addition at: K = 0.01 cm ⁻¹ 0.000 ... 1.999 µS/cm K = 0.1 cm ⁻¹ 0.00 ... 19.99 µS/cm Temperature: 23 ... 221 °F (-5.0 ... +105.0 °C) Salinity: 0.0 ... 70.0 TDS: 0 ... 1999 mg/l Spec. resistivity: 0.000 ... 1999 MΩcm
Accuracy (±1 digit)	pH: ±0.01 pH mV: ±0.3 mV, ±1 mV	±0.5 % of value	±0.5 % of value
Temperature compensation	Automatic 23 ... 221 °F (-5 ... +105.0 °C) with TFK 325 or built-in temperature sensor manual -4 ... 266 °F (-20 °C ... +130 °C)	Automatic via IMT compensation from 32 ... 104 °F (0 ... 40 °C)	Linear and non-linear function for ultrapure water and natural water to EN 27 888, can be switched off
Calibration	MultiCal® 2-point ConCal® 1-/2-point	With OxiCal®-SL	Calibration with 0.01 mol KCl

	pH/Cond 740, Multi 740	Multi 740	pH/Cond 740, Multi 740
	pH measurement	Oxygen measurement	Conductivity measurement
Range/Resolution	pH: -2.000 ... +20.000 pH -2.00 ... +20.00 pH mV: -999.9 ... +999.9 mV -2000 ... +2000 mV ISE: Concentration: 0.01 ... 2000 mg/l Temp.: 23 ... 221 °F (-5.0 ... +105.0 °C)	O ₂ concentration: 0.00 ... 20.00 mg/l 0.0 ... 90 mg/l* O ₂ saturation: 0.0 ... 200.0% 0 ... 600%* O ₂ partial press.: 0.0 ... 200.0 mbar 0 ... 1250 mbar Temp.: 32 ... 122 °F (0.0 ... 50.0 °C) <i>*depends on oxygen sensor and medium</i>	0.0 µS/cm... 2000 mS/cm in 5 ranges or AutoRange, in addition at: K = 0.01 cm ⁻¹ 0.000 ... 2.000 µS/cm K = 0.1 cm ⁻¹ 0.00 ... 20.00 µS/cm Temperature: 23 ... 221 °F (-5.0 ... +105.0 °C) Salinity: 0.0 ... 70.0 TDS: 0 ... 2000 mg/l Spec. resistivity: 0.000 ... 2000 MΩcm
Accuracy (±1 digit)	pH: ±0.004 pH ±0.01 pH mV: ±0.2 mV, ±1 mV	±0.5 % of value	±0.5 % of value
Temperature compensation	Automatic 23 ... 221 °F (-5 ... +105.0 °C) with TFK 325 or built-in temperature sensor manual -4 ... 266 °F (-20 °C ... +130 °C)	Automatic via IMT compensation from 32 ... 104 °F (0 ... 40 °C)	Linear and non-linear functions for ultrapure water and natural waters to EN 27 888, can be switched off
Calibration	MultiCal® 2-, 3- and 5-point ISECal 2- and 3-point ConCal® 1-/2-point	With OxiCal®-SL	Calibration with 0.01 mol/l KCl

Ordering Information

□ with BNC plug ▲ with DIN plug

inoLab® Multi-Parameter SETs – with wide-range power supply 100-240 VAC (50/60 Hz) included		□ Order No.	▲ Order No.
inoLab® Multi 720	Simple and reliable multi-parameter meter for routine work, including Sentix® 41/42, CellOx® 325, TetraCon® 325 and accessories	1F10-211711	1F10-111211
inoLab® Multi 740	Flexible and powerful – the intelligent pH, oxygen and conductivity measuring station consisting of multifunction box and universal terminal, with PC software/connection cable, Sentix® 81/82, CellOx® 325, TetraCon® 325 and accessories	1F30-211911	1F30-111411
Passive multifunction box	(not included in Multi 720 Set)		109 810
Other multi-parameter SETs see brochure "Product Details"			




Multi-Meters

Laboratory Multi-Parameter Instruments

inoLab® pH/ION/Cond 750

- 2 galvanically isolated pH inputs
- Multiple conductivity functions
- Menu-driven with backlit graphics display

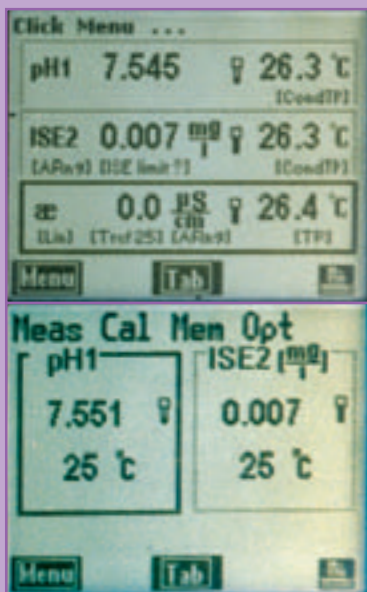

Display with menus

The Ultimate Measurement Technology by WTW:

This ultimate multiparameter meter with its two galvanically separated pH inputs allows independent and simultaneous measurements of pH, ORP or ion concentrations. Calibrations with up to 7 standard solutions can be carried out to achieve precise measurements across the entire characteristic curve of an ISE electrode. The wide range of calibration procedures allows for even the non-linear portion of the calibration curve to be accurately defined. In addition to direct potentiometric methods the following advanced methods are also included:

- Known addition/subtraction
- Addition/subtraction of samples
- Double known addition
- Blank correction
- Known addition with blank correction
- Reference measurement



Laboratory Multi-Parameter Instruments

In addition to pH and Ion capability, and with features such as the ability to measure salinity, TDS, specific conductance, resistance and temperature coefficients, this versatile meter can perform all required conductivity measurements. Temperature coefficients are pre-programmed for NaOH, HCl, KCl, and NaCl. More over that individual (non-) linear coefficients can be fed into the instrument. Additional features include the measurement of sample concentrations by conductivity; two dimensional temperature concentration field to determine accurate reference conductivities. Comprehensive additional functions such as measurement data management, PC-connection via MultiLab® pilot, GLP-conform calibration and measurement data acquisition and a bidirectional RS 232 interface permit seamless integration into modern laboratory routines.



IP 43



cETLus

3 Year Warranty

Technical Data inoLab®

	pH/ION/Cond 750	pH/ION/Cond 750
	pH/ION measurement	Conductivity measurement
Range/Resolution	pH: -2 ... 20.000 pH -2.00 ... 20.00 pH mV: -999.9 ... +999.9 mV -2000 ... +2000 mV Conc.: 0.01 mg/l ... 200 g/l Temp.: 23 ... 221 °F (-5.0 ... +105.0 °C)	0.0 µS/cm ... 2000 mS/cm in 5 ranges in AutoRange mode additional: 0.00 µS/cm ... 20.00 µS/cm (K=0.1 cm ⁻¹) 0.000 µS/cm ... 2.000 µS/cm (K=0.01 cm ⁻¹) Temp.: 23 ... 221 °F (-5.0 °C ... +105.0 °C) Salinity: 0.0 ... 70.0 TDS: 0 ... 2000 mg/l Spec. resistivity: 0.00 ... 2000 MOhm
Accuracy (±1 digit)	pH: ± 0.004 pH, ± 0.01 pH mV: ± 0.2 mV, ± 1 mV	Conductivity: ± 0.5 % of value
Temperature compensation	Automatic 23 ... 221 °F (-5 ... +105.0 °C) Manual -4 ... 266 °F (-20 ... +130 °C) NTC30 KOhm: ±0.1 K Pt 1000: ±0.1 K	23 ... 212 °F (-5.0 ... 100 °C) • Linear and non-linear functions for ultrapure water and natural waters to EN 27 888 • Lineare compensation from 0.01 % ... 3.00 %/K • Compensation can be switched off Special functions: non-linear temperature coefficients, pre-programmed coefficients, concentration determination, non-linear reference conductivity
Reference temp.	—	68 °F/77 °F (20 °C/25 °C) selectable
Cell constants	—	Fixed 0.01 cm ⁻¹ , Freely selectable 0.090 ... 0.110 cm ⁻¹ , 0.250 ... 25.000 cm ⁻¹ Calibrable 0.450 ... 0.500 cm ⁻¹

Ordering Information

□ with BNC plug ▲ with DIN plug

inoLab® Multi-Parameter SETs – with wide-range power supply 100-240 VAC (50/60 Hz) included		□ Order No.	▲ Order No.
inoLab® pH/ION/Cond 750 Flexible and powerful – the intelligent pH, ION and conductivity measuring station consisting of multifunction box, with PC software/connection cable, Sentix® 81/82, TetraCon® 325 and accessories		1K30-211901	1K30-111401
Passive multifunction box (not included in Multi 720 Set)			109 810
Other multi-parameter SETs see brochure "Product Details"			

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.

Parameter

pH

ORP

ISE

Oxygen (D.O.)

Conductivity

Multi-parameter

BOD/Respiration

Photometers

Turbidity

Colony Counter

Software/Printers



Multi-Meters

Portable Multi-Parameter Instruments



ProfiLine Multi 197i

The new WTW ProfiLine Multi 197i is both hose-proof (IP 66) and submersible (IP 67). With its recorder output, real-time clock and 500 data file datalogger standard, these rugged meters conform to all GLP requirements. It allows the simultaneous connection of pH, conductivity and oxygen sensors. The parameter to be measured is set in the display via the "M" function key and can then be measured or stored.

They are equipped with a carrying/support handle and carrying strap as standard.

The Multi 197i has a built-in preamplifier and is therefore suitable in combination with the WTW depth armatures for single-parameter operation at depths down to 330 ft (100 m).

Up to three depth armatures can be simultaneously connected using the adapter ADA/TA 197 pH.

New: Powerful NiMH rechargeable batteries.

- Robust, shockproof
- Waterproof, submersible
- The multi-talent for depth measurements

IP 66
IP 67


CE

cETLUS

3 Year
Warranty

Measurement of Oxygen, pH and Conductivity at Depth Profiles:

Depth armatures with integrated temperature measurement sensors, up to 330 ft (100 m) of cable with a waterproof plug (IP 67), VA 1.4571 steel armoring and screw-on protective hood, pressure resistant to max. 10 bar, suitable for small boreholes (2" diameter).

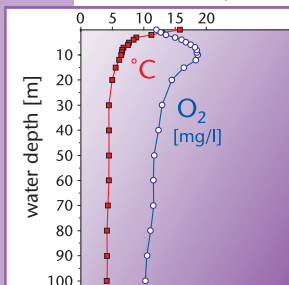

*f.l.t.r. oxygen depth armature TA 197 Oxi and battery-powered stirrer BR 325, pH depth armature TA 197 pH, 4-electrode depth measuring cell TA 197 LF
Please refer to price list for ordering information.*

Technical Data

	pH measurement	Oxygen measurement	Conductivity measurement
Range/ Resolution	pH: -2.00 ... +19.99 mV: -1999 ... +1999	O ₂ concentration 0.00 ... 19.99 mg/l 0.0 ... 90.0 mg/l* O ₂ saturation: 0.00 ... 19.99 % 0.0 ... 600 %* <i>*depends on oxygen sensor and medium</i>	1 μS/cm ... 500 mS/cm in 4 ranges Salinity: 0.0 ... 70.0
Accuracy (±1 digit)	pH: ± 0.01 pH, mV: ± 1 mV	±0.5 % of value	±1 % of value
Temperature compensation	Automatic 23 ... 221 °F (-5 ... +105.0 °C) Manual -4 ... 266 °F (-20 ... +130 °C)	Automatic via IMT Compensation from 32 ... 104 °F (0 ... 40 °C)	Non-linear function for ultrapure and natural waters to EN 27 888
Reference temp.	—	—	68/77 °F (20/25 °C) selectable
Calibration	1-2 point calibration with technical buffers	Automatic calibration	Automatic calibration

Ordering Information

Portable Multi-Parameter Instrument		Order No.
ProfiLine Multi 197i	Robust, waterproof, submersible multi-parameter instrument	3F30-110
Sensors and accessories see brochure "Product Details"		

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.




Multi-Meters

Handheld Multi-Parameter Instruments

Multi 350i



- Multi-functional, High degree of Accuracy
- Flexible
- All parameters simultaneously displayed

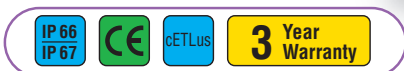

Multi 350i – compact Precision without Compromises

pH, mV, ISE, Oxygen, Conductivity: The new Multi 350i can measure all of these parameters. If desired, pH, oxygen, conductivity and temperature can be measured simultaneously: In the Laboratory using the **combined conductivity/oxygen sensor ConOx**, in the field with the multiparameter probe **MPP 350**. All current WTW pH, combination ISE's, oxygen and conductivity sensors can be additionally connected.

High resolution, high precision, simple, menu-driven operation. Even in poor lighting conditions the backlit graphics display provides clearly readable values. A data-logger, memory for 1,800 data sets and a real-time clock support GLP requirements.

Includes built-in NiMH rechargeable battery for up to 1,000 hours of continuous measurements and appropriate line adaptor.





ConOx

ConOx: Conductivity and Oxygen Measurement with fully automatic Salinity Correction.

A slender and handy sensor for three parameters: conductivity, dissolved oxygen and temperature. The oxygen module can be easily removed and works according to the time-proven galvanic principle. Immediately ready for use, little maintenance required, for all water analysis applications in laboratories and in the field. Including OxiCal® CX calibration vessel.

- Slender
- Handy
- Conductivity, dissolved oxygen and temperature at the same time

The conductivity sensor features the time-proven 4-electrode system. If desired it can also be used without the oxygen module. Precise, easy to clean, capable of accurate measurements even when used in normally electrode fouling samples. Available with cable lengths 4.9 ft, 9.8 ft, 19.7 ft, 32.8 ft, 49.2 ft, 65.6 ft (1.5 m, 3 m, 6 m, 10 m, 15 m, 20 m).

Technical Data

Electrode material	Graphite
Shaft material	Epoxy/POM
Shaft length	145 mm
Cell constant	K=0.475 cm ⁻¹
Diameter	0.60 in. (15.3 mm)
Range	1 µS/cm ... 2 S/cm
Temperature range	32...122 °F(0...50 °C)
Oxygen sensor	Galvanic sensor

Working time: 6 months with 1 electrolyte filling, zero current free



Multi-Parameter Probe MPP 350

An all-new Multi-Parameter Probe as the perfect Probe for use with Multi 350i:

With a diameter of only 1.6 in (41.5 mm) and a length of 11.42 in (29 cm), this probe is usable in a wide range of applications. Wherever simultaneous measurement of pH, dissolved oxygen and conductivity is required, this probe is the ideal supplement to the Multi 350i.

Suitable for measurements in lakes, rivers, seawater and brackish water, in boreholes down to a maximum depth of 330 ft. (100 m), as well in groundwater and spring water. The special pH sensor SensoLyt® MPP-A (please order separately) for measurements in water provides reproducible measurement values, even at low conductivity levels.

For cable lengths greater than 20 ft. (6 m) a battery operated stirrer is included. The conductivity measuring cell with proven 4-pole-measurement technology has a measuring range of 1 µS/cm to 2 S/cm. The MPP 350 is available with 8 different cable lengths up to 330 ft (100 m). Accessories include specialized calibration vessels. Two field sets are available including case, Multi 350i meter, accessories and MPP either with 20 ft. (6 m) or 81 ft. (25 m) cable length.

- pH, conductivity, dissolved oxygen and temperature at the same time
- For all areas of application in surface waters and 2 inch boreholes
- Depth measurement up to 330 ft (100 m)

Techn. Data

Range	pH: 4 ... 12 O ₂ : 0 ... 600% LF: 1 µS/cm ... 2 S/cm Temp.: 32...122 °F (0...50 °C)
Dimensions, Weight	Diameter 41,5 mm, Length approx. 290 to 410 mm (depends on special accessories), approx. 700 g
Materials	POM, Stainless steel 1.4571 (additional weight), PVC (Cable)



Handheld Multi-Parameter Instruments

Technical Data

	pH measurement	Oxygen measurement	Conductivity measurement
Range/Resolution	pH: -2 ... 20.000 -2.00 ... 20.00 mV: -999.9 ... +999.9 -2000 ... +2000 Conc.: 0.01 ... 2000 mg/l Temp.: 23 ... 221 °F (0.0 ... 105.0 °C)	O ₂ conc.: 0.00 ... 20.00 mg/l (19.9 mg/l*) 0.0 ... 90.0 mg/l (90 mg/l*) O ₂ saturation: 0.00 ... 200.0% (200%*) 0.0 ... 600 % O ₂ part. press.: 0.0 ... 200.0 mbar (200 mbar*) 0.0 ... 1250 mbar Temp.: 32 ... 122 °F (0.0 ... 50.0 °C) <i>*if DurOx® oxygen sensor is used</i>	0.0 µS/cm ... 2000 mS/cm in 5 ranges in AutoRange mode additional: 0.00 µS/cm ... 20.00 µS/cm (K=0.1 cm ⁻¹) 0.000 µS/cm ... 2.000 µS/cm (K=0.01 cm ⁻¹) Temp.: 23 ... 221 °F (-5.0 ... +105.0 °C) Salinity: 0.0 ... 70.0 TDS: 0 ... 2000 mg/l Spec. resistivity: 0.00 ... 2000 MOhm
Accuracy (±1 digit)	pH: ± 0.004 pH, ± 0.01 pH mV: ± 0.2 mV, ± 1 mV	O ₂ conc.: ±0.5 % of value O ₂ saturation: ± 0.5 % of value O ₂ part. press.: ±0.5 % of value	Cond: ±0.5 % of value
Temperature compensation	Automatic 23 ... 221 °F (-5 ... +105.0 °C) Manual -4 ... 266 °F (-20 ... +130 °C) NTC 30 kOhm: ± 0.1 K Pt 1000: ± 0.1 K	32 ... 122 °F (0 °C ... 50 °C) (on ambient temperature 41 ... 86 °F / 5 ... 30 °C) <2 % at 32 ... 104 °F (0 ... +40 °C) Temperature: ±0.1 K	23 ... 212 °F (-5.0 ... 100 °C) • Linear and non-linear function for ultrapure and natural waters to EN 27 888 • Linear comp. from 0.01 % ... 3.00 %/K • Compensation can be switched off NTC 30 kOhm: ± 0.1 K Pt 1000: ± 0.1 K
Air pressure compensation	—	Automatic with built-in pressure sensor (500 ... 1100 hPa)	—
Salinity correction	—	Automatic or manual	—
Ref. temperature	—	—	68 °F/77 °F (20 °C/25 °C) selectable
Call constants	—	—	Fixed 0.01 cm ⁻¹ , Freely selectable 0.090 ... 0.110 cm ⁻¹ , 0.250 ... 25.000 cm ⁻¹ Calibrable 0.450 ... 0.500 cm ⁻¹ , 0.800 ... 1.200 cm ⁻¹

Ordering Information

Handheld Multi-Parameter SET		Order No.
Multi 350i SET	Robust and waterproof pocket multi-parameter instrument with data logger and serial interface, including SenTix® 41-3 and ConOx-3, NiMH batteries and battery charger, PC communication package, professional case and accessories	2F40-114B0E
ConOx-3	Combined conductivity-oxygen-probe with 9.8 ft (3 m) cable and accessories	401 010
MPP 350-3	pH/oxygen/conductivity sensor without pH electrode, with 9.8 ft (3 m) cable and accessories	401 100
SensoLyt® MPP-A	armored pH electrode for MPP 350	401 152
SensoLyt® MPP-A Pt	armored ORP electrode for MPP 350	401 152
A 325/S	Stainless steel armoring for ConOx and CelloX®	903 831
SK-325	Protective hood suitable for A 325/S	201 580



- Waterproof
- Robust
- GLP conform

Multi 340i

The rugged, versatile Multimeter

The hose-proof (IP 66) instrument with battery or optional line adaptor also meets the requirements of IP 67 and is optimally suited for use in the field, in laboratories or at production sites.

Simultaneous connection of a pH/ORP electrode and an oxygen sensor or conductivity cell allows up to three parameters (including temperature) to be measured at the same time.

Multi 340i SET 1

- Multi-parameter instrument **Multi 340i**
- Professional case with built-in measuring set-up, two STH 320 stands, two beakers, SM 325 protective armoring and carrying strap with two cases
- Line adaptor power supply, calibration and maintenance supplies, operating instructions

additional features include:

- Handy
- Up to 2500 hours continuous operation
- Easy to use
- Complete in set

IP 66
IP 67



cETLus

3 Year
Warranty

pH/Oxi 340i, pH/Cond 340i

Multi-Parameter Instruments pH/Oxi 340i and pH/Cond 340i

WTW handheld multi-parameter instruments stand for precise multi-parameter measuring technology. The **pH/Oxi 340i** for the determination of pH, oxygen and temperature and the **pH/Cond 340i** for the determination of pH, conductivity and temperature are favorably-priced alternatives to the single parameter instruments for applications which require the measurement of several parameters. The instruments are hose-proof and also meet the requirements of IP 67. They are extremely robust and optimally suited for use in the field, in laboratories or at production sites.

additional features include:

- Handy
- Up to 2500 hours continuous operation
- Easy to use
- Complete in set

pH/Oxi 340i SET

Kit includes:

Professional case with sample beakers included, pH/Oxi 340i, pH electrode and oxygen sensor as chosen, STH 320 stand and calibration and maintenance supplies

pH/Cond 340i SET

Kit includes:

Professional case with sample beakers included, pH/Cond 340i, pH electrode and conductivity cell as chosen, STH 320 stand and calibration and maintenance supplies

IP 66
IP 67



cETLus

3 Year
Warranty





Handheld Multi-Parameter Instruments

Technical Data

	pH/Oxi 340i, pH/Cond 340i, Multi 340i	pH/Oxi 340i, Multi 340i	pH/Cond 340i, Multi 340i
	pH measurement	Oxygen measurement	Conductivity measurement
Range/Resolution	pH: -2.00 ... +19.99 pH mV: -1999 ... +1999	O ₂ concentration: 0.00 ... 19.99 mg/l 0.0 ... 90.0 mg/l* O ₂ saturation: 0.00 ... 19.99 % 0.0 ... 600 %* <i>*depends on oxygen sensor and medium</i>	1 µS/cm ... 500 mS/cm in 4 ranges Salinity: 0.0 ... 70.0
Accuracy (±1 digit)	pH: ± 0.01 pH, mV: ± 1 mV	±0.5 % of value	±1 % of value
Temperature compensation	Automatic 23 ... 221 °F (-5 ... +105.0 °C) Manual -4 ... 266 °F (-20 ... +130 °C)	Automatisch via IMT compensation from 32 ... 104 °F (0 ... 40 °C)	Non-linear function for ultrapure and natural waters to EN 27 888
Ref. temperature	—	—	68/77 °F (20/25 °C) selectable
Calibration	1-2 point calibration with technical buffers	Automatic calibration	Automatic calibration

Protective armoring



for safe on-site use

- ① **SM 325** Shock-absorbing, rubber protective armoring with support handle and sensor cable management
- ② **TG/ML** Sleeve set, suitable for SM 325 protective armoring, consisting of 2 sensor sleeves, holding device and additional carrying strap for field use. Can also be used for storing the sensor
- ③ **FM/ML** Field armoring, specially designed for rough use on-site and in industry, extremely robust and shock-resistant thanks to protective armoring with 2 sensor sleeves, carrying handle and additional carrying strap with holding device, sensor cable management and folding support for laboratory measurements

Ordering Information

Handheld Multi-Parameter Instrument SETs		Order No.
pH/Oxi 340i	Robust and waterproof handheld multi-parameter instrument with datalogger and serial interface for battery operation, including Sentix® 41-3, CellOx® 325-3, professional case and accessories	2D30-101B20
pH/Cond 340i	Robust and waterproof handheld multi-parameter instrument with datalogger and serial interface for battery operation, including Sentix® 41-3, TetraCon® 325-3, professional case and accessories	2E30-101B02
Multi 340i	Robust and waterproof handheld multi-parameter instrument with datalogger and serial interface for battery operation, including Sentix® 41-3, CellOx® 325-3, TetraCon® 325-3, professional case and accessories	2F30-104B22
Universal wide-range power supply 100 V - 240 V, 50 - 60 Hz; for 340i series		902 867
Other electrode/sensors in SET see brochure "Product Details"		

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.



BOD measurements

Respiration

measurement

Biological Oxygen Demand

BOD measurement according to EN 1899-1 and EN 1899-2 and for self-checks

Biological Oxygen Demand (BOD) is an important parameter in water resource management. BOD is a parameter used to measure the quality of water and treatment results in wastewater. In addition, BOD analysis potential is used in the planning and design wastewater treatment facilities.

In routine use BOD determination is used to check the wastewater in the inflow and discharge of wastewater treatment plants. Depending on the measurement site and type of wastewater the BOD value can lie between a few mg/l and several thousand mg/l. Several methods are available for carrying out the measurement.

WTW offers various measuring systems for these methods.

In “**dilution BOD**” the oxygen content of a sample is measured with an oxygen sensor before and after an incubation period of 5 days. The difference between the measurements is the BOD₅ value; this is the **official EPA method**.

In “**BOD self-checks**” with the respirometer the reduction in oxygen causes a definite pressure difference which can be measured by a pressure sensor. This method is very easy to carry out and is a practical method.

Also both methods are very different, the measurements correlate for discharge analysis in municipal wastewater treatment facilities. Both methods requires the samples to be kept at 68 °F (20 °C) for 5 days. WTW offers a wide range of temperature controlled incubators.

Depletion/Respiration

In the course of growing environmental consciousness, microbiological degradability tests have become increasingly important. These can be, for example, soil surveys from waste sites or environmental impact surveys for new chemical substances. The necessary respiration measurements for anaerobic or aerobic degradation can be easily carried out using the OxiTop®-C systems with excellent analysis. WTW offers a wide range of application specific packages with appropriate sample vessels.

User reports and application reports on this topic can be obtained from the WTW homepage (www.wtw.com).

BOD/Depletion/Respiration



Oxi 197i



inoLab® BSB/BOD 740
with StirrOx® G

“Dilution BOD”

according to EN 1899-1/EN 1899-2; official EPA method

- with inoLab® BSB/BOD 740

see p. 64

with easy-to-use analysis program, with PC control.

- with ProfiLine Oxi 197i

see p. 65

Recommended electrode: self-stirring oxygen sensor StirrOx® G



OxiTop® Control



OxiTop® IS 12

“BOD self-check measurement”

worldwide approved method according to the self-check regulations

- OxiTop®

see p. 68

Simple routine measurement, mercury-free pressure measurement

- OxiTop® Control

see p. 69

Routine, standard and special measurement, with automatic sample management

Depletion/Respiration

OxiTop® Control OC 110 for special measurements

see p. 70/74

- Respiration
- Biogas determination
- Soil respiration
- Biodegradability



Biogas
determination



Soil respiration

Accessories/Incubators

Upgrading and general Accessories

see p. 72

Incubators/Thermostat Cabinets

see p. 78



BOD measurements

measurement

"Dilution BOD"

according to DIN EN 1899-1 and DIN EN 1899-2; official EPA method

With inoLab® BSB/BOD 740

flexible and powerful

This laboratory oxygen meter has been specially developed for BOD_n measurements. BOD measurements are determined by regulation EN 1899-1 and is **EPA approved**. You can store up to 7 of your own routines for frequently occurring dilution ratios. A maximum of 30 measuring samples each with 18 dilutions allow the management of up to 540 diluted samples. The inoLab® BSB/BOD 740 can also be used as a conventional high-end oxygen meter (*For technical data such as inoLab® Oxi 740 refer to page 33*). Additional memory and editing options are available when operated using the MultiLab® pilot. The entire measurement and sample management can thus be easily handled via PC.

In combination with StirrOx® G with its automatic start/stop function the inoLab® BSB/BOD 740 is the ideal measuring system for routine oxygen measurement in the BOD₅ determination according to DIN EN 1899-1 and DIN EN 1899-2.


- Simple sample management
- Automatic BOD calculation
- EPA approved

Special features:

- BOD/depletion
- Determination of the biochemical oxygen demand according to DIN EN 1899-1
- Determination of oxygen depletion according to DIN 38 414 P6
- Up to 5 samples for dilution water
- Up to 30 measuring samples
- Up to 18 dilutions per measuring sample
- Up to 7 routines can be stored
- Adjustable incubation time, 5 ... 30 days

Additional features when using the MultiLab® pilot:

- Management of an arbitrary number of samples
- Max. of 32 dilutions per measuring sample
- Max. of 32 dilution waters (blank solutions)
- Adjustable incubation period, 1 to 32 days
- Allocation of names for dilution waters, samples, diluted samples and routines (max. of 255 characters per name)
- Warning indication for BOD values that are too high or too low
- Calculations by mouse click
- Automatic protocols

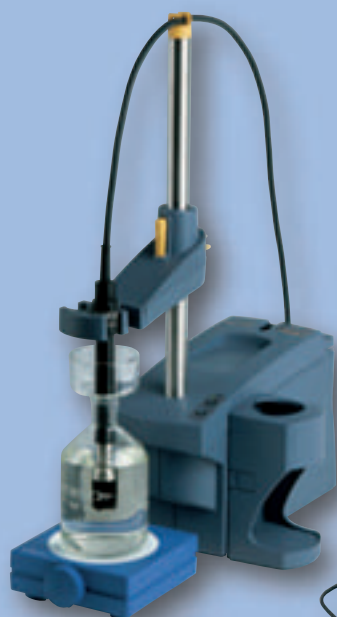

IP 43



GETLUS

3 Year Warranty



“Dilution BOD”

With ProfiLine Oxi 197i

Laboratory oxygen meter ProfiLine Oxi 197i with self-stirring oxygen sensor StirrOx® G.

Technical Data see page 34

- EPA approved method
- Accurate
- Battery and line power operation

Parameter
pH
ORP
ISE
Oxygen (D.O.)
Conductivity
Multi-parameter
BOD/Respiration
Photometers
Turbidity
Colony Counter
Software/Printers



StirrOx® G

Self-stirring oxygen sensor – simultaneous stirring and measurement

- Single-handed operation for series measurements
- Constant flow for high reproducibility
- Immediately ready for measuring – no polarization period required
- Extremely low self-consumption of oxygen – only 0.008 µg h⁻¹ (mg/l)⁻¹
- Zero-current free – no zero point calibration necessary
- With calibration and storage vessel OxiCal®-ST as standard
- Membrane life – Up to 6 months
- Temperature compensation with 2 built-in sensors
- Membrane leakage monitoring – damaged membranes are indicated

Ordering Information

BOD measurement		Order No.
inoLab® BSB/BOD 740P	High-end oxygen/BOD meter including terminal with built-in printer, active multifunction box, with self-stirring oxygen sensor StirrOx® G, wide range power pack and accessories	1H31-0114
ProfiLine Oxi 197i	ProfiLine oxygen meter, extremely robust, hose-water proof (IP 66), RS 232 digital output, operation from mains supply or rechargeable batteries, with wide-range power supply with connection for self-stirring oxygen sensor StirrOx® G and CellOx® 325	3B30-010
StirrOx® G	Self-stirring oxygen sensor for oxygen determination in Karlsruhe bottles, with OxiCal®-ST calibration and storage vessel and accessory case with spare parts and maintenance supplies	201 425



“BOD self-check measurement” / Respiration/ Biogas Determination

With OxiTop® and OxiTop® Control

Mercury-free measurement

Biochemical oxygen demand BOD determination is still one of the most important parameters in water resource management. It can be used to evaluate the impact of biodegradable substances in waters and wastewater. With its OxiTop® systems, WTW offers a unique, modular and mercury-free instrument system. It is not only suitable for BOD determination, but also for measuring biodegradability and depletion.

The advantages of OxiTop® and OxiTop® Control: simple operation, improved controllability and non-toxicity, and measuring ranges of up to 400 000 mg/l BOD (with OxiTop® Control OC 110). As the measured pressure is automatically converted the values can be directly read as mg/l BOD.

- Undiluted samples
- AutoTemp function for delayed start of cold samples
- Non-volatile memory of measured values

Application Range

	OxiTop®	OxiTop® Control OC 100	OxiTop® 110
Application	BOD routine	BOD routine, BOD standard	BOD routine, standard and BOD special, Respiration/Dilution, Soil respiration, Biodegradability, Biogas determination
BOD range	0 – 4,000 mg/l	0 – 4,000 mg/l	0 – 400,000 mg/l
Measured value memory	5 days	0.5 h – 99 days	0.5 h – 99 days
Pressure mode	—	—	Pressure p 500 – 1.350 hPa
Sample volume	Fixed	Fixed	Definable



BOD measurement

OxiTop® Complete Sets for 6 or 12 Measuring Vessels

The complete packages have been assembled so that they contain everything necessary to carry out the measurements. The make up of each package depends on the application and varies by number of vessels, controllers and utensils for sample preparation.

Special stirring platforms were developed in order to maintain a constant temperature and guarantee optimum oxygen distribution in the sample. These stirrer platforms have space for either 6 or 12 standard bottles or 6 large vessels for special applications.

Applicable systems

- **BOD**
OxiTop® IS 6 / IS 12
OxiTop® Control 6/12
- **Soil respiration**
OxiTop® Control B6M / B6
- **OECD / aerobic applications**
OxiTop® Control A6 / A12
OxiTop® Control S6 / S12
- **Biogas determination**
OxiTop® Control AN 6 / AN 12
- **Microbial applications**
OxiTop® Control AN 6 / AN 12
OxiTop® Control A6 / A12

Composition of complete packages



Accessories	OxiTop®		OxiTop® Control			
	IS 6 / IS 12	6 / 12	B6 / B6M / B6M 2.5	A6 / A12	S6 / S12	AN6 / AN12
Vessel with measuring head connection	Amber bottle, 510 ml with rubber sleeve	Amber bottle, 510 ml with rubber sleeve	Duran bottle 500 ml / 1,0 l vessel / 2.5 l vessel; with adapter	1000 ml vessel/ 250 ml vessel with adapter	Amber bottle, 510 ml with rubber sleeve	1000 ml vessel/ 250 ml vessel
Number	6 / 12	6 / 12	6 / 6 / 6	6 / 12	6 / 12	6 / 12
Measuring heads	OxiTop®	OxiTop®-C	OxiTop®-C	OxiTop®-C	OxiTop®-C	OxiTop®-C
Stirrer	IS 6/IS 12	IS 6/IS 12	—	IS 6-Var/IS 12	IS 6/IS 12	IS 6-Var/IS 12
Controller	—	OC 100	OC 110	OC 110	OC 110	OC 110
Software + Cable	—	—	●	●	●	●
CO ₂ absorbent	●	●	●	●	●	●
Nitrification inhibitor	●	●	—	●	●	●
Overflow measuring flask	164 / 432 ml	164 / 432 ml	—	—	—	—
Stirrer bars	6/12	6/12	—	6/12	6/12	6/12
Stirrer bar remover	●	●	—	●	●	●
Blocks of chart paper	●	●	—	—	—	—

see page

68

69

75

76

76

77



BOD self-check measurement

OxiTop® IS 6, IS 12

- High-precision
- 5-day automatic storage of measured values
- Mobile
- Extendable

Technical Data

	OxiTop® measuring head
Measuring principle	Manometric with pressure sensor
Quantity measured	BOD _n
Measuring range	0 ... 40 digits (display units) equals 0 ... 40 / 80 / 200 / 400 / 800 / 2000 / 4000 mg/L BOD
Display accuracy	±1 digit (± 3.55 hPa)
Operating range	500 - 1100 hPa
Memory	For BOD ₅ : 1 per day
Ambient temperature	Storage: -13 ... 149 °F (-25 ... +65 °C) Operations: 41 ... 122 °F (+5 ... +50 °C)
Dimensions	H: 2.7 in (69 mm), Ø 2.8 in (70 mm)

Complete packages for 6 or 12 simultaneous measurements

Measurement using OxiTop® is based on pressure measurement in a closed system: microorganisms in the sample consume the oxygen and form CO₂. This is absorbed by NaOH, creating a vacuum which can be read directly as a measured value in mg/l BOD.

The used sample volume regulates the amount of oxygen available for a complete BOD. Measurement ranges of up to 4,000 mg/l can be measured using different volumes.

The OxiTop® heads (green and yellow for differentiation of inflow/outflow) have an **AutoTemp function**: if the sample temperature is too cold, the start of measurement is automatically delayed (by at least 1 hour) until a constant temperature has been reached.

Apart from the **automatic** storage of 5 measured values (1 value per day), further measured values can be read at all times during or after the period of 5 days, which permits the tracking of check values or measurements over longer periods.



OxiTop® IS 12

Ordering Information

		Order No.
OxiTop® IS 6	Complete package, ready for use, for 6 simultaneous measurements, with IS 6 Inductive Stirring System, for mains operation 230 V / 50/60 Hz and 6 OxiTop® measuring systems, including accessories	208 210
OxiTop® IS 12-6	Complete package, ready for use, for 6 simultaneous measurements (extendable to 12 simultaneous measurements), with IS 12 Inductive Stirring System, for mains operation 230 V / 50/60 Hz and 6 OxiTop® measuring systems, including accessories	208 212
OxiTop® IS 12	Complete package, ready for use, for 12 simultaneous measurements, with IS 12 Inductive Stirring System, for mains operation 230 V / 50/60 Hz and 12 OxiTop® measuring systems, including accessories	208 211

Note: versions for 120 VAC/60 Hz see brochure "Product Details"



BOD measurement

BOD self-check measurement – for a larger number of samples with simple sample management

OxiTop® Control 6, OxiTop® Control 12



OxiTop® Control 12

- Controller-driven
- Simultaneous measurement of up to 100 samples
- Statistical evaluation
- Automatic sample ID
- Worldwide approved method

Complete package for 6 or 12 simultaneous measurements

OxiTop® Control is the logical further development of the successful OxiTop® system using software-controlled functions and infrared interface to communicate with a handy controller, the OC 100. This system enables the simultaneous and grouped start, management, storage and tracking of 100 measuring heads via the controller and tracked on a large display with graphic evaluation. Data can be transferred to the PC for evaluation and documentation via the **AK-540/B** cable (order no. 902 842) and the communication program **Achat OC** (order no. 208 990).

Controller OC 110 in combination with the OxiTop® Control S6 / S12 is ideal for users with other applications apart from BOD (see page 76).

Check sampling progress!

The data can be called up at any time, even during sampling, thus enabling checking of the samples for errors. The display of the progress curve allows immediate detection of irregularities and interferences, such as a BOD value set too high for the volume used or undesired nitrification. Corrections can thus be made at an early stage.





The Controller OC 100 and OC 110



common features

- Simultaneous sample management with option of grouping up to 100 OxiTop®-C measuring heads.
- Data call-up of one parallel sample with statistical evaluation and as individual data.
- Automatic calculation and graphical display of BOD value.
- Data transfer even through glass doors.
- Protocol and documentation of data via Achat OC communication program in combination with a PC
- GLP and AQS with inspection intervals for calibration with the OxiTop® PM calibration tablets
(see page 73: Accessories)

IP54



UL
CUL

1 Year
Warranty

OxiTop® OC 100

Application Range/Technical Data

	OxiTop® Control OC 100	OxiTop® Control OC 110
BOD routine	Individual samples up to 4,000 mg/l	Individual samples up to 4,000 mg/l
BOD standard	Parallel samples with statistical evaluation up to 4,000 mg/l	Parallel samples with statistical evaluation up to 4,000 mg/l
BOD special	—	Freely definable volumes, 0.5 h – 99 days, up to 400,000 mg/l BOD
Soil respiration	—	freely definable volume determination
OECD / Aerobic applications	—	freely definable volume determination
Biogas determination	—	Pressure p 500 - 1350 hPa 10 intermediate values
Data sets per measurement	180 ... 360 (depending on duration)	
Measurement period	0.5 h ... 99 days	
Power supply	3 mignon (AA); alkaline 1.5 V	
Interface	IR (infrared); RS 232 for communication with PC	
Ambient temperature	Storage: -13 °F ... 149 °F (-25 °C ... +65 °C), Operations: 41 °F ... 104 °F (+5 °C ... +40 °C)	
Dimensions	1.7 x 3.9 x 7.9 in (45 x 100 x 200 mm) (H x W x D)	
Weight	Approx. 390 g	



OxiTop®-C Measuring Head



- Instead of the display and keys, the OxiTop®-C measuring head has an infrared interface with which it communicates with Controller OC 100 or OC 110. By “pointing” the controller at an OxiTop®-C measuring head it can be identified and started, data can be called up or deleted and sampling progress can be displayed.
- Each measuring head has its own identification number; this means that manual identification of samples is no longer necessary, even for parallel samples. In addition, statistical evaluations can be easily performed for parallel samples.
- The OxiTop®-C measuring heads have an AutoTemp function for delaying the start of samples which are too cold by up to 4 hours. This mode can be deactivated for BOD standard.
- The measuring heads can store up to 360 data sets. Data are automatically stored in the corresponding interval according to the interval period set (0.5 h to 99 days).
- The built-in pressure sensor can register differences in pressure ranging from 500 to 1,350 hPa.

Technical Data

	OxiTop®-C measuring head
Measuring principle	Manometric with pressure sensor
Measurement of	BOD _n
Pressure range	500 - 1350 hPa
Accuracy	±1 % of value ±1 hPa
Resolution	1 hPa (corresponds to 0.7% of BOD _n measuring range)
Power supply	lithium batteries (280 mAh), 2 x CR2430
Ambient temperature	Storage: -13 ... 149 °F (-25 ... +65 °C) Operation: 41 ... 122 °F (+5 ... +50 °C)
Dimensions	H: 2.8 in (70 mm), Ø 2.8 in (70 mm)

Ordering Information

	Order No.
OxiTop® Control 6 Complete package, ready for use, for 6 simultaneous measurements, with Controller OC 100 and IS 6 Inductive Stirring System, for mains operation 230 V / 50/60 Hz and 6 OxiTop®-C measuring systems, including 6 sample bottles, 6 rubber sleeves, 6 stirrer bars and other accessories	208 201
OxiTop® Control 12 Complete package, ready for use, for 12 simultaneous measurements, with Controller OC 100 and IS 12 Inductive Stirring System, for mains operation 230 V / 50/60 Hz and 12 OxiTop®-C measuring systems, including 12 sample bottles, 12 rubber sleeves, 12 stirrer bars and other accessories	208 204
OxiTop® Control S6/S12 Complete package with Controller OC110 and software	see page 76

Note: versions for 120 VAC/60 Hz see brochure “Product Details”



BOD measurement System Extensions and general Accessories

OxiTop® Measuring Heads and SETs for Retrofitting



In order to meet the growing demand and the extension of potential applications, OxiTop® and OxiTop®-C systems are also available as individual items in different combinations, such as:

- Individual measuring heads OxiTop®/OxiTop®-C
- A set of two OxiTop® heads (yellow and green).
- Upgrade sets for a further 6 positions with 6 heads each and flasks, sleeves and stirring bars, as well as the stirring platform.

Please refer to the brochure "Product Details" for a precise listing of all available components.



Stirrers for BOD measurement

Technical Data

	IS 6	IS 12	IS 6-Var
No. of stirring positions	6	12	6
Stirrer speed	Program-controlled 180 ... 450 min ⁻¹		
Ambient temperature	Storage: -13 °F ... 149 °F (-25 °C ... +65 °C) Operation: 41 °F ... 104 °F (+5 °C ... +40 °C)		
Dimensions (H x W x D in mm)	67 x 265 x 181	67 x 350 x 266	67 x 265 x 181

Stirrers IS 6 and IS 12 have been specially developed for BOD measurement with the OxiTop® system. Software-controlled speed regulation prevents the magnetic stirrer bar from getting caught or wobbling.

The speed is selected so that an optimal gas exchange with the sample takes place. The stirrer is maintenance-free and non-wearing as it contains no moving parts.

The IS 6-Var model has been specially developed for use with large measuring vessels and has space for 6 measuring vessels. Its outer dimensions are identical to those of the IS 12.



Stirrer IS 6 and IS 12

IS 6-Var



BOD measurement

General Accessories

BOD measurement

Testing Aids for the OxiTop® system for Quality Control

Two testing aids are available for monitoring measurement and checking system leakage, which can be called up during a corresponding time interval using the AQA function in the controller.

- **OxiTop® PM**

These calibration tablets simulate a complete BOD and perform quantitative monitoring of measurement (approx. 308 mg/l, batch-dependent) as well as checks for leakage over the entire period.

- **OxiTop® PT**

This testing aid performs a “quick” check for under-pressure and leakage. The OxiTop® contains the pressure table required for the individual place of installation. OxiTop®-C automatically includes these values.

Additional Accessories

- **Storage racks**

For safe storage of OxiTop® measuring systems and OxiTop®-C measuring heads, for 6 measuring heads each.



- **Marking rings**

For identification of BOD bottles for OxiTop® instruments.

- **Overflow measuring flasks in different standard sizes for OxiTop®**

Apart from the supplied 164 ml and 432 ml overflow measuring flasks contained in the standard scope of supply, the following sizes are also available: 22.7 ml, 43.5 ml, 97 ml, 250 ml, 365 ml.

Ordering Information for accessories and spare parts see brochure “Product Details”.



Depletion



Depletion/Respiration with OxiTop® Control OC 110

The investigation and monitoring of biological cleaning treatment processes are becoming more and more important with respect to the environmental requirements such as wastewater treatment, soil remediation and waste treatment.

Biological tests are often in the foreground as well as the usual physical-chemical measuring methods. In order to determine the biodegradability of foodstuffs, pollutants, harmful substances or waste substances and the microbial activity so-called respiration (depletion) measurements are often carried out. In these measurements the respiration of the organisms is determined under defined conditions as the oxygen uptake or release of carbon dioxide.

Measurements are carried out via closed systems using the OxiTop®-C in combination with the OC 110 controller. Depending on the field of use, specially adapted measuring vessels, which are all equipped with the necessary connection thread and some of which are autoclavable, come into operation. For this task there are various complete packages available which come equipped with all required utensils.

For the incubation of larger measuring vessels, WTW offers the TS 1006-i thermostat cabinet as well as the IS 6-VAR stirrer platform, which has been specially designed for vessels with large base diameter.

	Applications and Procedures	Measuring
Soil respiration	Soil analysis/biodegradability of pollutants: laboratory method according to DIN 19 737	Aerobic using CO ₂ absorption, quantitative CO ₂ determination possible
Biodegradability	Determination according to OECD 301 F / DIN EN 29 408 / ISO 9408	Aerobic using CO ₂ absorption
Biogas determination	Determination of anaerobic degradation processes	Anaerobic, determination of CO ₂ + Methane
Microbiology	Growth and stress investigations: determination of the respiration rate	Aerobic, warning pressure possible



Determination of soil respiration

Laboratory method for determining the microbial soil respiration according to DIN 19 737.

OxiTop® Control B6 / BM6

Soil respiration measurements are used for forecasting, surveying and checking remediation work, for biodegradability measurements of substances (pesticides, fungicides, fertilizers, etc.) as well as for carrying out toxicity tests.

This determination is possible with the OxiTop® Control System and special, practically-tested measuring vessels in a very accurate, simple and favorably-priced way.

The expenditure for personnel and apparatus is considerably reduced when compared with conventional methods.

- Simple and precise
- Cost-efficient
- Optimum measuring vessels for subsequent quantitative determination of CO₂



Example of application using MG/... measuring vessels



Example of application using PF/45... sample vessels

Soil respiration measurements can be carried out in 2 different types of vessel.

For actively respiring soils with strong CO₂ development the MG 1.0 measuring vessel is recommended: its large opening (approx. 3.9 in / 100 mm dia.) means that it can be easily used with large-volume CO₂ absorber vessels for afterwards quantitative CO₂ determination.

Ordering Information

Model	Complete soil respiration package	Order No.
OxiTop® Control BM6	Package for soil respiration (aerobic) with 6 MG 1.0 measuring vessels, 1000 ml, with stopper adapters for OxiTop®-C	208 232
OxiTop® Control B6	Package for soil respiration (aerobic) with 6 PF 45/500 sample vessels, 500 ml, Duran and 6 OxiTop® AD/SK adapters, autoclavable	208 230



Determination of biodegradability

Laboratory procedures for determination of biodegradability according to DIN EN 29 408 / ISO 9408 / OECD 301 F

OxiTop® Control A6 / A12 and S6 / S12



The determination of the biodegradability should be checked before “new” chemicals are used for the first time, not only for environmental reasons but also to minimize disposal charges.

The sample and a blank are stirred at a constant temperature for 28 days in closed bottles.

The CO₂ produced is removed from the gas space by means of an absorber so that the resulting negative pressure is a measure of the biodegradability.

The continuous recording of the values in the OxiTop®-C means that the required documentation can be guaranteed in an optimal manner.

The measuring bottles and adapters can be autoclaved at 249.8 °F (121 °C).

Ordering Information

Model	Complete OECD packages	Order No.
OxiTop® Control A6	Package for aerobic applications with 6 x 1000 ml measuring units	208 220
OxiTop® Control A12	Package for aerobic applications with 12 x 250 ml measuring units	208 222
OxiTop® Control S6	Package for aerobic applications with 6 x 510 ml measuring units	208 196
OxiTop® Control S12	Package for aerobic applications with 12 x 510 ml measuring units	208 198
Note: versions for 115 VAC/50/60 Hz see brochure “Product Details”		



Biogas determination

Determination of anaerobic degradation processes: biogas determination

OxiTop® Control AN6 / AN12



Anaerobic degradation processes take place in the absence of oxygen. In order to be able to fill the gas space above the sample with inert gas the measuring bottle has a septum sealed nozzle. When anaerobic degradation has taken place the dissolved CO₂ can be driven off and then removed from the gas space by means of a CO₂ absorber. The resulting pressure difference is proportional to the CO₂ concentration, the remaining overpressure is proportional to the methane concentration.

The degradation process can be comfortably followed in the "pressure" operating mode.

Determination of the Respiration Rate

Microbiological growth and stress investigations: determination of the respiration rate (aerobic/anaerobic measuring operation)



OxiTop® Control AN6 / AN12 and A6 / A12

The use of special measuring bottles with a septum sealed nozzle allows the interference-free addition of substrates and solutions.

Pressure alterations could indicate a reduction in oxygen concentration, for

example, which could make the addition of oxygen or air necessary (possibly other gases as well).

It is possible to set a "warning pressure" or a pressure limit so that the operator can carry out manipulations on the system.

The momentary pressure can be stored so that the manipulation is fully documented. The recording of measured values (max. 10 values) permits long-term measurement.

Ordering Information

Model	Complete packages for microbiology	Order No.
OxiTop® Control AN6	Package for aerobic or anaerobic applications with 6 x 1000 ml measuring units	208 225
OxiTop® Control AN12	Package for aerobic or anaerobic applications with 12 x 250 ml measuring units	208 227
Complete packages for aerobic measurements		
OxiTop® Control A6	Package for aerobic applications with 6 x 1000 ml measuring units	208 220
OxiTop® Control A12	Package for aerobic applications with 12 x 250 ml measuring units	208 222
Note: versions for 115 VAC/50/60 Hz see brochure "Product Details"		



BOD Incubators measurement

OxiTop® Box

- Compact
- Precise
- Uniform temperature distribution



CE 18 Months Warranty

Example of an application:
 OxiTop® Box with OxiTop® Control 12

Thermostat box with forced air circulation for 68 °F (tolerance 67.1 - 68.9 °F / 20 ±0.5 °C)

OxiTop® Box is a benchtop model with hinged clear-view cover which can accommodate a maximum of either 12 OxiTop® simultaneous measurements or 20 Karlsruhe bottles.

The chamber is equipped with a connection for an IS 6 or IS 12 stirrer.

A special compartment is provided for thermostating 6 methylene blue samples.

The box is made of non-corrosive materials and the compressor is CFC-free.

A cross-flow fan ensures uniform temperature distribution.

The box has an automatic defrosting system with condensate evaporation.

Technical Data

Temperature control	68 °F (tolerance 67.1 - 68.9 °F) (20 °C ±0.5 °C)
Ambient temperature	Storage: -13 ... +122 °F (25 °C ... +50 °C) Operation: +50 ... 89.6 °F (+10 °C ... +32 °C)
Power consumption	200 W
Dimensions (H x W x D)	14.76 x 16.73 x 23.62 in (375 x 425 x 600 mm)
Weight	Approx. 30 kg

Ordering Information

BOD thermostat boxes		Order No.
OxiTop® Box	BOD OxiTop® Box, thermostat box with temperature-controlled forced ventilation for 230 V 50 Hz mains supply	208 432
Note: versions for 115 V / 60 Hz see brochure "Product Details"		



Incubators

Thermostat Cabinets



- Versatile
- Powerful
- Inexpensive

To incubate samples at a constant, desired temperature during the reaction period, a thermostat cabinet is necessary. WTW offers thermostat cabinets in various sizes with a variably adjustable temperature range of 50 °F - 104 °F (10 °C - 40 °C) and a power supply of 230 V/50 Hz. Temperature accuracy lies at ±1 °C deviation from the set temperature.

As the samples must be additionally stirred the thermostat cabinets are fitted with internal sockets to provide the stirrers with electricity. 2 – 4 shelves are available, according to the thermostat cabinet size, thus enabling simultaneous thermostating of up to 48 standard BOD samples, or equipping with 4 IS 12 or IS 6-Var stirrer platforms.

The largest model TS 1006-i is especially suited for special applications, as the compartment height between the 4 shelves leaves enough space for 1.5 l vessels or flasks with side nozzles.

The sizes TS 606/2-i aTS 606/4-i are available with transparent insulating glass doors and especially suited for use in combination with the OxiTop® Control system, as data can be called up through the closed glass door. This has the advantage that temperature fluctuations caused by opening the door can be avoided.

Technical Data

	TS 606/2-i	TS 606/3-i	TS 606/4-i	TS 1006-i
Shelves	2	3	4	4 widely spaced
Number of samples	2 x 12 BOD Standard	3 x 12 BOD Standard	4 x 12 BOD Standard	4 x 12 BOD Standard 4 x 6 special vessels
Glass door	Optional	—	Optional	—
Temp. control range	50 °F ... 104 °F (+10 °C ... +40 °C) ±1 K; Adjustment interval: 1 °C			
Ambient temperature	Operation: 50 °F ... 89.6 °F (+10 °C ... +32 °C) (Climate class SN); Storage: -13 °F ... 149 °F (-25 °C ... +65 °C)			
Gross contents	180 l	260 l	360 l	500 l
Dimensions outside: (H x B x D in mm) inside:	850 x 602 x 600 734 x 513 x 433	1215 x 602 x 600 1047 x 513 x 433	1589 x 602 x 600 1418 x 513 x 433	1515 x 755 x 715 1338 x 646 x 516
Weight	37 kg	45 kg	50 kg	72 kg

Ordering Information

BOD thermostat cabinets – only available for 230 V/50 Hz		Order No.
TS 606/2-i	Thermostat cabinet for 2 BOD OxiTop® systems	208 380
TS 606/3-i	Thermostat cabinet for 3 BOD OxiTop® systems	208 382
TS 606/4-i	Thermostat cabinet for 4 BOD OxiTop® systems	208 383
TS 1006-i	Thermostat cabinet for 4 BOD OxiTop® systems	208 385
Other thermostat cabinet see brochure "Product Details"		

Straightforward measuring!

Depending on your application, WTW offers the appropriate photometer and the appropriate test sets. Photometers and tests are perfectly matched to each other. The methods for all test sets are stored in the photometers.

... precision
for all-
purpose
use

Cuvette tests without bar code

Powder tests

Small but nice –
the powder tests
for field use

p. 92



pHotoFlex

p. 86

Cases/sets

The mobile lab
for field use

p. 88



... utmost precision
for lab use



photoLab®
p. 84



Reagents/ accessories

p. 92 / 101

Thermoreactors

The convenient and safe way
to digestions

p. 90



LabStation

The small lab solution:
pHotoFlex plus
LabStation

p. 88





A System of Instrument plus Reagent:

The Photometric Determination



A photometer is used in wastewater, drinking water and environmental analysis to determine the concentration of dissolved substances. For this purpose, specific reagents are added to the dissolved substance to convert it into a measurable color. The color results from the absorption of a certain proportion (wavelengths) of the white light, and hence the measuring is done at the wavelength with the largest absorption. Within a linear range the degree of light absorption can be translated into concentration.

Thus, photometers and test sets make up a system whose components are matched optimally. Modern photometers provide method data for every test. All necessary settings, e.g. wavelength, conversion factors, blank value (= the own color of a solution), etc. are set automatically. When calibrating user defined methods or new reagents, the values for the test are measured, put in and stored as a method.

The characteristics of the test sets are different for each instrument, because the optics are different. Due to the difference in optics, it covers different measurement ranges for the same parameter measured on different photometers.





Portable and Accurate:

The pHotoFlex and photoLab® series

Photometry

In order to choose the appropriate instrument, the following should be considered:

Mobile measuring

With pHotoFlex and pHotoFlex Turb

For fast and accurate measurements in the field consider these features:

- low power consumption
- robust
- mobile
- precise

These requirements are met by a special optical system working with a combination of LED and filters. The robustness of the portable pHotoFlex instruments is due to the low warming and to the longer lifespan of the LEDs utilized. With two cuvette sizes, these photometers allow for all common tests and offer a wide measuring range. The optional LabStation allows the user to conveniently download results collected in the field to a computer back at the lab.

Measuring in laboratory environment

With photoLab® S6/S12 and photoLab® Spektral

Highest demands are considered to be the basis for research, routine measurements and “calculation charges”. Therefore, the instruments have to offer:

- AQS/IOK
- accurate measuring
- wide measuring ranges
- comfort features, e.g. test and cuvette recognition

A complex optical system and a short warm up time guarantee constant measuring conditions. The constant power supply allows the use of bar-codes. The optical system and rectangular cuvettes up to 50 mm allow wide measuring ranges reaching up to trace elements analysis. The largely constant temperature in the lab allows extensive presettings for the methods, thereby providing a higher user comfort.

What are the common features of both series?

- Proven quality, adapted to the respective use
- Highest accuracy corresponding to the used optical system
- A large selection of cuvettes and outstanding instrument features for a simple use of the cuvettes.

Application Photometers

Application range	Portable Photometers		Laboratory Photometers		
	pHotoFlex	pHotoFlex Turb	photoLab® S6	photoLab® S12	photoLab® Spektral
Application areas	Environmental monitoring, water treatment, beverage industry, wine industry, process control, multi-parameter applications for photometry, pH and turbidity.		Routine measurements in wastewater and drinking water, optional field use	Routine measurements in wastewater and drinking water, comprehensive laboratory tests, optional field use	Routine measurements in wastewater and drinking water, professional instrument for comprehensive laboratory tasks in the VIS sector
Wavelengths	6 wavelengths: 436, 517, 557, 594, 610, 690 nm	6 wavelengths: 436, 517, 557, 594, 610, 690, 860 nm	6 wavelengths: 340, 445, 525, 550, 605, 690 nm	12 wavelengths: 340, 410, 445, 500, 525, 550, 565, 605, 620, 665, 690, 820 nm	330 nm – 850 nm stepless
Optical system	LED with filters		Filter/Reference beam	Filter/Reference beam	Zeiss spectrometer module
Special functions	pH measurement optional: LabStation for easy operation and laboratory evaluation with software and barcode reader support	pH measurement, Turbidity (IR)	AQA/IQC	AQA/IQC, Kinetics	AQA/IQC, Kinetics, Absorption spectrum; incl. PC software for easy data transfer
User-defined methods	100		No	50	100
Cuvettes	Round: 16 mm (height: 91 – 104 mm), 28 mm		Round 16 mm	Round and rectangular 10, 20, 50 mm	Round and rectangular 10, 20, 50 mm



photoLab[®] Series

The photoLab[®] Series:

Precision Laboratory Photometers for Every Demand



- AQA/IQC, multistage
- Automatic cuvette identification
- Built-in barcode reader

All the 3 proven models of the photoLab[®] series offer the highest possible convenience plus stringent quality assurance: Open the top cover, insert the cuvette, take the reading!

- Automatic test identification with barcodes
- Automatic cuvette identification
- Automatic self-check
- Quality assurance functions for tests and for the instrument
Password, intervals for instrument check and parameter, test with standards
- Use of rapid tests in reaction cuvettes



Photometry

photoLab® S6

Standard instrument with 6 wavelengths, for all routine determinations in reaction cuvettes (16 mm), especially for wastewater and drinking water.

photoLab® S12

Advanced instrument with 12 wavelengths for test sets in round and in rectangular cuvettes, for wide measuring ranges and low concentrations. In addition, 50 user-defined methods and Kinetics measurements are possible.

photoLab® Spektral

High-grade variable spectral photometer with Zeiss optics for all routine and special tasks in the VIS range: Test sets for round and rectangular cuvettes, Kinetics measurements, recording of absorption and transmission spectrums, as well as 100 user-defined methods with freely selectable wavelength between 330 and 850 nm. Including software Multi/ACHAT II for a comfortable data management and convenient calibration of user-defined methods.

Technical Data

Model	photoLab® S6 and S6-A	photoLab® S12 and S12-A	photoLab® Spektral
Type	Filter Photometer	Filter Photometer	Spectrophotometer with photodiode array technology
Photodiode array for	6 wavelengths	12 wavelengths	–
Wavelengths, nm	340, 445, 525, 550, 605, 690	340, 410, 445, 500, 525, 550, 565, 605, 620, 665, 690, 820	Range 330 ... 850, freely adjustable
User-defined methods	–	50	100
Auto-zero adjustment	Yes	Yes	Yes
AutoSelect-function	Yes	Yes	Yes
Cuvette recognition	Yes	Yes	Yes
Cuvette type	Round	Round, 10 mm, 20 mm and 50 mm	Round, 10 mm, 20 mm and 50 mm
Data storage and time	500 data sets with date and time	1000 data sets with date and time	1000 data sets with date and time
Essential functions	Concentration, absorption and transmission measurement, AQA/IQC, RS 232 interface	Concentration, absorption and transmission measurement, AQA/IQC, Kinetics, RS 232 interface	Concentration, absorption and transmission measurement, AQA/IQC, Kinetics, Absorption spectrum (Abs. +%T), RS 232 interface
Operation with rechargeable batteries (optional)	1 working day, total discharge protection, maintenance charging during mains operation	1 working day, total discharge protection, maintenance charging during mains operation	–
Test marks	CE, UL, CUL	CE, UL, CUL	CE, UL, CUL
Warranty	2 years	2 years	2 years

Ordering Information

Model		Order No.
photoLab® S6	Mains operated version, 230 V European standard plug	250 013
photoLab® S6-A	Version with rechargeable batteries, 230 V European standard plug	250 022
photoLab® S12	Mains operated version, 230 V European standard plug	250 024
photoLab® S12-A	Version with rechargeable batteries, 230 V European standard plug	250 026
photoLab® Spektral	230 V/115 V plug-in transformer with 4 plug adaptors	250 028

Note: versions for other mains supplies/countries on request



pHotoFlex Series

pHotoFlex:

The Portable Photometers

NEW

- Complete Field Case
 - Photometry
 - pH
 - Turbidity
- Undetachable adapter
- Simple push button operation



IP 67



CETLUS

2 Years Warranty



Field case with tray

The new portable photometers offer decisive advantages for field use: They are handy, have a low power consumption and a lot of extra features!

- Smart adapter solution for the use of different cuvettes
- Backlit display with automatic switch-off
- User guidance via display for easy operation without handbook reading
- Large selection of test sets for all requirements
- Method and software update via internet
- Integrated pH measurement with automatic temperature compensation
- Turbidity measurement with infrared light source according to DIN 27027/ISO 7027
- 100 program storage places for user-defined routine measurements
- Battery operation with 4 AA batteries for approx. 3000 measurements
- Optional LabStation with convenient data management for laboratory use
- Optional rechargeable battery set (alternative to LabStation)

The smart adapter solution

An ingenious flap mechanism has been developed to integrate the cuvette adapter undetachably in the instrument. Just slide the magnetic lid up and insert the 28 mm round cuvette. Or, simply flip up the adapter and use a 28 mm round cuvette. The cuvette can have a height of 91 to 104 mm and allows the use of different test sets.



pHotoFlex – portable photometer with pH

The portable photometer pHotoFlex demonstrates its capability with complex tasks in environmental and process monitoring at changing locations.

pHotoFlex features a highly robust optical system, thus being optimally suited for mobile applications under changing conditions. The LEDs plus filters for 6 wavelengths have a remarkably low power consumption and deliver accurate measuring results. The intuitive menu guidance allows a smooth operation even without studying instruction booklets. Dilution functions and timer ease work in special cases.

pH function

The integrated pH function allows measurements of pH 0 ... 16 with automatic buffer recognition (TEC/NIST). Temperature compensation is automatic within the permitted range of 23 ... 212 °F (– 5 ... 100 °C). WTW's MultiCal®-routine allows the automatic calibration with up to 3 calibration points. WTW offers a large selection of pH sensors as an optional accessory: For field use, the maintenance-free SenTix® 41 is recommended, whereas for precision measurements in the laboratory, the SenTix® 81 glass electrode should be used. The electrodes are described in detail in the pH measuring chapter (from p. 19 onward).

NEW

- 100 methods available
- Integrated pH measurement
- Color measurements



with pH sensor SenTix® 41

pHotoFlex Turb – Total Capability



The pHotoFlex Turb is analogous to the pHotoFlex, but has additionally an infrared (IR) light source for nephelometric turbidity measurement (90°), according to the requirements of DIN 27027/ISO 7027. The calibration with the supplied AMCO® standards can be documented and output via RS232 like the measured data.

With the optional LabStation and the LSdata software, all data can be displayed and processed on the computer monitor conveniently and according to GLP (see p. 88).

additionally:

- Turbidity measurement according to DIN 27027/ISO 7027
- 0-1100 NTU/FNU
- Calibration kit (0.02-10-1000 NTU)



pHotoFlex series in a convenient field case!



NEW

- The "in-field laboratory"
- Integrated tray
- Convenient

A small lab for in-field use. The integrated tray with placings for the instrument, cuvettes, measuring beaker and a stand for the pH electrode is particularly practical. Complete sets with:

- pH electrode SenTix® 41 for all pHotoFlex models
- 1 variable pipette with 5 ml volume for all pHotoFlex models
- Calibration standards for pHotoFlex Turb and Turb 430 IR/T
- Many useful accessories: empty cuvettes, buffer solutions with pH 4.01 and 7.00, PC cable AK Labor 540 B, stand for the pH electrode, cleaning tissues, screwdriver for battery change
- Space for other accessories

LabStation plus LSdata –

The smart way of data management!

NEW

The LabStation upgrades the portable pHotoFlex and Turb 430 models to make it a small laboratory solution. With the new software package LSdata the measured data can be processed on a PC conveniently and according to GLP standards. The software is included in the delivery of the LabStation.



- Data export from the instrument to the PC according to GLP and with password protection
- Subsequent processing in Excel format, e.g. for clear documentation of individual sampling points
- Generation, administration and matching between instrument and PC of user-defined methods via dialogue window
- Calculation of calibration curve for user-defined methods

The LabStation also serves as charging station for the rechargeable battery set included in the delivery package. Alternatively a rechargeable battery set for pHotoFlex and Turb 430 models is available separately.



pHotoFlex Turb and Accessories

Accessories

Making your work easier

The rechargeable battery set RB Flex/430

Upgrades pHotoFlex/Turb 430 to rechargeable battery set: The set consists of a rechargeable battery pack with universal power supply. For the connection to a PC via RS 232, the AK540/B cable (Ordering No. 902 842) is required.



LabStation LS Flex/430

The LabStation pHotoFlex LS upgrades the models pHotoFlex and pHotoFlex Turb as well as the turbidity meters Turb 430 IR/T (p. 108) to a small laboratory instrument.

Because of the constant environmental conditions and permanent power supply, the test sets can be run more conveniently via barcode and without new zeroing. Barcodes are offered in the analysis descriptions and on the WTW website (www.WTW.com).

LSdata is an outstanding solution for the management of stored measurement data and for user defined methods too! The delivery includes LSdata, the rechargeable battery set RB Flex/430 and the connection cable AK Labor. The LabStation serves as a charging station for the rechargeable battery set as well.

Technical Data

Model	pHotoFlex	pHotoFlex Turb
Light source	LED	LED
Wavelengths nm	436, 517, 557, 594, 610, 690	436, 517, 557, 594, 610, 690 + 860
User-defined methods	100	100
Timer	3	3
Data storage	1000 data sets	1000 data sets
pH	0-16	0-16
Turbidity	—	0-1100 NTU/FNU
Accuracy	Photometry <2 nm wavelength accuracy, 0.005 abs. reproducibility ±0.01 pH pH Turbidity (NTU/ FNU) —	< 2nm wavelength accuracy, 0.005 abs. reproducibility ±0.01 pH 0.01 NTU/FNU or ±2% of the measured value
Auto-zero adjustment/calibration:	Photometry With start of new method, with LabStation once a day pH / Turbidity 3 point	With start of new method, with LabStation once a day 3 point
Interface	RS 232, USB via adapter (optional)	RS 232, USB via adapter (optional)
Measuring parameters	Photometry, pH	Photometry, pH, Turbidity
Battery	Type AA batteries 4x1.5 V, for approx. 3000 measurements	Type AA batteries 4x1.5 V, for approx. 3000 measurements
Rechargeable battery	Optional: rechargeable battery or LabStation	Optional: rechargeable battery or LabStation
Test marks	cETLus	cETLus
Warranty	2 years	2 years

Ordering Information

	pHotoFlex	Order No.
pHotoFlex	Portable photometer with pH	251 100
pHotoFlex Turb	Portable photometer with pH and turbidity	251 110
pHotoFlex/SET	Portable universal LED filter photometer in a field case with tray to hold instrument and accessories	251 200
pHotoFlex Turb/SET	Portable universal LED filter photometer with integrated turbidity measurement and pH functions in a field case with tray to hold instrument, calibration standard kit and accessories	251 210
FC pHotoFlex/Turb 430	Field case with tray to hold instrument, for all pHotoFlex and Turb 430 models	251 304
LS Flex/430	LabStation for all pHotoFlex and Turb 430 models with LSdata software, rechargeable battery and universal mains adapter	251 301
RB Flex/430	Rechargeable battery for all pHotoFlex models and Turb 430 IR/T with universal plug	251 300

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.



Photometers

Thermoreactors



CR 2200



CR 3200



CR 4200



Thermoreactors for COD and all other thermal Digestion Processes

Thermoreactors are required for the determination of COD, total nitrogen or total phosphorus. They ensure a complete digestion of the sample, as they maintain the necessary high reaction temperature throughout the whole of the defined period.

In each of the thermoreactors from WTW the most important temperatures and digestion times are stored in 7, easily selectable digestion programs. In addition to these 7 fixed standard programs, CR 3200 and CR 4200 thermoreactors allow you to store 8 of your own user-defined programs. Suitable for 16 mm cuvettes.

- Programs for routine tests

NEW Rapid digestion for COD

- Quality assurance with separate sensor

New programs for COD

For the COD digestion, the user can now select among 3 programs: 289.4 °F (148°C), 302 °F (150°C, according to US EPA) for 120 minutes and, at the request of many customers, a **rapid digestion** for 20 minutes at 298.4 °F (148 °C). This timespan has proven sufficient for many purposes in practice.

Safety Precautions

All WTW thermoreactors optimize the heat transmission between the heating block and cuvettes as well as their superior safety. Apart from the built-in safety hood, which prevents chemicals from being splashed about should a cuvette break and the contact protection for the heating block surface, all reactors have timer functions. All reactors display when the reaction temperature is reached.



Thermoreactors

The right Instrument for the right Test!

CR 2200

is ideal for anyone who needs to perform routine water analysis tests with small sample amounts, as 7 programs are available for digestion of 12 sample cuvettes at 212, 248 and 298.4 °F (100, 120, 148 and 150 °C).

CR 3200

In addition, you can program the CR 3200 to carry out 8 of your individual digestions at freely selectable temperatures up to 338 °F (170 °C).

CR 4200

is the right choice for anyone who needs to perform multiple tests simultaneously, such as COD (298.4 °F/148 °C) and total-N (248 °F/120 °C), as the two thermoblocks for 12 cuvettes each can be controlled separately. It also has memory for 8 of your own user-defined programs with free temperature selection up to 338 °F (170 °C).

Quality Assurance:

Quality assurance is constantly increasing in importance, even in the operational analysis sector. The CR 3200 and CR 4200 thermoreactors are both equipped with the external temperature sensor TFK CR (Order No. 250 100) as a testing aid. This temperature sensor can be plugged into the interface in place of a cuvette and the set and actual temperatures can be outputted either to a printer or a PC. This means that the function can not only be monitored, but also documented.

Application Areas and Technical Data

	CR 2200	CR 3200	CR 4200
Application Areas	Routine measurements, wastewater	Routine and specialized tasks in wastewater and in laboratories	Routine and specialized tasks in wastewater and in laboratories
Number of samples, max.:	1 x12	2 x 12 same program	2 x 12, different programs
7 pre-stored programs:	212 °F (100 °C) 60 min, 248 °F (120 °C) with 30 min, 60 min, 120 min, 298.4 °F (148 °C) 120 min, 20 min 302 °F (150 °C) 120 min	212 °F (100 °C) 60 min, 248 °F (120 °C) with 30 min, 60 min, 120 min, 298.4 °F (148 °C) 120 min, 20 min 302 °F (150 °C) 120 min	212 °F (100 °C) 60 min, 248 °F (120 °C) with 30 min, 60 min, 120 min, 298.4 °F (148 °C) 120 min, 20 min 302 °F (150 °C) 120 min
Own programs	–	8 freely selectable 77-338 °F (25-170 °C)	8 freely selectable 77-338 °F (25-170 °C)
Control accuracy	±1 °C ±1 digit		
Safety class	I to DIN VDE 0700 part 1/11.90		
Instrument safety	EN 61010, UL 3101, CAN/CSA C22.2-1010; EN 61010-2-010, IEC-CAN/CSA C22.2-1010.2.010		
Dimensions	W: 10.08 in (256 mm); H: 7.28 in (185 mm), open: 11.42 in (290 mm); D: 12.4 in (315 mm)		

Ordering Information

Model		Order No.
CR 2200	Reactor (230 VAC with Europlug) for COD and other thermal digestions. For up to 12 reaction cuvettes. (Regional power supply available on demand)	1P21-1
CR 3200	Reactor (230 VAC with Europlug) for COD and other thermal digestions. For up to 2x12 reaction cuvettes. (Regional power supply available on demand)	1P22-1
CR 4200	Reactor (230 VAC with Europlug) for COD and other thermal digestions. For up to 2x12 reaction cuvettes in two separately controllable heating blocks. (Regional power supply available on demand)	1P23-1



Photometers

Reagents from A – Z

Reagents –

The right Test for every Application



- Convenient
- Precise
- Assured quality by AQA/IQC

A wide choice of tests is available for routine analysis in different applications. Depending on the optical system and the wavelength employed, photometer and test set make up a matched system with different specific advantages.

For use with portable photometers, test sets should first of all be straightforward. The low consumption LED optics allow the use of easy-to-use and cost-effective test sets, e.g. powder tests. In the laboratory, instruments with bar code and utmost optical sensitivity suggest the use of high-precision tests with bar code reader, certificate and quality assurance support.

WTW continues to expand our reagent offering. Not only are new tests developed, but the usability of tests with different instruments is continuously expanded. Due to the different photometer optics, one and the same test may yield different measuring ranges in different instruments. LED photometers usually have smaller measuring ranges for the same test.

Taking measurements correctly

Every concentration determination is accurate only within the linear absorption range. At the limits of the measuring range you have to expect deviations within the given tolerance. Therefore it may be well worth to repeat the measurement using a test set with a better suited measuring range.

Test Types Overview

Type	Cuvette test	Reagents test	Powder test
Certificate	With certificate (●) for optimum precision Without certificate (TC) for very good precision	With certificate (■) for optimum precision	Without certificate (TP), precise
Test identification	Bar code and/or method selection	Bar code and/or method selection	Method selection, bar code optional
Advantages:	Reaction cuvette with bar code or method selection, 16 mm: Sample adding, inserting, measuring and reading at minimum work, QA support for assured results	Large measuring range, recording of low concentrations in rectangular cuvettes up to 50 mm, QA support for assured results	Compact, straight forward procedure few paraphernalia
Application area:	Laboratory, infrequent work or very large sample throughput	Laboratory, low concentrations, cost-effective routine work with large sample throughput	Mobile measurements, screening and monitoring tasks



Reagents

Model	Measuring Range (Specification max.)	Cuvette (mm)	ml*	Order No.	No.	CC	SW	photoLab®				pHotoFlex
								S6	S12	Spektral		
Acid Capacity up to pH 4.3												
● 01762/1	0.20-8.00 mmol/l 10 - 400 CaCO ₃	round 0.2, 1.0, 5.0	5	252 059	90	-	-	●	●	●	●	
● 01762/2*	0.20-8.00 mmol/l 10 - 400 CaCO ₃	round 0.2, 1.0, 5.0	16		450	-	-	●	●	●	●	
Ag Silver												
■ 14831	0.25 - 3.00 mg/l Ag (total-Ag: 100 °C or 120 °C, 1 h) Digestion reagents are contained in the test set	10, 20	10	250 448	100	-	-	-	●	●	-	
Al Aluminium												
● 00594*	0.02 - 0.50 mg/l Al	round	6.0 + 0.250		25	-	-	●	●	●	-	
■ 14825	0.020 - 1.20 mg/l Al	10, 20, 50, 28	5	250 425	300	✓	✓	-	●	●	●	
TP AI-1 TP	0.00 - 0.25 mg/l Al	28	20	251 400	100	-	-	-	-	-	●	
Ammonium:												
see NH ₄												
Antimony:												
Please ask for application leaflets												
AOX												
● 00675	0.05-2.50 mg/l AOX	round	100	252 023	25	-	-	●	●	●	-	
Arsenic												
■ 01747	0.001 - 0.100 mg/l As	10, 20, 16, 28	350	252 063	30	-	-	-	●	●	●	
<i>also required: AS absorption tube</i>				252 066	1							
Ascorbic acid:												
Please ask for application leaflets												
Au Gold												
● 14821	0.5 - 12.0 mg/l Au	10, 16	2	250 436	80	✓	✓	-	●	●	●	
B Boron												
■ 14839	0.050 - 0.800 mg/l B	10, 20	5	250 427	60	-	-	-	●	●	-	
● 00826	0.05 - 2.00 mg/l B	round	4	252 041	25	-	✓	-	●	●	-	
Br₂ Bromine												
■ 00605	0.020 - 10 mg/l Br ₂	10, 20, 50	10	252 014	200	-	-	-	●	●	-	
C₂H₅OH Alcohol (to be discontinued in course of 2006)												
● 14965	0.40 - 5.00 g/l Alcohol	round	0.2	252 031	25	-	-	●	●	●	-	
C₆H₅OH Phenol												
■ 00856	0.002 - 0.100 mg/l C ₆ H ₅ OH 0.025 - 5.000 mg/l C ₆ H ₅ OH	20 10, 20, 50	200 10	252 058	50 250	-	✓	-	●	●	-	
● 14551	0.10 - 2.50 mg/l C ₆ H ₅ OH	round	10	250 412	25	-	✓	-	●	●	-	
Ca Calcium												
■ 14815	5 - 160 mg/l Ca	10, 20, 16, 28	0.1	250 428	100	-	✓	-	●	●	●	
● 00858	10 - 250 mg/l Ca	round	1	252 047	25	-	-	●	●	●	-	
Cd Cadmium												
● 14834	0.025 - 1.000 mg/l Cd	round	5	250 314	25	✓	-	●	●	●	●	
■ 01745	0.002- 0.500 mg/l Cd	10, 20, 50	10	252 051	55	-	-	-	●	●	-	
Chromium plating bath:												
See reagent-free tests												
● = Reaction cuvettes tests;			TC* = Cuvette test;			CC = CombiCheck test;			ml* = Sample volume;			
■ = Reagent tests;			TP* = Powder pillows;			SW = Sea water;			* = available Q3/2006			



Photometers

Reagents from A-Z

	Model	Measuring Range (Specification max.)	Cuvette (mm)	ml*	Order No.	No.	CC	SW	S6	S12	Spektral	photoFlex
Cl Chloride												
●	14730	5 - 125 mg/l Cl	round	1	250 353	25	✓	✓	●	●	●	●
■	14897	2.5 - 250 mg/l Cl	10	1, 5	250 491	100	✓	✓	-	●	●	●
Cl₂ Chlor												
●	00595	0.03 - 6.00 Cl ₂	round	5	250 419	200	-	-	●	●	●	●
●	00597	0.03 - 6.00 Cl ₂	round	5	250 420	200	-	-	●	●	●	●
■	00598/1	0.010 - 6.00 Cl ₂	10, 20, 50	10	252 010	1200	-	-	-	●	●	-
■	00598/2	0.010 - 6.00 Cl ₂	10, 20, 50	10	252 011	200	-	-	-	●	●	-
■	00599	0.010 - 6.00 Cl ₂	10, 20, 50	10	252 012	200	-	-	-	●	●	-
■	00602/1	0.010 - 6.00 Cl ₂	10, 20, 50	10	252 013	200	-	-	-	●	●	-
■	00602/2	0.010 - 6.00 Cl ₂	10, 20, 50	10	252 055	1200	-	-	-	●	●	-
■	14828	replaced by 00598, 00599, 00602										
■	14732	replaced by ClO ₂ 00608 and Ozone 00607										
TP	Cl-1 TP	0 - 2.00 mg/l Cl ₂ , free	round, 28	10	251 401	100	-	-	-	-	-	●
TP	Cl-2 TP	0.00 - 5.00 mg/l Cl ₂ , free	round, 28	25	251 402	100	-	-	-	-	-	●
TP	Cl-3 TP	0.00 - 2.00 mg/l Cl ₂ , total	28	25	251 414	100	-	-	-	-	-	●
ClO₂ Chlorine dioxide												
■	00608	0.020 - 10.00 mg/l ClO ₂	10, 20, 50, 28	10	252 017	150	-	-	-	●	●	●
■	14732	replaced by ClO ₂ 00608 and Ozone 00607										
ClO₂ Chlorine dioxide/Chlorine/Ozone												
■	14732	replaced by ClO ₂ 00608 and Ozone 00607										
CN Cyanid Cyanid (free and easily liberatable cyanide)												
●	14561	0.010 - 0.500 mg/l CN	round	5	250 344	25	-	-	●	●	●	●
■	09701	0.002 - 0.500 mg/l CN	10, 20, 50	5, 10	250 492	100	-	-	-	●	●	-
Copper plating bath:												
See reagent-free test												
Cr Chromate (chromium VI and total chromium)												
●	14552	0.05 - 2.00 mg/l Cr	round	10	250 341	25	-	✓	●	●	●	●
■	14758	0.01 - 3.00 mg/l Cr	10, 20, 50	5	250 433	250	-	✓	-	●	●	-
CrO₃ Chromium plating bath:												
See reagent-free tests												
Cu Copper												
●	14553	0.05 - 8.00 mg/l Cu	round	5	250 408	25	-	✓	●	●	●	●
■	14767	0.02 - 6.00 mg/l Cu	10, 20, 50, 16, 28	10	250 441	250	-	✓	-	●	●	●
TP	Cu-1 TP	0.00 - 5.00 mg/l Cu	28	10	251 403	100	-	-	-	-	-	●
Cu plating bath:												
See reagent-free tests												
Detergents:												
See Surfactants: anionic, cationic, nonionic												
F Fluoride												
●	14557	0.10 - 1.5 mg/l F	round	5	250 365	25	-	✓	-	●	●	●
■	14598	0.10 - 20.0 mg/l F	10	5 resp. 0.5	252 048	100	-	-	-	●	●	-

● = Reaction cuvettes tests;
■ = Reagent tests;

TC* = Cuvette test;
TP* = Powder pillows;

CC = CombiCheck test;
SW = Sea water;

ml* = Sample volume;
* = available Q3/2006



Reagents

	Model	Measuring Range (Specification max.)	Cuvette (mm)	ml*	Order No.	No.	photoLab®					
							CC	SW	S6	S12	Spektral	pPhotoFlex
Fe Iron												
●	14549	0.05 - 4.00 mg/l Fe	round	5	250 349	25	✓	✓	●	●	●	●
●	14896	1.0 - 50.0 mg/l Fe	round	1	250 361	25	-	-	●	●	●	-
■	14761/1	0.005 - 5.00 mg/l Fe	10, 20, 50, 16, 28	5	250 435	1000	✓	✓	-	●	●	●
■	14761/2	0.005 - 5.00 mg/l Fe	10, 20, 50, 16, 28	5	250 439	250	✓	✓	-	●	●	●
■	00796	0.010 - 5.00 mg/l Fe	10, 20, 50	8	252 042	150	✓	✓	-	●	●	-
TP	Fe-1 TP	0.00 - 1.80 mg/l Fe	28	10	251 404	100	-	-	-	-	-	●
TP	Fe-2 TP	0.00 - 3.00 mg/l Fe	28	10	251 405	100	-	-	-	-	-	●
Formaldehyde:												
See HCHO												
Halogens (total):												
See Cl ₂ , Br ₂ , I ₂ , ClO ₂ , O ₃												
Hazen:												
See reagent-free tests: Coloration												
H₂O₂ Hydrogen peroxide												
●	14731	2.0 - 20.0 mg/l H ₂ O ₂	round	10	250 402	25	-	✓	-	●	●	-
■	18789*	0.10 - 6.00 mg/l H ₂ O ₂	10, 20	8.0		approx. 100	-	✓	-	●	●	-
HCHO Formaldehyde												
●	14500	0.10 - 8.00 mg/l HCHO	round	2	250 406	25	-	-	●	●	●	-
■	14678	0.02 - 8.00 mg/l HCHO	10, 20, 50	3	250 331	100	-	-	-	●	●	-
Heavy metals:												
See lead, cadmium, chromium												
I₂ Iodine												
■	00606	0.050 - 10.00 mg/l I ₂	10, 20, 50	10	252 015	200	-	-	-	●	●	-
Iodine number:												
See reagent-free tests: Coloration												
K Potassium												
●	14562	5.0 - 50.0 mg/l K	round	2	250 407	25	-	✓	●	●	●	●
●	00615	30 - 300 mg/l K	round	0.5	252 020	25	-	✓	●	●	●	-
Mg Magnesium												
●	00815	5.0 - 75.0 mg/l Mg	round	1	252 043	25	-	✓	●	●	●	●
Mn Manganese												
■	01739	0.005 - 2.000 mg/l Mn	10, 20, 50	8	252 056	250	-	-	-	●	●	-
■	14770	0.01 - 10.0 mg/l Mn	10, 20, 50	5	250 442	500	✓	✓	-	●	●	●
●	00816	0.10 - 5.00 mg/l Mn	round	7	252 035	25	✓	-	●	●	●	●
TP	Mn-1 TP	0.0 - 20.0 mg/l Mn	round, 28	10	251 406	100	-	-	-	-	-	●
Mo Molybdenum												
●	00860	0.02 - 1.00 mg/l Mo	round	10	252 040	25	-	-	-	●	●	-
TP	Mo-1 TP	0.0 - 35.0 mg/l Mo	round, 28	10	251 407	100	-	-	-	-	-	●
Monochloramine												
■	01632	0.05 - 10.0 mg/l Cl ₂	10, 20, 50		252 057	150	-	-	-	●	●	-

● = Reaction cuvettes tests; TC* = Cuvette test; CC = CombiCheck test; ml* = Sample volume;
 ■ = Reagent tests; TP* = Powder pillows; SW = Sea water; * = available Q3/2006

Parameter
pH
ORP
ISE
Oxygen (D.O.)
Conductivity
Multi-parameter
BOD/Respiration
Photometers
Turbidity
Colony Counter
Software/Printers



Photometers

Reagents from A-Z

Model	Measuring Range (Specification max.)	Cuvette (mm)	ml*	Order No.	No.	photoLab®					
						CC	SW	S6	S12	Spektral	photoFlex
N₂H₄ Hydrazine											
09711	0.005 - 2.00 mg/l N ₂ H ₄	10, 20, 50	5	250 493	100	-	-	-	●	●	-
Na Sodium											
00885	10 - 300 mg/l Na	round	0.5	252 044	25	-	-	●	●	●	●
NH₄ Ammonium											
14739	0.010 - 2.000 mg/l NH ₄ -N 0.01 - 2.60 mg/l NH ₄ ⁺	round	5	250 495	25	✓	-	●	●	●	-
A5/25	0.20 - 8.00 mg/l NH ₄ -N 0.26 - 10.3 mg/l NH ₄ ⁺	round	1	250 323	25	✓	✓	●	●	●	●
14544	0.5 - 16.0 mg/l NH ₄ -N 0.6 - 20.6 mg/l NH ₄	round	0.5	250 329	25	✓	✓	●	●	●	●
14559	4.0 - 80.0 mg/l NH ₄ -N 5.2 - 103.0 mg/l NH ₄	round	0.1	250 424	25	✓	✓	●	●	●	-
14752	0.010 - 3.00 mg/l NH ₄ -N 0.013 - 3.86 mg/l NH ₄	10, 20, 50, 16, 28	5	250 426	500	✓	✓	-	●	●	●
00683	2.0 - 150 mg/l NH ₄ -N 2.6 - 193 mg/l NH ₄	10	0.1, 0.2	252 027	100	✓	✓	-	●	●	-
TP NH ₄ -1 TP	0.00 - 0.50 mg/l NH ₄ -N 0.00 - 0.64 mg/l NH ₄	28	10	251 408	100	-	-	-	-	-	●
TC NH ₄ -2 TC (LR)	0.00 - 2.50 mg/l NH ₄ -N 0.00 - 3.20 mg/l NH ₄	round	2	251 997	50	-	-	-	-	-	●
TC NH ₄ -3 TC (HR)	0 - 50 mg/l NH ₄ -N 0 - 64 mg/l NH ₄	round	0.1	251 998	50	-	-	-	-	-	●
Ni Nickel											
14554	0.10 - 6.00 mg/l Ni	round	5	250 409	25	✓	-	●	●	●	-
14785	0.02 - 5.00 mg/l Ni	10, 20, 50	5	250 443	250	✓	-	-	●	●	-
Nickel plating bath:											
See reagent-free tests											
Nitrogen (total):											
See N _{Total}											
NO₂ Nitrite											
N4/25	0.020 - 0.600 mg/l NO ₂ -N 0.05 - 2.00 mg/l NO ₂	round	4	250 343	25	-	✓	●	●	●	●
00609*	1.0 - 90.0 mg/l NO ₂ -N 3.28 - 295.2 mg/l NO ₂	16	8.0		25	-	✓	●	●	●	-
14776/1	0.005 - 1.000 mg/l NO ₂ -N 0.016 - 3.29 mg/l NO ₂	10, 20, 50	5	250 445	1000	-	✓	-	●	●	●
14776/2	0.005 - 1.000 mg/l NO ₂ -N 0.016 - 3.29 mg/l NO ₂	10, 20, 50	5	250 440	335	-	✓	-	●	●	●
TP NO ₂ -1 TP	0.00 - 0.33 mg/l NO ₂ -N 0.00 - 1.08 mg/l NO ₂	round, 28	10	251 409	100	-	-	-	-	-	●
TC NO ₂ -2 TC	0.03 - 0.60 mg/l NO ₂ -N (LR) 0.10 - 1.97 mg/l NO ₂ (LR)	round, 16	2	251 994	24	-	-	-	-	-	●
	0.30 - 3.00 mg/l NO ₂ -N (HR) 0.99 - 9.85 mg/l NO ₂	round, 16	0.5								

● = Reaction cuvettes tests;
■ = Reagent tests;

TC* = Cuvette test;
TP* = Powder pillows;

CC = CombiCheck test;
SW = Sea water;

ml* = Sample volume;
* = available Q3/2006



Reagents

	Model	Measuring Range (Specification max.)	Cuvette (mm)	ml*	Order No.	No.	photoLab®					pHotoFlex
							CC	SW	S6	S12	Spektral	
NO₃ Nitrat												
●	14556	0.10 - 3.00 mg/l NO ₃ -N 0.4 - 13.3 mg/l NO ₃	round	2	250 411	25	✓	✓	-	●	●	●
●	N1/25	0.5 - 23.0 mg/l NO ₃ -N 2 - 100 mg/l NO ₃	round	0.5	250 342	25	✓	-	●	●	●	-
●	14542	0.5 - 18.0 mg/l NO ₃ -N 2.2 - 79.7 mg/l NO ₃	round	1.5	250 410	25	✓	-	●	●	●	●
●	14764	1.0 - 50.0 mg/l NO ₃ -N 4 - 221 mg/l NO ₃	round	0.5	250 347	25	✓	-	●	●	●	-
●	00614	23 - 225 mg/l NO ₃ -N 102 - 996 mg/l NO ₃	round	0.1	252 019	25	-	-	●	●	●	-
■	14942	0.2 - 17.0 mg/l NO ₃ -N 0.9 - 75.3 mg/l NO ₃	10, 20, 50	1	250 422	50	✓	✓	-	●	●	●
■	14773	0.2 - 20.0 mg/l NO ₃ -N 0.9 - 88.5 mg/l NO ₃	10, 20	1.5, 3	250 444	100	✓	-	-	●	●	-
■	09713	0.1 - 25.0 mg/l NO ₃ -N 0.45 - 110.7 mg/l NO ₃	10, 20, 50	0.5	250 421	90	✓	-	-	●	●	-
TC	NO ₃ -1 TC	0 - 30.0 mg/l NO ₃ -N 0-133 mg/l NO ₃	round, 16	2	251 993	50	-	-	-	-	-	●
N_{Total} Total Nitrogen												
●	14537	0.5 - 15.0 mg/l N _{Total} (120 °C, 1 h)	round	10	250 358	25	✓	-	●	●	●	●
●	14763	10 - 150 mg/l N _{Total} (120 °C, 1 h)	round	1	250 494	25	✓	-	●	●	●	-
●	00613	0.5 - 15.0 mg/l N _{Total} (120 °C, 1 h)	round	10	252 018	25	✓	-	●	●	●	-
TC	N _{tot} 1 TC (LR)	0 - 25.0 mg/l N _{Total} (120°C, 30 min.)	16	2; 2	251 995	50	-	-	-	-	-	●
TC	N _{tot} 2 TC (HR)	5 - 150 mg/l N _{Total} (120°C, 30 min.)	16	0.5; 2	251 996	50	-	-	-	-	-	●
O₂ BOD Biochemical oxygen demand												
●	00687	0.5 - 3000 mg/l BOD	round	-	252 028	50	-	✓	●	●	●	-
<i>also required:</i> BOD nutrient salt 00688 (252 029) BOD SFR Oxygen reaction bottles 14663 (252 032) <i>available:</i> BOD standard (252 030)												
O₂ COD Chemical oxygen demand												
●	14560	4.0 - 40.0 mg/l COD (148 °C, 2 h)	round	3	250 303	25	✓	-	●	●	●	-
●	C1/25	15 - 160 mg/l COD (148 °C, 2 h)	round	2	250 302	25	✓	-	●	●	●	●
●	14895	15 - 300 mg/l COD (148 °C, 2 h)	round	2	250 359	25	✓	-	●	●	●	-
●	14690	50 - 500 mg/l COD (148 °C, 2 h)	round	2	250 304	25	✓	-	●	●	●	-
●	C2/25	25 - 1500 mg/l COD (148 °C, 2 h)	round	2	250 308	25	✓	-	●	●	●	●
●	14691	300 - 3500 mg/l COD (148 °C, 2 h)	round	2	250 351	25	✓	-	●	●	●	-
●	14555	500 -10000 mg/l COD (148 °C, 2 h)	round	1	250 309	25	✓	-	●	●	●	-
TC	COD1 TC (LR)	0 - 150 mg/l COD (148 °C, 2 h)	round	2	251 990	25	-	-	-	-	-	●
TC	COD2 TC (MR)	0 - 1500 mg/l COD (148 °C, 2 h)	round	2	251 991	25	-	-	-	-	-	●
TC	COD3 TC (HR)	0 - 15000 mg/l COD (148 °C, 2 h)	round	0.2	251 992	25	-	-	-	-	-	●
● = Reaction cuvettes test; TC* = Cuvette test; CC = CombiCheck test; ml* = Sample volume; ■ = Reagent tests; TP* = Powder pillows; SW = Sea water; * = available Q3/2006												

Parameter
pH
ORP
ISE
Oxygen (D.O.)
Conductivity
Multi-parameter
BOD/Respiration
Photometers
Turbidity
Colony Counter
Software/Printers



Reagents from A-Z

							photoLab®				
Model	Measuring Range (Specification max.)	Cuvette (mm)	ml*	Order No.	No.	CC	SW	S6	S12	Spektral	photoFlex
O₂ CSB COD Chemical oxygen demand (HG free)											
● 09772	10 - 150 mg/l COD (148 °C, 2h)	round	2	250 301	25	✓	-	●	●	●	●
● 09773	100 - 1500 mg/l COD (148 °C, 2h)	round	2	250 306	25	✓	-	●	●	●	●
O₂ Oxygen											
● 14694	0.5 - 12.0 mg/l O ₂	round		250 403	25	-	-	●	●	●	-
O₃ Ozone											
■ 00607/1	0.010 - 4.00 mg/l O ₃	10, 20, 50, 28	10	252 016	200	-	-	-	●	●	●
■ 00607/2	0.010 - 4.00 mg/l O ₃	10, 20, 50, 28	10	252 054	1200	-	-	-	●	●	●
■ 14732	replaced by ClO ₂ 00608 and Ozone 00607										
Organic Acids (volatile)											
● 01763	50-3000 mg/l	round		252 060	100	-	-	●	●	●	-
Pb Lead											
● 14833	0.10 - 5.00 mg/l Pb	round	5	250 313	25	✓	-	●	●	●	-
■ 09717	0.010 - 5.00 mg/l Pb	10, 50, 16, 28	8	252 034	50	✓	-	-	●	●	●
pH											
● 01744	pH 6.4 - 8.6	round	10	252 050	280	-	✓	●	●	●	-
Phenol:											
See C ₆ H ₅ OH Phenol											
PO₄ Phosphate											
● P4/25	0.05 - 1.50 mg/l PO ₄ -P 0.05 - 1.50 mg/l P _{Total} 0.20 - 4.50 mg/l PO ₄	round	4	250 366	25	✓	✓	●	●	●	●
● 14543	0.05 - 5.00 mg/l PO ₄ -P 0.05 - 5.00 mg/l P _{Total} 0.2 - 15.3 mg/l PO ₄	round	5	250 324	25	✓	✓	●	●	●	●
● P5/25	0.3 - 15.0 mg/l PO ₄ -P 0.3 - 15.0 mg/l P _{Total} 1.0 - 45.0 mg/l PO ₄	round	0.5	250 368	25	✓	✓	●	●	●	●
● 14546	0.5 - 25.0 mg/l PO ₄ -P 1.5 - 76.7 mg/l PO ₄	round	5	250 413	25	✓	✓	●	●	●	●
● 14729	0.5 - 25.0 mg/l PO ₄ -P 0.5 - 25.0 mg/l P _{Total} 1.5 - 76.7 mg/l PO ₄	round	1	250 334	25	✓	✓	●	●	●	●
● 00616	3.0 - 100.0 mg/l PO ₄ -P 10 - 307 mg/l PO ₄	round	0.2	252 021	25	-	✓	●	●	●	●
■ 14848	0.01 - 5.00 mg/l PO ₄ -P 0.03 - 15.3 mg/l PO ₄	10, 20, 50, 16, 28	5	250 446	420	✓	✓	-	●	●	●
■ 14842	0.5 - 30.0 mg/l PO ₄ -P 1.5 - 92.0 mg/l PO ₄	10, 20	5	250 447	400	-	✓	-	●	●	-
■ 00798	1.0 - 100 mg/l PO ₄ -P 3 - 307 mg/l PO ₄	10	8	252 045	100	-	✓	-	●	●	●
TP PO4-1 TP	0.00 - 0.80 mg/l PO ₄ -P 0.00 - 2.45 mg/l PO ₄	round, 28	10	251 410	100	-	-	-	-	-	●
TC PO4-2 TC	0.00 - 1.60 mg/l PO ₄ -P 0.00 - 4.91 mg/l PO ₄	round, 16	5	251 989	50	-	-	-	-	-	●
TC PO4-3 TC	0.00 - 1.10 mg/l PO ₄ -P 0.00 - 1.10 mg/l P _{Total} (digestion) 0.00 - 3.37 mg/l PO ₄	round, 16	5	251 988	50	-	-	-	-	-	●



Reagents

Model	Measuring Range (Specification max.)	Cuvette (mm)	ml*	Order No.	No.	photoLab®					
						CC	SW	S6	S12	Spektral	pHotoFlex
S Sulfide/Hydrogensulfide											
● 14779	0.02 - 1.50 mg/l S	10, 20, 50	5	250 450	220	-	-	-	●	●	-
Si Silicate/Silicic acid											
■ 14794	0.005- 5.00 mg/l Si	10, 20, 50	5	250 438	300	-	✓	-	●	●	●
■ 00857	0.5 - 500 mg/l Si	10	4/0.5	252 046	100	-	-	-	●	●	●
TP Si-1 TP (LR)	0.00 - 1.60 mg/l SiO ₂ 0.00 - 0.75 mg/l Si	round, 28	10	251 411	100	-	-	-	-	-	●
TP Si-2 TP (HR)	0.0 - 100.0 mg/l SiO ₂ 0.0 - 46.7 mg/l Si	round, 28	10	251 412	100	-	-	-	-	-	●
Sn Tin											
● 14622	0.10 - 2.50 mg/l Sn	round	5	250 401	25	-	✓	-	●	●	-
SO₃ Sulfite											
● 14394	1.0 - 20.0 mg/l SO ₃	round	3	250 416	25	-	-	-	●	●	-
■ 01746	1.0-60.0 mg/l SO ₃	10	2	252 053	150	-	-	-	●	●	-
SO₄ Sulfate											
● 14548	5 - 250 mg/l SO ₄	round	5	250 414	25	✓	✓	●	●	●	●
● 00617	50 - 500 mg/l SO ₄	round	2	252 022	25	✓	✓	●	●	●	-
● 14564	100 - 1000 mg/l SO ₄	round	1	250 415	25	✓	✓	●	●	●	-
■ 14791	25 - 300 mg/l SO ₄	10, 20	2.5	250 449	200	✓	-	-	●	●	-
TP SO4-1 TP	0 - 70 mg/l SO ₄	round, 28	10	251 413	100	-	-	-	-	-	●
Surfactants											
a-Ten (anionic) ● 14697	0.05 - 2.00 mg/l a-Ten	round	5	250 333	25	-	-	-	●	●	-
c-Ten (cationic) ● 01764	0.05 - 1.50 mg/l CTAB	round	5	252 062	25	-	-	-	●	●	-
n-Ten (nonionic) ● 01787	0.10 - 7.50 mg/l Triton X-100	round	4	252 061	25	-	-	-	●	●	-
TOC Total organic carbon											
● 14878	5.0 - 80.0 mg/l TOC	round	3	252 036	25	-	-	●	●	●	-
● 14879	50 - 800 mg/l TOC	round	3	252 037	25	-	-	●	●	●	-
<i>also required: TOC Screw Caps (252 038)</i>											
<i>available: TOC Standard Solution 1000 mg/l (250 499)</i>											
Total nitrogen:											
See N _{Total}											
Total phosphorus:											
See PO ₄ Phosphate											
Water hardness, total hardness											
● 00961	0.7 - 30.1 °d 5 - 215 mg/l Ca	round	1	252 039	25	-	-	●	●	●	●
Water hardness, RH Residual hardness											
● 14683	0.075 - 0.750 °d 0.50 - 5.00 mg/l Ca	round	4	250 404	25	-	-	●	●	●	-
Zn Zinc											
● 00861	0.025 - 1.000 mg/l Zn	round	2	252 049	25	-	-	●	●	●	●
● 14566	0.20 - 5.00 mg/l Zn	round	0.5	250 417	25	✓	-	●	●	●	-
■ 14832	0.05 - 2.50 mg/l Zn	10	5	250 451	90	-	-	-	●	●	-
<i>Necessary reagent: 06146 Zinc reagent 6</i>											
				250 452	180						

● = Reaction cuvettes tests; TC* = Cuvette test; CC = CombiCheck test; ml* = Sample volume;
 ■ = Reagent tests; TP* = Powder pillows; SW = Sea water; * = available Q3/2006

Parameter
pH
ORP
ISE
Oxygen (D.O.)
Conductivity
Multi-parameter
BOD/Respiration
Photometers
Turbidity
Colony Counter
Software/Printers



CombiCheck

CombiCheck solutions are ready-to-use multi-parameter standards. Each package contains a standard solution and an addition solution. Both solutions can be used directly **without dilution** for quality assurance.

- The standard solution is used to check the accuracy of the results for the complete system: procedure - analytical method - reagents - photometer.
- The addition solution is used to check sample-dependent influences by measuring the recovery rate and to determine the most suitable sample preparation method.

The maximum number of determinations which can be made with a **CombiCheck** standard solution depends on the test set used. With the addition solution 280 determinations are always possible.

Read also all test kit instructions!



Storage: 35.6 ... 46.4 °F (+2 ... +8 °C)

Model	Order No.
14676 CombiCheck 10	250 482

Parameter	Concentration	Suitable for test set model	Max. no. of determinations
Ammonium	4.00 mg/l NH ₄ -N	A5/25 14558	90
Chloride	25.0 mg/l Cl	14730	90
COD	80 mg/l COD	C1/25 14540	45 30
Nitrate	2.5 mg/l NO ₃ -N	14556 14773	45 60
Phosphate	0.80 mg/l PO ₄ -P	P4/25 14543 14848	22 18 9
Sulfate	100 mg/l SO ₄	14548 14791 00617	18 40 48

Model	Order No.
14675 CombiCheck 20	250 483

Parameter	Concentration	Suitable for test set model	Max. no. of determinations
Ammonium	12.0 mg/l NH ₄ -N	14544	180
Chloride	60 mg/l Cl	14730	90
COD	750 mg/l COD	C2/25 14541	45 30
Nitrate	9.0 mg/l NO ₃ -N	N1/25 14542 14563 14773 14942 09713	180 60 90 60 60 180
Phosphate	8.0 mg/l PO ₄ -P	P5/25 14729	180 90
Sulfate	500 mg/l SO ₄	14564	90

Model	Order No.
14677 CombiCheck 30	250 484

Parameter	Concentration	Suitable for test set model	Max. no. of determinations
Cadmium	0.500 mg/l Cd	14834	19
Copper	2.00 mg/l Cu	14553 14767	19 19
Iron	1.00 mg/l Fe	14549 14761 00796	19 9 12
Manganese	1.00 mg/l Mn	14770 00816	9 13

Model	Order No.
14692 CombiCheck 40	250 485

Parameter	Concentration	Suitable for test set model	Max. no. of determinations
Aluminium	0.75 mg/l Al	14825	19
Nickel	2.00 mg/l Ni	14554 14785	19 19
Lead	2.00 mg/l Pb	14833 09717	19 11
Zinc	2.00 mg/l Zn	14566	190

Model	Order No.
14695 CombiCheck 50	250 486

Parameter	Concentration	Suitable for test set model	Max. no. of determinations
Ammonium	1.00 mg/l NH ₄ -N	14739 14752	19 19
Nitrogen	5.0 mg/l N _{Total}	14537 00613	9 9
COD	20.0 mg/l COD	14560	32

Model	Order No.
14696 CombiCheck 60	250 487

Parameter	Concentration	Suitable for test set model	Max. no. of determinations
COD	250 mg/l COD	14690 14895	48 48
Chloride	125 mg/l Cl	14897	96

Model	Order No.
14689 CombiCheck 70	250 488

Parameter	Concentration	Suitable for test set model	Max. no. of determinations
Ammonium	50.0 mg/l NH ₄ -N	14559 00683	950 480
COD	5000 mg/l COD	14555	95
Nitrogen	50.0 mg/l N _{Total}	14763	95

Model	Order No.
14738 CombiCheck 80	250 489

Parameter	Concentration	Suitable for test set model	Max. no. of determinations
COD	1.500 mg/l COD	14691	48
Nitrate	25.0 mg/l NO ₃ -N	14764	190
Phosphate	15.0 mg/l PO ₄ -P	14729	95



Photometers

Accessories

Accessories Photometry

Standard Solutions

Parameter	Conc. in mg/l	Amount in ml	Model	Order No.
Aluminium	1000	500	19770	250 460
Ammonium	1000	500	19812	250 461
AOX	20	85 (8-16 Checks)	00680	252 026
Lead	1000	500	19776	250 462
Boron	1000	500	19500	250 463
BOD	210	10 bottles for 10 x 1l	00718	252 030
Cadmium	1000	500	19777	250 464
Calcium	1000	500	19778	250 465
Chloride	1000	500	19897	250 466
Chromium	1000	500	19779	250 467
Chromate	1000	500	19780	250 468
COD 160	100	30	KCSB 100	250 356
COD 1500	400	30	KCSB 400	250 357
Iron	1000	500	19781	250 469
Fluoride	1000	500	19814	250 470
Potassium	1000	500	70230	252 471
Silicic acid (Silicon)	1000	500	70236	252 472
Copper	1000	500	19786	250 473
Manganese	1000	500	19789	250 474
Nickel	1000	500	19792	250 475
Nitrate	1000	500	19811	250 476
Nitrite	1000	500	19899	250 477
Phosphate	1000	500	19898	250 478
Silver	1000	500	19797	250 479
Sulfate	1000	500	19813	250 480
TOC	1000	100	09017	250 499
Zinc	1000	500	19806	250 481

Standard solutions which limited stability, to be freshly prepared at regular intervals:

- Free chlorine
- Bound chlorine
- Formaldehyde
- Hydrazine
- Hydrogen peroxide
- Hydrogen sulfide
- Phenol
- Silicon
- Sulfide
- Sulfite
- Anionic surfactants

PhotoCheck

AQA/IQC: Comprehensive testing aid for optics and measurement linearity!

The stable colored solutions are used for checking the filter and the wavelength settings 445 nm/446 nm, 520 nm/525 nm as well as 690 nm. The correctness of the wavelength setting and the linearity of the extinction measurement are checked with 4 solutions per wavelength. The control is fast and comfortable, via a simple, menu-guided function. The traceability of this testing aid to international standards is guaranteed by checking the solutions in a reference photometer monitored with primary standards (NIST standards). These values are documented accordingly.

PipeCheck

Testing aid for the right pipetting volume!

The appropriate test solution is diluted with distilled water using the pipette to be checked and the extinction of the dilute solution is compared with that of a reference solution. Pipettes with a variation in volume of more than $\pm 2.5\%$ must be regarded as being faulty.



Reagent-free Tests

% transmittance

0 - 100 % T, 10, 20, 50 mm cuvette (self-absorption).

FAU turbidity

(EN ISO 7027) Determination of turbidity

Turbidity is caused in liquids by the presence of undissolved substances. For undissolved finely dispersed substances the turbidity can be measured by measuring the reduction in the intensity of a beam of light when passed through the liquid, or by measuring the intensity of the scattered radiation.

A formazin solution, which must be freshly prepared and is not commercially available, is used as a reference solution. According to EN ISO 7027, all instruments may be used which satisfy the following requirements: Incoming radiation at 860 nm. The results are given in FAU units (Formazin Attenuation Units) when the radiation passing through is measured.

Extinction

According to the Lambert-Beer law, the extinction $E = \epsilon(\lambda) \cdot c \cdot d$ is proportional to the concentration of substances contained in the water. The proportionality constant $\epsilon(\lambda)$ depends on the wavelength. These constants, and other data required for the determination of the solids in the water are stored in modern photometers as method data. The basic quantity measured is and remains the extinction.

Coloration

(EN ISO 7887: 1994)

If a layer of several meters of pure water is observed in transmitted light it appears to have a weakly blue coloration. This coloration can alter in the presence of contaminants to form a wide range of colorations. Natural waters are usually colored yellow-brown by contamination with iron or clay particles or humic matter.

(A green coloration can be produced by algae.) The "true" color of water is determined after filtration through a 0.45 μm filter

Normally most yellow-brown colored waters and the outflows of municipal sewage treatment plants can be measured at 436 nm. The outflows of industrial wastewater treatment plants show no sharp and distinctive extinction maxima. For the investigation of such water it is obligatory to measure at 436 nm (mercury line); the two other measuring wavelengths 525 nm and 620 nm can, depending on the filter used, vary slightly from these wavelengths. For discontinuous measurements the standard permits the use of filter photometers with a spectral bandwidth of < 20 nm for measurements at 436 nm, 525 nm and 620 nm. Thus, for example, instruments with 445 nm and 520 nm interference filters with a bandwidth of 10 nm are also suitable. For comparability with the standard, however, a spectrophotometer is required. The results are presented in m^{-1} together with the measuring wavelength, spectral bandwidth, water temperature and pH.

In some publications the result is given in DFZ (translucent coloration number); which is identical with the m^{-1} result.

(DIN ISO 6271: 19988)

Clear liquids, determination of the color number with the platinum-cobalt scale (Hazen color number, APHA color number). Spectrophotometers are mentioned as being suitable for measuring the stock solutions at 430 nm, 455 nm, 480 nm and 510 nm. According to the standard the measurement itself is carried out with a color comparator which allows a visual comparison.

Chrome-plating bath

Reagent-free measurement of the self-coloration of an electroplating bath. 5 ml of the sample are pipetted into a 100 ml volumetric flask, filled up to the mark with distilled water and well mixed. 4 ml of the diluted sample are pipetted into a 100 ml volumetric flask, filled up to the mark with distilled water and well mixed. 5 ml of the 1:500 dilution are placed in a screw-cap glass and 5 ml 40% sulfuric acid are added. The glass is sealed and the contents well mixed. The solution is filled into a rectangular cuvette for the measurement.

Nickel-plating bath

Reagent-free measurement of the self-coloration of an electroplating bath. 5 ml of the sample are pipetted into a round cuvette and 5 ml 40% sulfuric acid are added. The cuvette is sealed and the contents mixed. The solution is filled into a rectangular cuvette for the measurement.

Copper-plating bath

Reagent-free measurement of the self-coloration of an electroplating bath. 25 ml of the sample are pipetted into a 100 ml volumetric flask, filled up to the mark with distilled water and well mixed. 5 ml of the diluted sample are placed in a screw-cap glass and 5 ml 40% sulfuric acid are added. The glass is sealed and the contents well mixed. The solution is filled into a rectangular cuvette for the measurement.



Turbidity

Quality Control using Turbidity Measurements

- AMCO®-standards
- AQA functions
- US EPA

Typical turbidity values for various liquids

Liquid	NTU
Deionized water	0.02
Drinking water	0.02 ... 0.5
Spring water	0.05 ... 10
Wastewater (untreated)	70 ... 2000
White water (paper industry)	60 ... 800

Please note:

As floating and moving particles are measured in turbidity, slight measurement deviations are possible. In order to achieve results which are as representative as possible, attention should be paid to the following:

- samples should be measured immediately, as particle otherwise settle.
- constant lamp operating temperature.
- condensation on samples should be avoided.
- the position of the standards should be marked to exclude the influence of glass inhomogeneities.

Turbidity measurements are of extreme importance in quality monitoring in water, wastewater, beverage production, electroplating and petrochemical applications.

Light passing through liquid which contains undissolved solids, such as algae, mud, microbes and other insoluble particles, is both absorbed and scattered. Turbidity increases with the amount of undissolved solids present in the sample. However, the shape, size and composition of the particles also influence the degree of turbidity. Turbidity has been determined by simply measuring light passing through the sample. Measuring the **scattered light at an angle of 90°** has proved to be a more accurate method particularly at lower measuring ranges. Instruments that use this method are also referred to as **nephelometers**.

Turbidity or nephelometers instruments differ by the light source they utilize. Infrared units (IR-LED) with a wavelength of 860 nm are required for methods: ISO 7027/ DIN EN 27027 (EN ISO 7027). Standard methods specifies the use of units that use white light by a tungsten wide-band lamp for water and wastewater analysis.

Which light source – infrared (IR) or white light (tungsten)?

An infrared light source minimizes or even eliminates the influence of coloration in a solution, because there is practically no absorption at a wavelength of 860 nm. The detection sensitivity for small particles, on the other hand, is somewhat lower at this wavelength because of the generally lower light scattering of small particles.

White light has a higher sensitivity for small particles, however the inherent coloration of the solution has a stronger disturbing effect in this case.

The IR measurement is required by DIN ISO, the white light measuring by US EPA.

Nephelometric or transmittance measuring?





The nephelometric measuring with 90° scattered light is advantageous for the lower turbidity range. In contrast, the transmittance method is better suited for medium to high turbidities, where there are stronger light scattering and shadowing effects among the larger-sized particles. In this case, the diminution of light intensity gives a more accurate result than a 90° scattered light metering. Laboratory instruments for higher turbidity ranges therefore offer significantly more measuring options.



The Right Instrument for the right Use!

4 models to choose from:

2 portable models, each with either IR or tungsten light source, and
2 laboratory meters with IR or Tungsten light source:

Applications		Measuring range	Calibration	Special features
Applications	Light source			
 <p>Turb 355 T/IR</p> <p>Portable use for waste water, surface water and ground water applications</p>	Tungsten lamp/IR LED	0-1100 NTU/FNU	Automatic 1-3 point	Portable field meter
 <p>Turb 430 T/IR</p> <p>Portable use for all water testing applications incl. drinking water, wine industry, process control</p> <p>Laboratory use; optional for all applications up to 1100 NTU/FNU with LabStation</p>	Tungsten lamp/IR LED	0-1100 NTU/FNU	Automatic 3 point	Portable field meter Calibration interval Calibration documentation Storage for measured values Optional: LabStation for -Easy data evaluation with LSdata -Laboratory use -As a charging station
 <p>Turb 550/Turb 550 IR</p> <p>Routine meter for laboratories; drinking water</p>	Tungsten lamp/IR LED	0-1000 NTU/FNU	Automatic 1-3 point	AQA Flow-through measurement (unpressurized)
 <p>Turb 555/Turb 555</p> <p>Routine meter for precision measurements</p>	Tungsten lamp/IR LED	0-10000 NTU/FNU/FAU	Automatic 1-5 point	AQA complete with password protection, ratio method for the reduction of interferences; transmission, flow-through measurement (unpressurized/ up to 4 bar)



Turbidity

Laboratory Turbidity Meters

Turb 550 / Turb 550 IR



- AutoRange
- Automatic 1-3 point calibration
- Flow-through measurement

THE professional turbidity meter – Up to 1,000 NTU

Laboratory turbidity meters for nephelometric measurements with automatic 1-3-point calibration and calibration interval monitoring. Measuring range selection from 0.01 ... 1000 NTU is carried out automatically and for comparative measurements the current and previous values can be shown on the 2-line display.

Standard equipment includes instrument with built-in short operating instructions, 3 empty cuvettes and 3 standards (0.02 – 10.0 – 1000 NTU, AMCO® standards with approval for drinking water as primary standards according to US EPA and according to EN ISO 7027).

An unpressurized flow-through adapter is available for continuous measurements.



Technical Data

	Turb 550	Turb 550 IR	Turb 555	Turb 555 IR
Measuring principles	Nephelometric	Nephelometric	Nephelometric ratio methode transmission	Nephelometric ratio methode transmission
Light source	Tungsten lamp	IR-LED	Tungsten lamp	IR-LED
Measuring range	NTU 0 ... 1000 FNU – EBC – Nephelos – FAU –	0 ... 1000 0 ... 1000	0 ... 10000 – 0 ... 2450 0 ... 67000 –	0 ... 10000 0 ... 10000 0 ... 2450 – 0 ... 10000
Resolution	0.01 NTU from 0.00 ... 9.99 0.1 NTU from 10.0 ... 99.9 1 NTU from 100 ... 1000		0.0001 NTU from 0.0001 ... 9.9999 NTU 0.001 NTU from 10.000 ... 99.999 NTU 0.01 NTU from 100.00 ... 999.99 NTU 0.1 NTU from 1000.0 ... 9999.9 NTU	
Accuracy	±2% of value or ±0.01 NTU		0 ... 1000 NTU: ±2% of value or ±0.01 NTU 1000 ... 4000 NTU: ±5% of value 4000 ... 10000 NTU: ±10% of value	
Reproducibility	±1% of value or ±0.01 NTU			
Calibration	Automatic 1...3 point calibration		Automatic 1...5 point calibration	
Response time	< 3 seconds		< 6 seconds	
Cuvettes	1.1 x 2.76 in (28 x 70 mm) round cuvette, 25 ml sample volume			
AQA functions	Calibration interval monitoring Calibration protocol		Calibration interval monitoring Calibration protocol Password-protected access to calibration and configuration time-controlled data transmission	
Operating temp.	50 ... 104 °F (+10 ... +40 °C)		32 ... 122 °F (0 ... +50 °C)	
Power supply	Plug-in power supply 100 - 240 VAC ±10% / 47 - 63 Hz			



Laboratory Turbidity Meters

Turb 555 / Turb 555 IR



The ADVANCED professional meter – Measuring range up to 10,000 NTU

High-precision laboratory turbidity meter with a wide measuring range of 0.0001 to 10000 NTU (automatic measuring range switching) for all turbidity measuring applications from ultrapure and drinking water measurements, through quality assurance in soft drinks and wastewater treatment.

The measuring system with its 4 detectors allows not only nephelometric (90° scatter) measurements and transmittance measurements, but also ratio measurements in which the influences of stray light and sample color are reduced.

Comprehensive AQA functions such as monitoring the calibration interval or password protection for calibration and setup access fulfil quality assurance requirements for measured values and are all also included in the documentation of the measurements.

Units come complete with all accessories required for accurate measurements.

Continuous flow-through measurements are possible up to a pressure of 4 bar with FLOW-THRU-TURB vessel.



- Measuring range 0.0001 to 10000 NTU with AutoRange function
- Automatic 1...5 point calibration
- Values displayed in
 - NTU
 - EBC
 - FNU, FAU (Turb 555 IR)
 - Nephelos (Turb 550)
- Flow-through measurement



Ordering Information

Model		Order No.
Turb 550	Laboratory turbidity meter with universal power supply 90 ... 250 V, 3 calibration standards 0.02 – 10.0 – 1000 NTU, 2 empty cuvettes	600 100
Turb 550 IR	Laboratory turbidity meter for measurements according to DIN EN 27 027, ISO 7027 (EN ISO 7027) universal power supply 90 ... 250 V, 3 calibration standards 0.02 – 10.0 – 1000 NTU, 2 empty cuvettes	600 110
Turb 555	High-end laboratory turbidity meter according to US EPA with universal power supply 90 ... 250 V, 4 calibration standards 0.02 – 10.0 – 100 – 1750 NTU, 3 empty cuvettes	600 200
Turb 555 IR	High-end laboratory turbidity meter according to DIN/ISO (EN ISO 7027) with universal power supply 90 ... 250 V, 4 calibration standards 0.02 – 10.0 – 100 – 1750 NTU, 3 empty cuvettes	600 210
Flow-through vessels, calibration standards and other accessories see brochure "Product Details"		



Turbidity Portable Turbidity Meters

Turb 430 IR / Turb 430 T

- Scattered light characteristics according to Pharmacopoeia 5.0
- Multifunctional LabStation
- GLP/AQA conform documentation



NEW



NEW: A turbidity measuring lab in the field – the new Turb 430IR/T sets



Lab accuracy & comfort in portable field instrument

With the new turbidimeters **Turb 430 T** and **Turb 430 IR**, the user now has the choice to perform nephelometric measurements at 90° scattered light according to the application and standard required.

The **Turb 430 IR** meets the DIN 27027/ISO7027 requirements, the **Turb 430 T** those of US EPA 180.1. The measuring range is from 0 to 1100 NTU/FNU and is identified automatically. Accurate measurements in the lower range, e.g. in drinking water are no problem!

Menu driven calibration with up to 3 points and all measurement function are easy for even the most inexperienced operator to perform accurate and precise measurements. The calibration is via an AMCO® standards set (0.02-10-1000 NTU). Up to 1000 data sets with ID nos. can be stored and subsequently output via the LabStation in combination with the powerful LSdata software. The quality of the measurement results is supported by adjustable calibration intervals with documentation. (Order no. 251301. see p. 3).

The Turb 430 is not only a field measuring instrument (especially with the practical field case), but also a “**small lab instrument**” for applications up to 1100 NTU/FNU and with optimum data management.

Optional: Field case and rechargeable battery set (see p. 4 and brochure “Product Details” p. 71).



Portable Turbidity Meters

Turb 355 T / Turb 355 IR

Small portable turbidimeter for control purposes

Battery-operated portable turbidity meter with Tungsten lamp according to US EPA or infrared LED (860 nm) for nephelometric measurements according to ISO 7027 / DIN/ EN 27 027 (EN ISO 7027): Handy, lightweight and easy to operate.

The Turb 355 T / IR comes in a handy carrying case. All necessary accessories (calibration standards 0,02 – 10,0 and 1000 NTU, empty cuvettes and batteries) are included. The instrument is powered by 4 AAA batteries.

0 - 1100 NTU/FNU

Easy operation



IP 67



2 Year Warranty

Technical Data

	Turb 430 IR / Turb 430 T	Turb 355 T / 355 IR
Measuring principles	Nephelometric (90° scatter)	Nephelometric (90° scatter)
Light source	IR-LED/Tungsten lamp	Tungsten lamp/IR-LED
Measuring range NTU FNU	0 ... 1100 / 0-1100 0 ... 1100	0 ... 1100 0 ... 1100
Resolution	0.01 from 0.00 ... 9.99 0.1 from 10 ... 99.90 1 from 100 ... 1100	0.01 NTU from 1 ... 9.99 0.1 NTU from 10.0 ... 99.9 1 NTU from 100 ... 1000
Accuracy	0.01 NTU or ±2 % of the measured value	±2% of the measured value or ±0,1 NTU last decimal place in range 1 ... 500 NTU ±3 % of the measured value in range 500 ... 1100 NTU
Reproducibility	<0.5% of the measured value or 0,01 NTU/FNU	±1% of the measured value or ±0,05 NTU/FNU
Calibration	Automatic 3 point calibration	Automatic 1...3 point calibration
Response time	Approx. 3 seconds (IR) / approx. 7 seconds (T)	14 seconds
Cuvettes	1.10x2.36 in. (28x60 mm), 20 ml sample volume	0.98x1.77 in (25x45 mm), 15 ml sample volume
Interface	RS 232, USB via adapter	
Special functions		
Calibration protocol	Yes	—
Storage of measured value	1000	—
RS 232	Yes	—
Date/Time	Yes	—
Data evaluation	Yes	—
Rechargeable battery	Optional	—
Operating temp.	32 ... 122 °F (0 ... +50 °C)	32 ... 122 °F (0 ... +50 °C)
Power supply	4 x AA batteries for approx 3,000 measurements	4 micro (AAA) alkaline manganese batteries suitable for more than 1,500 measurements

Ordering Information

Model		Order No.
Turb 355 IR	Portable turbidity meter according to ISO 7027 / DIN EN 27 027 (EN ISO 7027) in professional case with 4 micro (AAA) alkaline manganese batteries, 3 calibration standards 0.02 – 10.0 – 1000 NTU and 2 empty cuvettes	600 311
Turb 355 T	same as Turb 355 IR, but with tungsten lamp according to US EPA	600 312
Turb 430 IR	Portable turbidity measuring instrument for nephelometric measurements (90°) according to DIN EN 27027, incl. calibration kit (0.02 - 10 - 1000), 2 empty cuvettes, cleaning tissues, batteries (4 x AA), suited for drinking water. (Optional LabStation or rechargeable battery pack, see brochure "Product Details" p. 71)	600 320
Turb 430 T	Portable turbidimeter for nephelometric measurement (90°, tungsten) acc. US EPA 180.1, incl. calibration standard kit (0.02-10-1000 NTU) and accessories: 2 empty cuvettes (28 mm), cleaning tissues, batteries (4 x AA); suitable for drinking water. Optional: LabStation or rech.batt. pack or as set (see brochure "Product Details" p.71)	600 325

For information visit www.WTW.com for a customer care center near you or inside US: call WTW 800 645 5999.

Parameter

pH

ORP

ISE

Oxygen (D.O.)

Conductivity

Multi-parameter

BOD/Respiration

Photometers

Turbidity

Colony Counter

Software/Printers



Colony Counter

Colony Counting

BZG 30



- Intelligent counting technology
- Operator Comfort
- Safe
- Flexible

One and two and three...

The most time-consuming part of carrying out a microbial count is the actual counting process on Petri dishes. Colony counters facilitate this task and as a result are essential in every bacteriological laboratory. Their particular advantages are the easy, rapid and reliable counting of bacterial colonies with very sample handling.

The **colony counter BZG 30** is easy to use, and allows fatigue-free and safe work thanks to its problem-free counting mechanism: The surface on which the Petri dish is placed is pressure-sensitive; marking the colonies on the Petri dish with the marker pen activates the counter. The pressure sensitivity is adjustable for optimal adaptation to the operator.



Colony Counter

Intelligent counting technology

The colony counter **BZG 30** has an acoustic check signal and automatically compensates for the weights of different Petri dishes; the pressure sensor system also provides uniform sensitivity over the whole of the working field. Zero-voltage protection guards against power interruptions.

The ring-shaped lamp provides uniform working field illumination; the magnifier offers distortion-free optics and can be individually positioned on its extremely flexible arm.

Versatile work is guaranteed by selectable direct or indirect illumination. Direct illumination is particularly suitable for nutrient agars with a dark background; indirect illumination for lighter agars. Colony counter **BZG 30** has an additional counter input for counter marker **KS 30** and is supplied with an interchangeable Wolffhügel grid.

Technical Data

Counter	LED display (0 ... 999) with reset to 0
Standard magnifier	1.7-fold surface magnification, 100 mm dia.
Illumination	20 W ring lamp
Power supply	230 V ± 15%, 50 ... 60 Hz or 115 V ± 15%, 50 ... 60 Hz, as selected
Power consumption	21 W
Dimensions	300 x 325 x 90 mm (B x D x H) with magnifier arm
Weight	5.6 kg
Test certificate	CE
Warranty	1 year

Ordering Information

Colony Counter		Order No.
BZG 30	Colony counter with magnifier (1.7-fold, 100 mm dia.) complete with plug-in flexible arm, switchable background (light - dark) and Wolffhügel grid; for Petri dishes 70 mm and 100 mm dia. 230 V / 50 ... 60 Hz	803 314
BZG 30	Ditto, but for 115 V / 50/60 Hz	803 315
BZG/LUP 8	Surcharge for colony counter with special magnifier (2.3-fold magnification, 60 mm dia.) and plug-in flexible arm (replaces 1.7-fold magnifier)	803 319



Software Printers

MultiLab[®] pilot communication Software

WTW software – simple and comfortable

MultiLab[®] pilot, the communication program specially created for inoLab[®] 740 and 750. For these instruments MultiLab[®] pilot is not just limited to comfortable on-line data transfer, but can also control the active multifunction box or the inoLab[®] and 750.

MultiLab[®] pilot can also be used to comfortable data transfer with the inoLab[®] 730, handheld meters MultiLine P3, P4 and 340/340i and other instruments.

Digital on-line recorder

Export function for Word, Excel, Access and other programs

Free-of-charge download from the internet



New: Adapter for USB to serial interfaces

Minimum system requirements

Hardware:

- IBM-compatible PC with 486DX processor
- 8 MB RAM
- 20 MB free memory
- VGA/SVGA graphics board
- CD-ROM drive
- one free interface (RS 232) for each connected instrument

Software:

- Windows 95

MultiLab[®] pilot, the free-of-charge software package for all laboratory, handheld and field instruments, can be downloaded from the internet

“MultiLab[®] pilot” controls the measurement and carries out data transfer and storage. The data can be shown online as a table or graph or off-line transferred to a printer. The inoLab[®] 740 and 750 can be controlled active.


Well laid out



All display modules are arranged as independent windows which can be moved about as required. Each window has its own menu bar with which settings for the measurement can be made quickly and conveniently. Additional graphic symbols, on-line help texts and a large measurement display round out the picture.

Graphic

A comfortable digital recorder is used as a graphic support for data monitoring in real-time mode. Apart from zoom and scrolling functions, individual values can be displayed or limiting values laid down via the cursor.

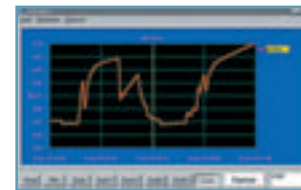

Interface cables for connecting a PC or an external printer (P3001)

Instrument	Model Suitable accessories	Instrument	Model Suitable accessories
Cond 197i	b ④,⑤	Oxi 340	b ④,⑤
Cond 340i	b ④,⑤	Oxi 340i	b ④,⑤
inoLab [®] 730	b ④,⑤	Oxi 538	u ④,⑤
inoLab [®] 735	b ④,⑤	Oxi 597	u ④,⑤
inoLab [®] 740	f ④,⑤,⑥,⑦	pH 197i	b ④,⑤
inoLab [®] 750	f ⑦,⑧,⑨	pH 197	u ④,⑤
inoLab [®] Level 2	b ④,⑤	pH 325	u ④,⑤
inoLab [®] Level 3	f ④,⑤,⑥,⑦	pH 340	b ④,⑤
LF 197	u ④,⑤	pH 340i	b ④,⑤
LF 325	u ④,⑤	pH 340/ION	b ④,⑤
LF 340	b ④,⑤	pH 538	u ④,⑤
LF 538	u ④,⑤	pH/Cond 340i	b ④,⑤
LF 597	u ④,⑤	pH/ION 340i	b ④,⑤
Multi 197i	b ④,⑤	pH/Oxi 340i	b ④,⑤
MultiLine [®] P3 pH/LF	b ④,⑤	photoLab [®] Series	b ⑥
MultiLine [®] P3 pH/Oxi	b ④,⑤	pHotoFlex Series	u ⑥
MultiLine [®] P4	b ④,⑤		
Multi 340i	b ④,⑤	④ = AK Labor	④ = AK 540 B
Multi 350i	b ④,⑤	⑤ = AK 340/B	⑤ = AK 325/S
Oxi 197i	b ④,⑤	⑥ = M-PC/5	⑥ = AK T-PC
Oxi 197	u ④,⑤	⑦ = AK T-P9 PIN/25 PIN	⑦ = AK T-R 2ST
Oxi 325	u ④,⑤	b = bidirectional	f = can be remote controlled
		u = unidirectional	



Communication Program ACHAT OC for BOD Measurement with OxiTop® Control

Stored measuring data can be transferred simple and comfortable from the OxiTop® Control system to a PC with the communication program ACHAT OC. These transferred data can be processed by PC or the Controller OC 100/OC 110.

Min. software requirements:


Windows 3.1 update options on request.

Necessary accessories:

Connection cable AK 540/B.

PC Software Multi/ACHAT II:

PC Communication with WTW Photometers photoLab®

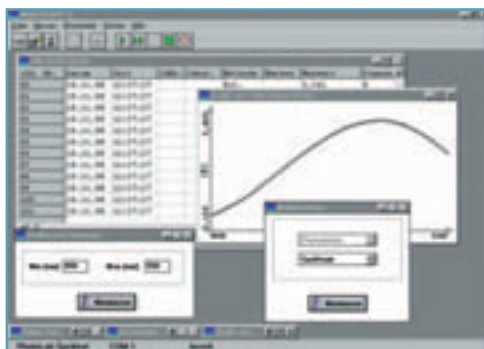

- dBASE format
- Sorting according to measurement parameter
- Windows program with context-sensitive menus and help programs

WTW offers the **Multi/ACHAT II** PC software program so that stored measuring data can be transferred simply and comfortably from the instrument to a PC for documentation and statistical evaluation.

PC communication with pHotoFlex photometers: LSdata see page 88

Ordering Information

	MultiLab® pilot	Order No.
KOM pilot	Communications package, consisting of 1 x MultiLab® pilot and 1 AK 340/B	902 915
AK M-PC	Adapter cable for connecting a PC to the multifunction box, incl. MultiLab® pilot	902 901
AK T-PC	Interface cable for connecting the inoLab® 740/Level 3 terminal to the PC	902 902
ADA USB/Ser	Adapter USB serial interface RS 232 (9-pin socket)	902 880
	Multi/ACHAT II	Order No.
Multi/ACHAT II	Software for PC under Windows, German and English	902 750
KOM Labor	Communications package, consisting of: 1 x Multi/ACHAT II and 1 AK Labor	902 754
AK Labor	Interface cable for connecting instrument to a PC via the RS 232 interface including 9/25-pin adapter and operating instructions	902 758
	ACHAT OC	Order No.
ACHAT OC	PC communication program for Controller OxiTop® OC 100 or OC 110 for further processing of measuring data	208 990
AK 540/B	Interface cable for connecting instrument to a PC via the RS 232 interface including 9/25-pin adapter and operating instructions	902 842
Additional accessories see brochure "Product Details"		



Software Printers

WTW Printers

WTW instruments with a serial interface can be connected directly to a PC (see software section) and can therefore use the PC printer for their printouts.

For protocol purposes a printer can also be connected directly.

For instruments with a serial interface WTW offers suitable printers and cables. As the transmission rate (baud rate) of most instruments is permanently fixed it may be necessary to set the printer to the suitable transmission rate.

Additional Information
see brochure "Product Details".



P 3001



LQ 300+

Technical Data

Model	P 3001	inoLab® printer	LQ 300+
Printing method	Thermal printer		24-pin matrix printer
Printing line	40 characters/line	-	80 characters at 10 cpi
Papierbreite	112 mm		182 mm ... 216 mm (single sheet, autom. feed)
Papierart	Thermal printer paper, normal quality: legible approx. 5 years high quality: legible at least 10 years		Normal paper 52.3 g/m ² ... 90 g/m ²
Size (W x D x H)	170 x 170 x 66 mm	-	366 x 275 x 141 mm
Weight	Approx. 1 kg	-	Approx. 4,3 kg
Power supply	230 V AC, 50 Hz rechargeable battery	-	220 V AC ... 240 V AC, 50 Hz ... 60 Hz
Ambient temperature - Operation - Storage	0 °C ... 40 °C -20 °C ... 55 °C		+5 °C ... 35 °C -20 °C ... 55 °C
Tested safety to			EN 60 950
Interface(s)	RS 232 (serial) Centronics (parallel)	-	RS 232 (serial) Centronics (parallel)

Ordering Information

Printer P 3001		Order no.
P 3001	Thermal matrix printer, mains and rechargeable battery operation. Paper width 112 mm, preset baud rate: 4800; preset characters/line: 40	250 045
AK 325/S	Interface cable to connect instrument to P 3001	902 837
AK 540/S	Interface cable to connect instrument to P 3001	902 843
Printer LQ 300+		Order no.
LQ 300+	Matrix printer, mains operation on 230 VAC. Standard paper (A4 or endless)	250 046
AK/LQ 300	Interface cable to connect instrument to LQ 300+ (only photoLab® series, not inoLab®, not 3xx(i) series)	250 746
Further connection cable and accessories see brochure "Product Details"		



Customer Service/Certificates

We are doing something...


to solve your quality assurance problems

If you make a measurement you make a mistake. This applies in particular to the calibration of instruments. In order to quantify this error the variation of the measured value from a national or international standard is given. In a transferred sense this is the "standard meter" of the quantity involved.

In chemical analysis reference materials are used. These substances are measured in appropriate metrological facilities. The uncertainty of a measurement for such a material is documented. Examples of facilities which can provide such qualified evidence are the National Institute of Standards (NIST, Gaithersburg, USA) and the Physikalisch-Technische Bundesanstalt (PTB, D-Braunschweig).

In further steps secondary, tertiary, etc. reference materials are prepared from the primary reference materials by comparative measurements. The uncertainty of each of these steps compared to the "standard meter" can be given; this takes into account the measuring agents and methods used. It is important that the calibration of a measuring system can be traced back to the particular standard in an unbroken chain with defined uncertainty.

In practice working reference buffer solutions are used; these are obtained by comparing them with primary or secondary material. WTW pH buffers meet these requirements. The individual uncertainty of the pH of a particular buffer solution is documented by a certificate.

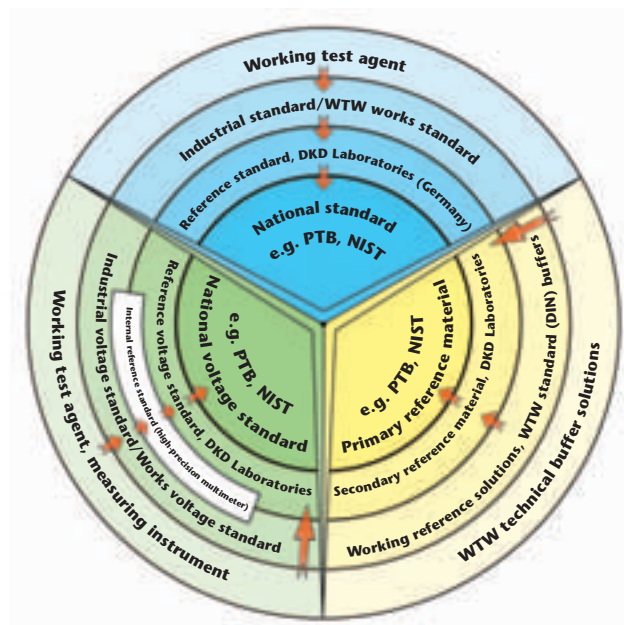



What we can offer you

Calibration of measuring systems for determination of conductivity, pH, and dissolved oxygen in aqueous media

No certification to DIN ISO 9000 without test agent monitoring

The perfect functioning of the test agent used is a constant requirement for the correctness and comparability of measured values. This is why it belongs to the basic rules of **quality assurance** and **Good Laboratory Practice** that the accuracy of each test agent is monitored at regular intervals after a defined period of use by carrying out a calibration. This task is faced by a continually increasing number of companies and laboratories who are trying to achieve or have already achieved certification of their QA system according to the DIN ISO 9000 series of standards.

Why you should make use of the manufacturer's technical competence

Proper calibration requires specially qualified personnel with particular knowledge of the individual instrument and the presence of suitable calibration facilities. This is why it is usually more efficient and economical to allow test

agent monitoring to be carried out by an external calibration laboratory or directly by the manufacturer.

WTW is available as a competent partner and can provide this service for all WTW measuring systems for the determination of conductivity, pH, and dissolved oxygen.

We have been certified to ISO 9001 since 1993 and are completely familiar with the requirements of the standard. Our calibration facilities are linked to national standards. Calibration agents for which no national standards exist are prepared in accordance with recognized national and international standard methods.

We carry out a works calibration and provide you with a calibration certificate.

If required we can also carry out test agent monitoring for our photometers and BOD measuring instruments. Please ask for our advice.

WTW offers different types of certificate

1. Certificate of Compliance

General certificate (without mentioning the serial number) which certifies that the product complies with the technical data given in the operating instructions. The certificate is not signed and is free of charge.

2. Manufacturer's Test Certificate

Individual certificate (mentioning the serial number) which states that the product has been tested and complies with the information given about accuracy on the certificate. Contains a passage about the regular calibration of the test agents used by us and their traceability to national and international standards. Can be produced by the customer as evidence for ISO 900 purposes.

Certificate for brand-new products:

These Certificates are added to all instruments. The certificate is not signed and is free of charge.





Customer Service/Certificates

CE Declaration of Conformity

Certifies that the product complies with the valid EC directives.

Certificates according to FDA regulations

Validation of instruments according to FDA regulations like IQOQPQ on request.

Manufacturer's certificate for calibration solutions

When ordering or within 3 months after purchasing we can supply a manufacturer's certificate for the pH buffer solutions and conductivity calibration solutions offered in our range of products; this certifies their controlled manufacture on the basis of national and international standards.

Calibration certificates for which there is a charge

Calibration certificate for an instrument


The measuring functions of the instrument are calibrated independently of the signal generator by using electrical standards.



Calibration certificate for a signal generator

For pH electrodes and conductivity cells the calibration is made by using calibration solutions. For oxygen sensors the slope is calibrated by using air saturated with water vapor and the zero current by using a zero solution or in pure nitrogen.

With pH electrodes and oxygen sensors a gradual alteration to the technical data occurs (aging). This is why they have to be calibrated by the operator at regular intervals; the procedure is described in the instruction manuals of the corresponding instrument.

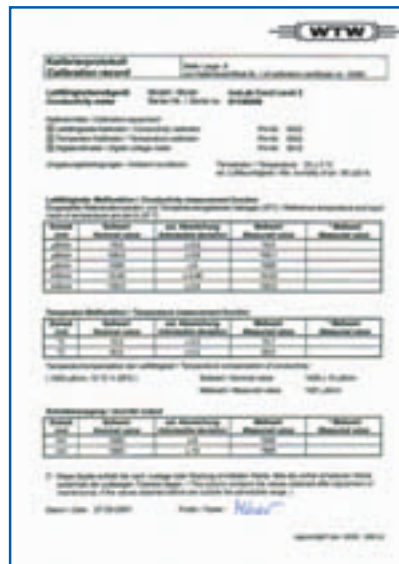

Certificate for used products:

Provided at the customer's request in association with a repair contract. Test data are listed in a protocol. The certificate is signed by our QM officer and will be invoiced.



Analog

Monitors
EcoLine/QuadroLine®
and sensors

- High accuracy and enhanced EMC performance using integrated pre-amplifier
- Integrated lightning protection
- EcoLine 170: Monitors for field installation
- QuadroLine® 296: Monitors for panel mounting 96 x 96 mm

D.O.
pH
Cond
Turb

NH_4

NO_3

NO_2

Analyzer

TresCon®/
TresCon® Uno

- TresCon®: Multi-parameter-analyzer for up to three parameters
- Self-calibrating systems: easy to use – easy to extend
- Also available as a compact single parameter unit



Digital

Multiparameter System

IQ SENSOR NET

System 182 (XT), 184 XT
and 2020 XT

TSS

NH₄

NO₃

COD

TOC

DOC

SAC

BOD

P Total

PO₄

- One system for all parameters
- High accuracy and enhanced EMC performance using – integrated pre-amplifier & digital processing
- Integrated lightning protection
- Sensor can be pre-calibrated in Lab
- Universal sensor connection – Standard for all digital sensors
- Easily expandable using 2-wire technology
- 0/4 ... 20 mA
RS 232, RS 485, PROFIBUS-DP, Modbus RTU



Are you interested?

Please order the new
WTW On-line Catalog



WTW

About us

Company highlights

- 1945** Company founded by **Dr. Karl E. Slevogt**
- 1948** Renamed to *Wissenschaftlich-Technische-Werkstätten (WTW)*
- 1954** Introduction of first WTW pH meter
- 1965** Introduction of first WTW oxygen meter
- 1976** Bavarian State Award for the **Combibox** compact multi-parameter system
- 1982** Introduction of the first zero-current-free (stable zero point) oxygen sensor for field measurements in the world
- 1983** Start of WTW's on-line measuring technology program
- 1986** First company to offer a 3-electrode oxygen sensor (**TriOxmatic®**) with automatic calibration on air (**OxiCal®**)
- 1987** First company to offer a 4-electrode conductivity sensor (**TetraCon®**) for portable water analysis
- 1993** First manufacturer of D.O., pH and conductivity measuring systems to be certified to ISO 9001
- 1995**
 - Introduction of the mercury-free **OxiTop®** system for manometric BOD determination
 - First company to offer monitors with built-in lightning protection
- 1997** New **photoLab®** laboratory photometers combine precision with outstanding ease of use
- 1998**
 - Introduction of the **PurCon®** sample preparation system as a replacement for conventional filtration systems
 - First WTW spectrophotometer
- 1999** "Innovations that make sense": the new laboratory instruments of the **inoLab®** family with their 3 function levels set new standards for the analytical measurement of pH, oxygen, conductivity and temperature as well as for ISE and multi-parameter measurements



Sean Donnelly, CEO WTW GmbH

There are not many companies around that can look back on 60 successful years. WTW is such a company, and one of the secrets of its enduring success is undoubtedly the permanent strive for top performance in the development and production of water analysis instruments. We are driven by our goal to offer complete solutions to make our customers work easier.

"To make good things even better" is one of the most difficult tasks one could imagine, but we believe there are always opportunities for improvement. This will be part of WTW's strategy for the next 60 or more years – take our word for it! Our customers and partners can look forward to new and innovative products in 2006 and beyond.

Laboratory & Field Instrumentation

Offering the world's most complete line of pH/ORP, D.O./BOD/Respirometry and Conductivity Instruments, Turbidity Meters, Photometers with reagents, WTW systems range from rugged waterproof, hand-held, and portable field meters to an integrated line of laboratory instruments, supplies and accessories that include the new and innovative inoLab™ systems and a wide range of technically advanced Multi-parameter instruments that represent the state-of-the-art in measurement systems. Our latest innovation is the 350i Multi-parameter handheld meter. This product delivers pH, Conductivity, Dissolved Oxygen, and temperature measurements in a portable, water-proof system with a compact multi-parameter probe.





About us

On-Line Instrumentation

The IQ SENSOR NET system is a flexible, digital based system providing from 1 up to 20 measuring points suited for – conventional instrumentation with analog outputs as well as for field bus instrumentation. The innovative digital sensors in this system represent the state of the art in process measuring systems. Our latest addition to the digital sensor offering is the NiCaVis™ product. This measurement system measures Nitrate and Carbon (COD), simultaneously in-situ.

The world's most complete line of pH/ORP, D.O., Conductivity and unique self cleaning Turbidity instrumentation. Complete systems, probes and accessories are available for measuring and monitoring water and wastewater.

TresCon™ Analyzers provide unparalleled modular flexibility and convenience for the on-line measurement and monitoring of up to three different parameters and uses the unique PurCon™ sample conditioning system for high accuracy and reproducibility in highly polluted wastewater samples.

The dependability, reliability, and versatility of WTW field proven Ammonia, Phosphate, Nitrite and Nitrate Analyzers, probes, and pH, ORP, D.O., and Conductivity systems and meters have established WTW products as industry standards world-wide.

WTW has built a solid reputation in its 60 year history by providing "best in class" products with unparalleled customer and technical support. WTW strives to deliver solutions to our customers measuring problems. Our Customer Care Centers are dedicated to ensuring each customer's individual success. WTW's extensive applications library coupled with knowledgeable applications specialist provides for rapid resolution to technical challenges.

With support facilities around the globe, the WTW manufacturing center, located just south of Munich, Germany, delivers quality technical instrumentation with continuous support. We are proud to present our product offering to you and look forward to serving your needs.



Company highlights

2000 Introduction of **TresCon®** – the modular analytical system for the continuous measurement of ammonium, nitrite, nitrate, phosphate

2001

- **IQ SENSOR NET** – the multi-parameter measuring system offers unlimited possibilities for on-line measurements
- The new **VisoTurb®** and **ViSolid®** turbidity and solid sensors with their revolutionary ultrasonic keeping clean system give "low-maintenance" a completely new meaning

2002

- **AmmoLyt® 700 IQ** enables reliable On-line "in-situ" measurement of Ammonium
- **PurCon® IS: Sample Preparation** – directly without pump

2003 **NitraLyt 700 IQ** is a perfect supplementary nutrient parameter (Nitrate) for On-line "in-situ" measurement

2004

- Multi-parameter handheld meter **Multi 350i** represents state-of the art in field applications
- **NitraVis®, CarboVis®** und **NiCaVis®** – spectral "in-situ" On-line sensors for combined Nitrate, Carbon and TSS measurement providing new solutions for waste water control

2005



- Portable photometers and turbidity meters for universal applications:
pHotoFlex/pHotoFlex Turb Turb 430 IR
- **IQ SENSOR NET System 182**
compact 2 channel transmitter as a supplement to the IQ SENSOR NET family
- **ECDO DO-Sensors**
Best for DO online measurement

2006 • **VARION** ammonium and nitrate multisensor with automatic compensation of interference ions



WTW

Contact Addresses



**Wissenschaftlich-Technische
Werkstätten GmbH**

Dr.-Karl-Slevogt-Strasse 1

D-82362 Weilheim

Germany

Tel: +49 (0) 881 183-0

+49 (0) 881 183-100

Fax: +49 (0) 881 183-420

E-Mail: Info@WTW.com


Internet: <http://www.WTW.com>

Looking for your local contact?

Here you can find your
"local WTW":

contact addresses, representatives,
distributors...

Click!

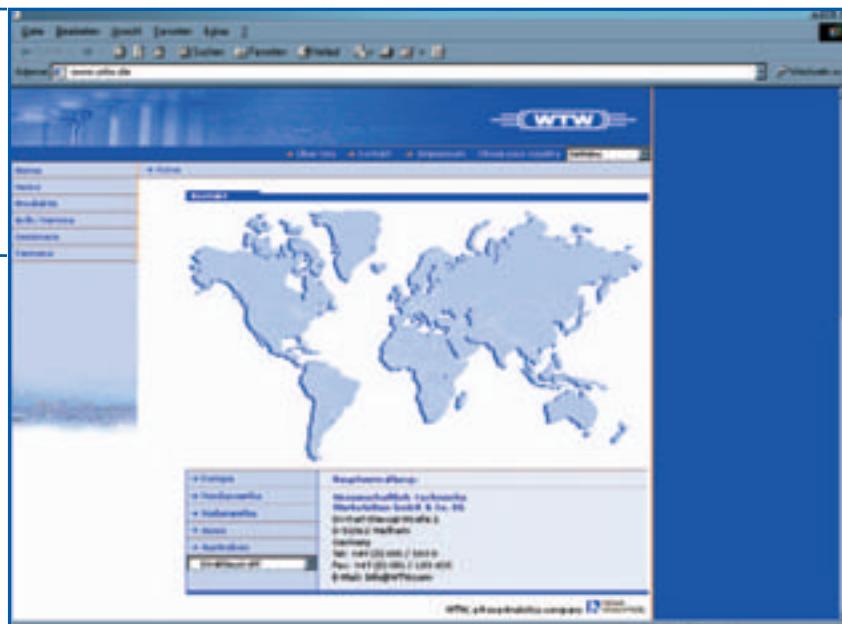




WTW
www.WTW.com

WTW Contact

Information around the Clock!

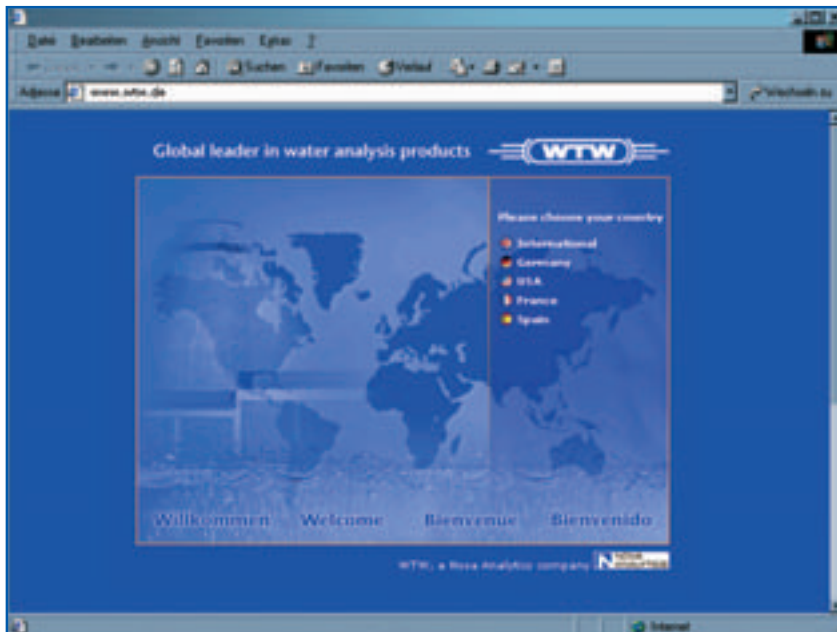

- Up-to-date
- Informative
- Multilingual

New Products

Click in – WTW presents its complete line of new products, innovative measuring and analytical instruments, helpful accessories, useful system extensions, special sets and much more –

24 hours a day..

Applications

WTW can provide you with a solution for all your measurement needs. In addition you will find tech tips, application notes,

and much more



Downloads

You need a Manual, Application Report or a WTW Certificate?

Have a look at our Download Area...



Important information!

General information

1. Special versions of instruments on request.
2. Accessories and spare parts for older models – please make separate inquiry.
3. In order to avoid our customers having to pay a surcharge for small-volume purchases, we supply our consumables in practical minimum ordering quantities.

Technical alterations

The technical description corresponds to the current products. Alterations because of technical improvements are possible.

Illustrations

We draw your attention to the fact that the illustrations are intended to clarify certain points. There may therefore be discrepancies between the illustrations and the written text.

Liability

We accept no responsibility for printing errors, writing errors or mistakes in the translation.

edition January 2006

Publisher



Wissenschaftlich-Technische Werkstätten GmbH

Dr.-Karl-Slevogt-Strasse 1

D-82362 Weilheim

Germany

Tel: +49 (0) 881 183-0
+49 (0) 881 183-100
Fax: +49 (0) 881 183-420
E-Mail: Info@WTW.com
Internet: <http://www.WTW.com>



Index

00594*	93	14695 CombiCheck 50	100
00595	94	14696 CombiCheck 60	100
00597	94	14729	99
00598/1	94	14730	98
00598/2	94	14731	94
00599	94	14732	95
00602/1	94	14738 CombiCheck 80	94, 98
00602/2	94	14739	100
00605	93	14752	96
00606	95	14758	96
00607/1	98	14761/1	94
00607/2	98	14761/2	94
00608	94	14763	95
00609*	96	14764	95
00613	97	14767	97
00614	97	14770	94
00615	95	14773	95
00616	98	14776/1	97
00617	99	14776/2	96
00675	93	14779	96
00680	101	14785	99
00683	96	14791	96
00687	97	14794	99
00718	101	14815	99
00796	95	14821	93
00798	98	14825	93
00815	95	14828	94
00816	95	14831	93
00826	93	14832	99
00856	93	14833	98
00857	99	14834	93
00858	93	14839	98
00860	95	14842	98
00861	99	14848	98
00885	96	14878	99
00961	99	14879	99
01632	95	14895	97
01739	95	14896	95
01744	98	14897	94
01745	93	14942	97
01746	99	14965	93
01747	93	18789*	95
01762/1	93	19500	101
01762/2*	93	19770	101
01763	98	19776	101
01764	99	19777	101
01787	99	19778	101
06146	99	19779	101
09017	101	19780	101
09701	94	19781	101
09711	96	19786	101
09713	97	19789	101
09717	98	19792	101
09772	98	19797	101
09773	98	19806	101
14394	99	19811	101
14500	95	19812	101
14537	97	19813	101
14542	97	19814	101
14543	98	19897	101
14544	96	19898	101
14546	98	19899	101
14548	99	70230	101
14549	95	70236	101
14551	93		
14552	94	A	
14553	94	A 325/S	page
14554	96	A5/25	59
14555	97	ACHAT OC	96
14556	97	ADA USB/Ser	113
14557	94	AK 325/S	113
14559	96	AK 540/B	114
14560	97	AK 540/S	114
14561	94	AK Labor	113
14562	95	AK M-PC	113
14564	99	AK T-PC	113
14566	99	AK/LQ 300	113
14598	94	Al-1 TP	114
14622	99	AS absorption tube	93
14675 CombiCheck 20	100		
14676 CombiCheck 10	100	B	
14677 CombiCheck 30	100	BOD nutrient salt 00688	page
14678	95	BOD SFR oxygen reaction bottles 14663	97
14683	99	BOD standard	97
14689 CombiCheck 70	100	BOD thermostat boxes	78
14690	97	BOD thermostat cabinets	79
14691	97	BSB/BOD 740P	65
14692 CombiCheck 40	100	BZG 30	111
14694	98	BZG/LUP 8	111

Index

100	C	page
100	C1/25	97
99	C2/25	97
98	Calibration agents	48
94	CellOx® 325	37
95	Cl-1 TP	94
94, 98	Cl-2 TP	94
100	Cl-3 TP	94
96	COD1 TC (LR)	97
96	COD2 TC (MR)	97
94	COD3 TC (HR)	97
95	Colony Counter	111
95	Cond 197i	42
97	Cond 315i	44
97	Cond 330i	44
94	Cond 340i	44
95	Cond 720	41
97	Cond 730	41
96	Cond 740P	41
96	Conductivity flow-through cell	47
99	ConOx-3	59
96	CR 2200	91
99	CR 3200	91
99	CR 4200	91
93	Cu-1 TP	94
93		
93	D	page
94	D 01/T	49
93	D 1/T	49
99	D 201	49
98	D 530	49
93	DurOx® 325-3	37
93		
98	E	page
98	E/SET	48
99	ELY/ORP/AG	23
99	EP/SET	48
97		
95	F	page
94	FC pHotoFlex/Turb 430	89
97	Fe-1 TP	95
93	Fe-2 TP	95
95		
101	H	page
101	Handheld Conductivity Meter SETs	44
101	Handheld Ionmeter	29
101	Handheld Multi-Parameter SETs	59, 61
101	Handheld Oxygen Meter SETs	36
101	Handheld pH Meter SETs	17
101		
101	I	page
101	inoLab® BSB/BOD 740P	65
101	inoLab® Cond 720	41
101	inoLab® Cond 730	41
101	inoLab® Cond 740P	41
101	inoLab® Laboratory Conductivity Meter SETs	41
101	inoLab® Laboratory Ionmeters	28
101	inoLab® Laboratory Oxygen Meter SETs	33
101	inoLab® Laboratory pH Meter SETs	15
101	inoLab® Multi-Parameter SETs	53, 55
101	inoLab® Multi 720	53
101	inoLab® Multi 740	53
101	inoLab® Oxi 730P	33
101	inoLab® Oxi 740	33
101	inoLab® pH 720	15
101	inoLab® pH 730	15
101	inoLab® pH 740P	15
page	inoLab® pH/ION/Cond 750	28, 55
59		
96		
113	K	page
113	KCSB 100	101
114	KCSB 400	101
113	Kit for measuring the conductivity according to USP 28	48
113	KOM Labor	113
113	KOM pilot	113
113	KS 100µS	48
114	KS 5µS	48
93		
93	L	page
page	Laboratory Conductivity Meter SETs	41
97	Laboratory Ionmeters	28
97	Laboratory Oxygen Meter SETs	33
97	Laboratory pH Meter SETs	15
78	LQ 300+	114
79	LR 325/001	47
65	LR 325/01	47
111	LR01 V	47
111	LS Flex/430	89



Index

M

Mn-1 TP	page 95
Mo-1 TP	95
MPP 350-3	59
Multi 197i	56
Multi 340i	61
Multi 350i SET	59
Multi 720	53
Multi 740	53
Multi/ACHAT II	113
MultiLab® pilot	113

N

N1/25	page 97
N4/25	96
NH4-1 TP	96
NH4-2 TC (LR)	96
NH4-3 TC (HR)	96
NO2-1 TP	96
NO2-2 TC	96
NO3-1 TC	97
Ntot1 TC (LR)	97
Ntot2 TC (HR)	97

O

Oxi 197i	page 34, 65
Oxi 315i	36
Oxi 330i	36
Oxi 340i	36
Oxi 730P	33
Oxi 740	33
OxiTop® Box	78
OxiTop® Control 12	71
OxiTop® Control 6	71
OxiTop® Control A12	76, 77
OxiTop® Control A6	76, 77
OxiTop® Control AN12	77
OxiTop® Control AN6	77
OxiTop® Control B6	75
OxiTop® Control BM6	75
OxiTop® Control S12	76
OxiTop® Control S6	76
OxiTop® Control S6/S12	71
OxiTop® IS 12	68
OxiTop® IS 12-6	68
OxiTop® IS 6	68
Oxygen sensors	37

P

P 3001	page 114
P4/25	98
P5/25	98
Passive multifunction box	15, 41, 53, 55
pH 197i	16
pH 315i	17
pH 330i	17
pH 340i	17
pH 720	15
pH 730	15
pH 740P	15
pH/Cond 340i	61
pH/ION 340i	29
pH/ION 735P	28
pH/ION 740P	28
pH/ION/Cond 750	28, 55
pH/Oxi 340i	61
pHotoFlex	89
pHotoFlex/SET	89
pHotoFlex Turb	89
pHotoFlex Turb/SET	89
photoLab® S12	85
photoLab® S12-A	85
photoLab® S6	85
photoLab® S6-A	85
photoLab® Spektral	85
PO4-1 TP	98
PO4-2 TC	98
PO4-3 TC	98
Portable pH Meter	16
Portable Conductivity Meter	42
Portable Multi-Parameter Instrument	56
Portable Oxygen Meter	34
Printer LQ 300+	114
Printer P 3001	114
ProfiLine Cond 197i	42
ProfiLine Multi 197i	56
ProfiLine Oxi 197i	34, 65
ProfiLine pH 197i	16

R

RB Flex/430	page 89
RH 28	23

S

SensoLyt® MPP-A	page 59
SensoLyt® MPP-A Pt	59
SenTix® 20	19
SenTix® 21	19
SenTix® 21-3	19
SenTix® 22	19
SenTix® 41	19
SenTix® 41-3	19
SenTix® 42	19
SenTix® 51	19
SenTix® 52	19
SenTix® 60	19
SenTix® 61	19
SenTix® 62	19
SenTix® 81	19
SenTix® 82	19
SenTix® 91	19
SenTix® 92	19
SenTix® Ag	23
SenTix® Au	23
SenTix® B	19
SenTix® FET-B	20
SenTix® FET-D	20
SenTix® H	20
SenTix® HW	20
SenTix® HWS	19
SenTix® L	19
SenTix® Mic	20
SenTix® Mic-B	19
SenTix® Mic-D	19
SenTix® pH	19
SenTix® PtR	23
SenTix® R	19
SenTix® RJS	19
SenTix® SP	20
SenTix® Sur	20
SenTix® V	20
SenTix® ORP	23
Si-1 TP (LR)	99
Si-2 TP (HR)	99
SK-325	59
SO4-1 TP	99
SORT/RH	23
Special conductivity cell	47
Standard conductivity cell	47
StirrOx® G	37, 65

T

Testing and maintenance supplies for	page
ORP measurements	23
TetraCon® 325	47
TetraCon® 325/S	47
TetraCon® DU/T	47
TetraCon® V	47
TOC Srew Caps	99
TOC Standard Solution 1000 mg/l	99
Trace conductivity cell	47
TS 1006-i	79
TS 606/2-i	79
TS 606/3-i	79
TS 606/4-i	79
Turb 355 IR	109
Turb 355 T	109
Turb 430 IR	109
Turb 430 T	109
Turb 550	107
Turb 550 IR	107
Turb 555	107
Turb 555 IR	107

U

Ultrapure water conductivity cell	page 47
Universal wide-range power supply 100 V-240 V	
50-60 Hz; for 340i series	17, 29, 36, 44, 61
USP Kit 1	48
USP Kit 2	48

V

VARIO	page 18
VARIO Cond	45
VARIO Cond Set	45
VARIO Set	18