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Plankton Screen

Field of application

The AquaCare plankton screen is made for filtering algae or zooplankton cultures. Other applications should be discussed with AquaCare before.

Safety instructions

Never filter sharp particles (e.g. glass dust).
Never clean with sharp tools.
Never put heavy things onto the screen.
Never use oxidizing reagents.

Handling

Pore the suspension carefully and even through the screen. Screen with pore sizes smaller than 50 µm should be filled with more water. Alternatively you may stroke with your hand at the bottom side. The water will flow due to the capillary forces.

You may wash out the filtered plankton with a smooth water jet.

Maintenance

Clean the screen directly after use by washing with clean water (R.O. water oder deionized wa-

ter) in opposite direction. If only tap water is available dry the screen after cleaning with a clean cloths. Otherwise lime crusts will block the screen. – For other cleaning methods ask AquaCare.

Store the plankton screen dark and dust-free.

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Material of body:PMMA (Sperpex)
Material of screen:PA (Polyamide)
Diameter: 200 mm
Pore size: 6...200 µm (see side of screen)
Open filter area: 5...54% (see side of screen)

Warranty

You have 24 months warranty on all AquaCare units excepts spare parts like pump bearings and rotors. You have no warranty if parts are broken by violent.

AquaCare is not responsible for consequential damages.

www.aquacare.de